MINISPENSOR

Mikroliter-Dispenser
Microliter Dispenser

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Content

1 Safety- and user instructions	11
1.1 General safety instructions	11
1.2 Before operating the dispenser	12
2 General product description	13
3 Functional characteristics	14
3.1 Dispensing	14
3.2 Function and ventilation	14
3.3 Volume adjustment	14
4. Cleaning and maintenance	15
4.1 Preparation	15
4.2 Disassembly	15
4.3 Cleaning and maintenance process	16
4.4 Sterilization	16
6 Spare part list	17

1 Safety- and user instructions

1.1 General safety instructions

Highest safety precautions should be used when dispensing corrosive, poisonous, radioactive or hazardous chemicals.

- Observe general safety regulations (e.g. wear safety goggles, protective clothing, etc).
- Carefully follow the operation manual.
- Use the instrument only with regard to resistance of material and for the purpose for which it is intended. In case of doubt concerning the suitability please contact the manufacturer.
- Always check the dispenser for leaks. Make sure that the suction and ejection cannula are firmly in position before starting to use the instrument.
- NEVER use force! Breakage of any part will lead to hazardous exposition to the user/ operator as well as others.
- We only consider ourselves responsible in regard to safety, reliability and performance of the MINISPENSOR if:
 - The dispenser has been used in accordance to the instruction manual
 - Only the parts supplied by the manufacturer have been used

MINISPENSOR

1.2 Before operating the dispenser

Check if dispenser has not been damaged during in transit.

The Minispensor is suitable for nearly all liquids and dispensing methods.

EXCEPTIONS:

- Solutions of hydrofluoric acid or of substances which form crystals, contain or form solids, as well as solutions of substances which are catalyzed by platinum and iridium alloy. Inorganic oxidizable solutions (e.g. biuret reagent) may precipitate metallic oxides, therefore clean the dispenser daily.
- The temperature of the instrument and the reagent should not exceed 40 °C (104 °F).
- The user has to ensure the compatibility of the dispenser and its chemical resistance for himself.

Select the appropriate length of the cannula according to the bottle which will be used and plug it on tightly. Screw the ejection cannula on firmly.

Do not use the plunger before the dispenser has been assembled completely!

The Minispensor fits to the corresponding 100 ml bottle, also supplied with an adapter A28/32; for more adapters, see spare part list

Warning: Avoid personal injury from chemicals. To protect yourself always wear eye protection and other appropriate safety equipment and clothing. Please follow all safety instructions and observe the operating procedures.

Safe and proper operation and function is only guaranteed with the ejection cannula as originally supplied. Never use damaged or deformed cannulas. Do not use any other type of tubing. Inspect the tubes for leakage regularly and replace immediately damaged or deformed ones.

2 General product description

The **MINISPENSOR** is ideal for long pipetting series where you need a fixed micro-volume. In contrast to fixed volume microliter pipettes, there are no pipette tips required. There are four color-coded models available, with two definite volumes in a range of 50 μl to 2 000 μl . You can simply change the dispensing volume to 50% by turning the push button.

Dispensing system

The direct displacement system of the **MINISPENSOR** is for universal use and ideal for crystalline solutions.

Volume setting

Every **MINISPENSOR** is able to dispense two definite volumes. They are easy to set: Just push and turn the button at 180°.

One handed dispensing

The **MINISPENSOR** can easily be used with just one hand.

3 Functional characteristics

3.1 Dispensing

Which fluid do you wish to dispense?

Be sure to observe all safety instructions and operating limitations.

Always place a collecting vessel under the ejection tube. The ejection cannula must face away from operator or any other persons at all times.

Warning: For your own personal safety:

Never press down the plunger as long as there is no vessel to collect the liquid.

CAUTION: If the valve is damaged, reagent may drop out of the ejection cannula!

3.2 Function and ventilation

In contrast to other dispensers, Minispensor is the only dispenser which enables a effortless use. Long series of work can easily be done. This new developed dispenser can be operated with only three fingers. The plunger has to be pushed down with the thumb (just like with an automatic pipette), the selected volume will then be obtained through a spring.

Before the first use the dispenser has to be vented. The air caught in the supply pumps can be removed with 2-3 pumps.

3.3 Volume adjustment

The Minispensor is supplied with a volume adapter, which is divided in 2 fix-volumina.

By rotating the pressure knob you reach a second volume raster, which shows 50% of the total lift.

The two different volumina are named on the volume adapter and the arrow on the plunger, must be lined up with the 2nd arrow, showing on the volume adapter.

4 Cleaning and maintenance

4.1 Preparation

1. Place the complete dispenser (with bottle) in a sink.
2. Unscrew the device from the bottle so that the suction tube will not be immersed in the liquid.
3. Tap the tube carefully against the inner side of the bottle. This allows the remaining reagent to disperse.
4. Remove dispenser from the bottle and rinse with distilled water.
5. It is now possible to disassemble the dispenser. See instructions “Disassembling”.

Warning: Avoid personal injury from chemicals. The instruments, suction and ejection tube may contain reagent. Always face them away from your body. Wear eye protection and protective clothing!

4.2 Disassembly

CAUTION! The dispenser must be completely empty.

1. Screw of the volume adapter (1) of the valve head (2).
2. You now have the pressure knob (function (3)) and volume adapter with the spring (4) in your hand. The PTFE-plunger is fixed to the function button (5).
3. The glass cylinder (6) is attached into the valve head, which you can easily clean now.
4. For assembly just add all components back together again in reverse order.

Caution! The upper part must be screwed on the valve head, otherwise it can leak.

Warning: Avoid personal injury from chemicals. The instruments, suction and ejection tube may contain reagent. Always face them away from your body. Wear eye protection and protective clothing!

Please note:

Never press the plunger downwards with force. Never raise or depress it too fast. Always leave the plunger down and empty after use.

MINISPENSOR

4.3 Cleaning and maintenance process

As you can see on the sectional drawing, MINISPENSOR can be disassembled in seconds. Therefore the dispenser is optimal to clean and, if necessary, it is easy to change spare parts.

To ensure a proper performance and a long life-span, the dispenser should be cleaned in the following cases:

- When changing reagents
- Before storing it over a long period of time
- Daily after you have dispensed:
 - Substances which form crystals
 - Alkaline solutions
 - Organic solutions
 - anorganic oxidizable solutions (fe. Biuret-reagent).

Prevention of clogs of valves during prolonged intervals

When the dispenser is screwed onto the bottle make sure that the valves are surrounded by liquid. Once unscrewed you can maintain an easy running of the valves, if you rinse it with distilled water and/or laboratory cleanser. Afterwards, rinse the dispenser with alcohol.

In order to ensure accuracy, the device should be examined with double-distilled water on semi-micro balance every once a while. Pay attention to the temperature dependence (1 ml water 20° C = 0,998 g)!


4.4 Sterilization

After removal of the suction tube and ejection cannula the dispenser can be steam-sterilized at 121 °C and maximum 2 bar. Place the instrument in an upright position on a cloth. Avoid any contact with hot metal surfaces.

Caution: Slowly cool down the device. After sterilization, let the device cool down to room temperature before using it again. Check all parts and tubing for leakage after every sterilization. In case of deformation return the Minispensor

It is also possible to sterilize the dispenser chemically with alcohol, formaldehyde etc.

5 Spare part list

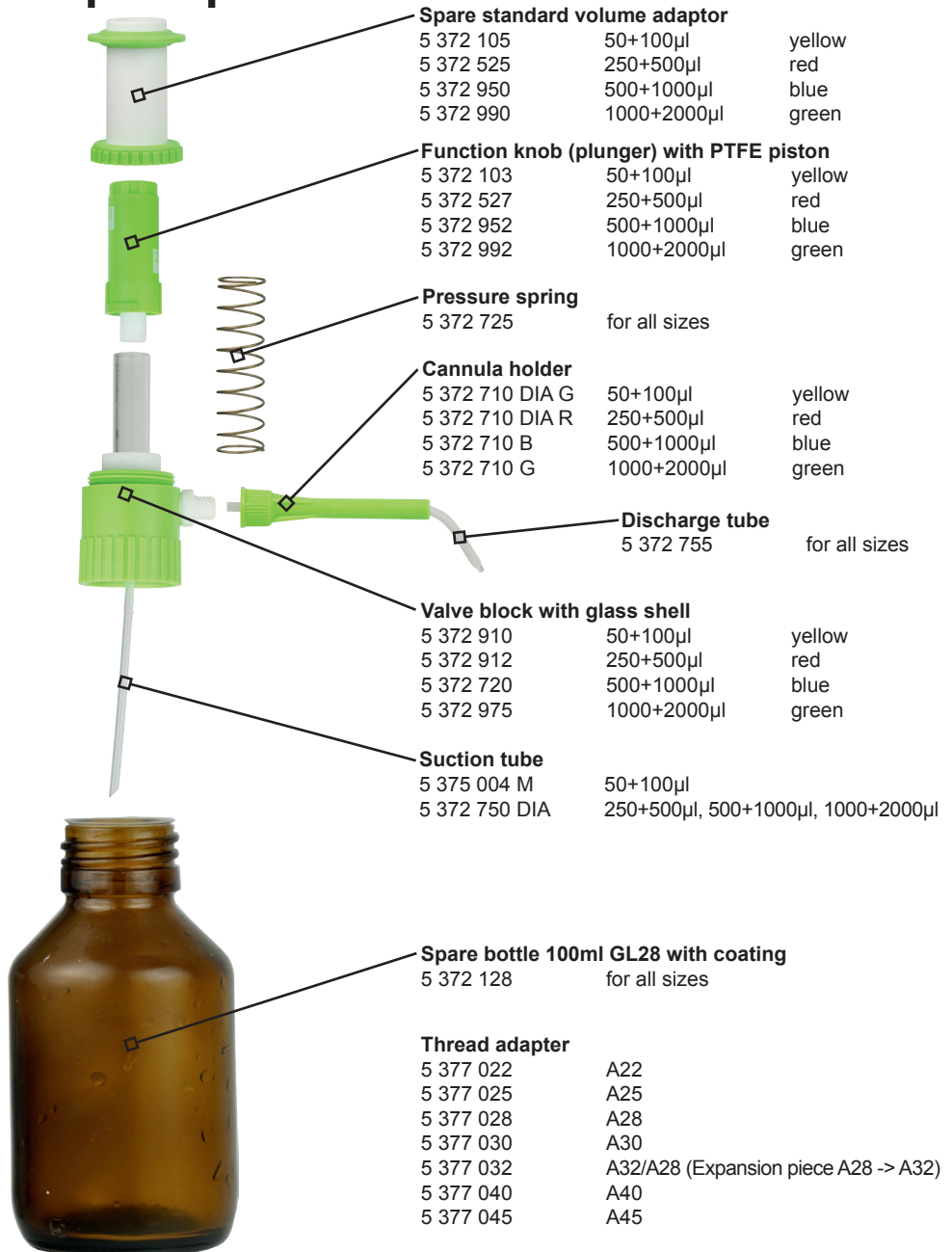Spare standard volume adaptor

5 372 105	50+100µl	yellow
5 372 525	250+500µl	red
5 372 950	500+1000µl	blue
5 372 990	1000+2000µl	green

Function knob (plunger) with PTFE piston

5 372 103	50+100µl	yellow
5 372 527	250+500µl	red
5 372 952	500+1000µl	blue
5 372 992	1000+2000µl	green

Pressure spring

5 372 725	for all sizes
-----------	---------------

Cannula holder

5 372 710 DIA G	50+100µl	yellow
5 372 710 DIA R	250+500µl	red
5 372 710 B	500+1000µl	blue
5 372 710 G	1000+2000µl	green

Discharge tube

5 372 755	for all sizes
-----------	---------------

Valve block with glass shell

5 372 910	50+100µl	yellow
5 372 912	250+500µl	red
5 372 720	500+1000µl	blue
5 372 975	1000+2000µl	green

Suction tube

5 375 004 M	50+100µl
5 372 750 DIA	250+500µl, 500+1000µl, 1000+2000µl

Spare bottle 100ml GL28 with coating

5 372 128	for all sizes
-----------	---------------

Thread adapter

5 377 022	A22
5 377 025	A25
5 377 028	A28
5 377 030	A30
5 377 032	A32/A28 (Expansion piece A28 -> A32)
5 377 040	A40
5 377 045	A45

Notizen / Notes

Notizen / Notes

Liquid Handling

MINISPENSOR

witeg Labortechnik GmbH
Am Bildacker 16
97877 Wertheim
Germany

Liquid Handling

DE-M

www.witeg.de

Phone.: +49 9342 / 9301-0

Fax: +49 9342 / 9301-77

www.witeg.de

info@witeg.de

5 371 XXX_2016-06-16