


Assembly


1. Locate Stand in a place you want to install. <Figure 3>
2. Join Rod in the stand. <Figure 4>
3. Combine Clamp and turn it to clockwise to clamp. <Figure 5>
4. Insert support rod into near Stirrer and fasten by L-wrench. <Figure 6>
5. Put Support bar on Clamp and fasten Clamp's handle to clockwise direction. <Figure 7>
6. Insert Impeller into the bottom of Drill Chuck. <Figure 7>
7. Impeller Length: up to 3~4cm from top of unit, if not the Impeller will shaking when the unit is stirring. <Figure 8>
8. After that, fasten Drill Chuck Key such as Figure 8 <Figure 8>
9. Completed. <Figure 9>




< Figure 3> Stand (ST120)

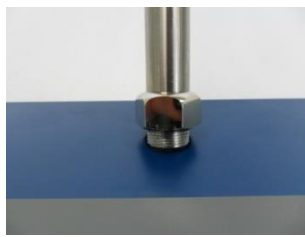


<Figure 4> Stainless Steel Rod

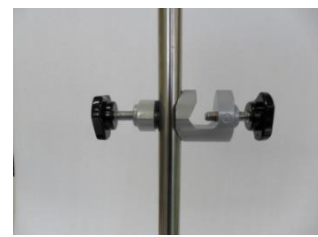

< Figure 5> Clamp (CL200)

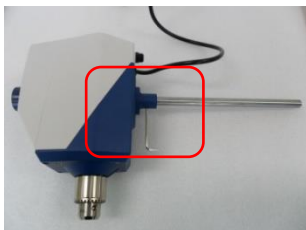

< Figure 6> Stirrer, support Rod and L-Wrench

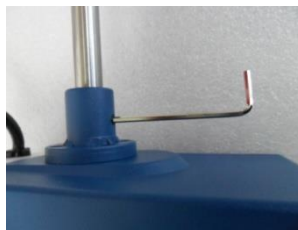

< Figure 7> Stirrer, Support Rod and Clamp



< Figure 8> Drill Chuck, Impeller and Chuck Key



< Figure 9> Completed