BML

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Contents

1 Introduction	12
2 Safety Instructions	12
2.1 Description of symbols	12
2.2 Instructions	13
3 Packing contents.....	14
4 Product description	14
4.1 Key features	14
4.2 General survey.....	15
4.3 Operating surface	15
4.4 Term description	16
5 Positioning the device and commissioning.....	16
6 Job description.....	17
6.1 Quick start	17
6.2 Create Program	17
6.3 Start Program	18
6.4 Storage function	19
7 Maintenance and cleaning.....	19
8 Troubleshooting Guide	19
9 Technical data.....	20
10 Certificate.....	21

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Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you have to follow the safety instructions of the second chapter.

Please keep this operating manual in a save place for future reference and hand it out if you consign the device to a third person.

Witeg reserves to modify the capsule or make amendments for the improvement of quality and efficiency without prior notice.

2 Safety Instructions

2.1 Description of symbols

**WARNING**

Information for avoiding injuries or fatal accidents.

**ATTENTION**

Safety instruction to avoid a damage of the product.

**ATTENTION**

Information about technical parameters.

**Environment**

Information about the operating environment.

**WARNING of explosions**

Safety instruction to avoid explosions.

**ATTENTION to high temperatures**

Instructions to avoid injuries caused by high temperatures.

2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Connect the device only to the suitable power supply. Please have a look to the identification plate.
- Use the device only in electric circuits with ground fault circuit interrupter.
- Check the power supply cord before you stick it into the electrical outlet.
- Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.
- Do not touch the device with wet hands during the usage because this might cause an electrical shock.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.

- Pay attention if you open the door, hot air is coming out.
- Do not put heat-sensitive items at the device.
- Warn all attendees if you leave the device unattended.
- Only use adequate sized vessels and avoid an over boil of liquids.
- Do not clean the scorching device.

- Reparations and amendments should only made by **witeg** authorized people. Contraventions exclude the warranty.

- Use this device only if you read the whole safety instructions.
- This device should only used by persons who are familiarized with safety instructions made for laboratories.
- Pay attention that no liquid reach the controller or inward of the capsule.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective work clothing during the usage.
- All circuits and electrotechnical components are **witeg** owned technology. In fact of this it is necessary that only **witeg** authorized personal repair the device.
- Pull out the power cable if you didn't want to use the device for a longer time.

- It is important that the device stands up on a horizontal and firm subsurface to avoid vibrations.
- If you set the device in a site with vibrations, make sure that the device can not slip away.
- The device could get damaged if the vibrations are too strong.
- To avoid heat congestion or fires make sure that there is enough space for air circulations.
- Avoid direct solar radiation.
- Do not use the device near highly combustible materials.
- Do not set the device outdoors.
- It is superiorly to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- To set the device in wet or dusty places could cause overheating, short circuits or fires.

- Do not use this device in area with potentially explosive atmosphere.

- Do not put your hands in the running device.

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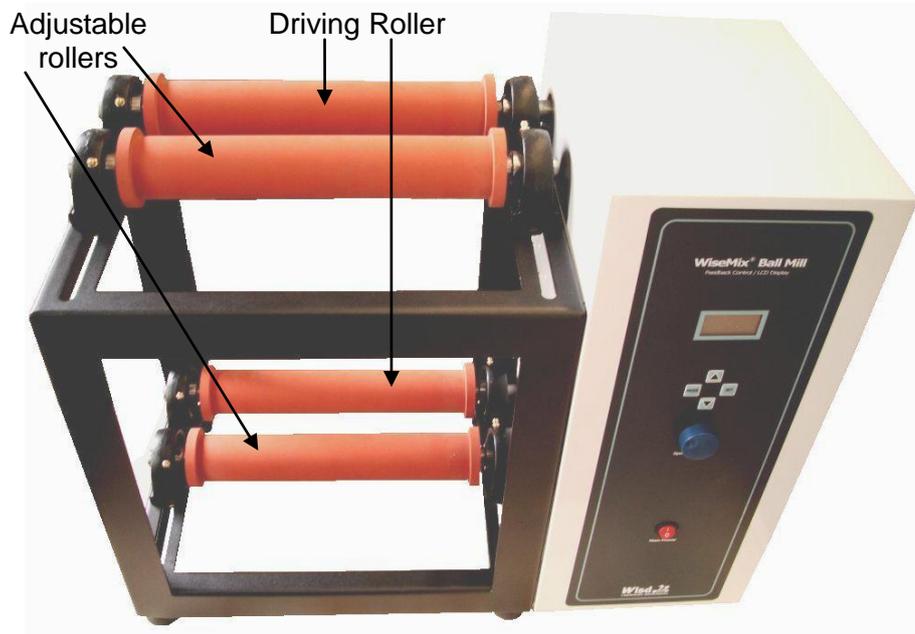
Key features are as follows:


- Simplest control by Jog-Shuttle knob.

- Fast commissioning.
- Long life by simply and solid construction.
- Digital LCD with Back-Light.
- Programmable: 6 programs a 10 steps.

4.2 General survey


Grease nipple

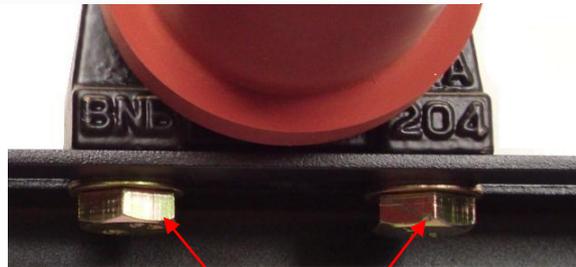Fixing screws for adjustable rollers

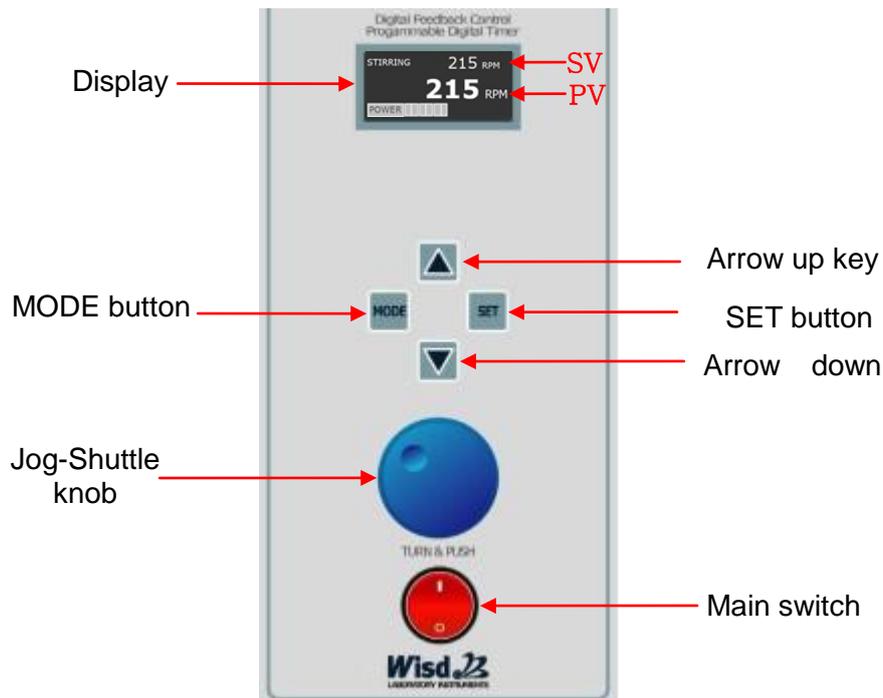4.3 Operating surface

4.4 Term description

SV: set value (upper line, in key lock menu items no SV)

PV: present value, measured by the device (lower line, in key lock menu item also upper line)

Main switch

- Switches the device on and off.

Jog-Shuttle knob

- Change SV by turning.
- Push the knob for start and stop or probate set values.

MODE button

- Display reset
- Change menu item

SET button

- Required for start, stop, storage, and menu change.

Arrow keys

- Required for selection of programs and steps.

5 Positioning the device and commissioning

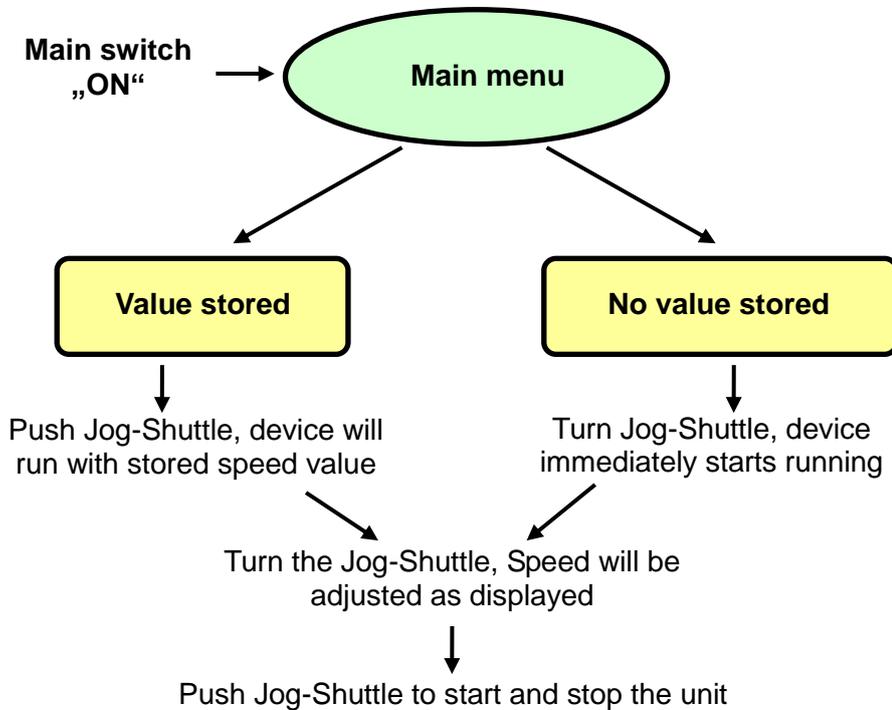
1. Follow especially the safety instructions marked with following symbol:
2. Remove the package carefully.
3. If there are any unexpected transport damages, please inform your forwarder and your customer service.
4. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatised.
5. Make sure that you follow all safety instructions.

6. Adjust the space between the rollers depending on the pod size. Therefore you need a 17 mm screw-wrench. Pay attention that the rollers are parallel to each other.
7. Connect the device with the power outlet.
8. Switch the device on.
9. Set your values and start the process (look at job description).

6 Job description


6.1 Quick start

Working without timer and program function.


6.2 Create Program

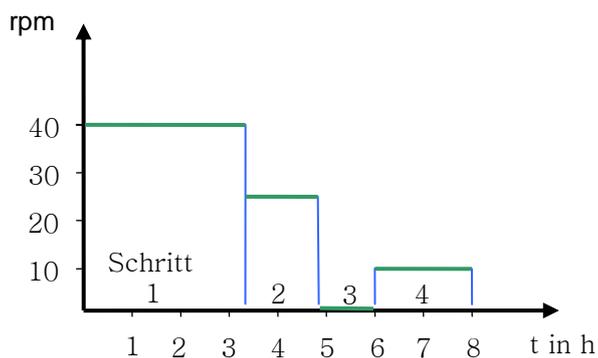
The program function provides automatic operation with different speed levels and breaks. You can create 6 programs with 10 steps each:

The following picture shows the menu structure. The descriptions in the diagram explain how to create a program.

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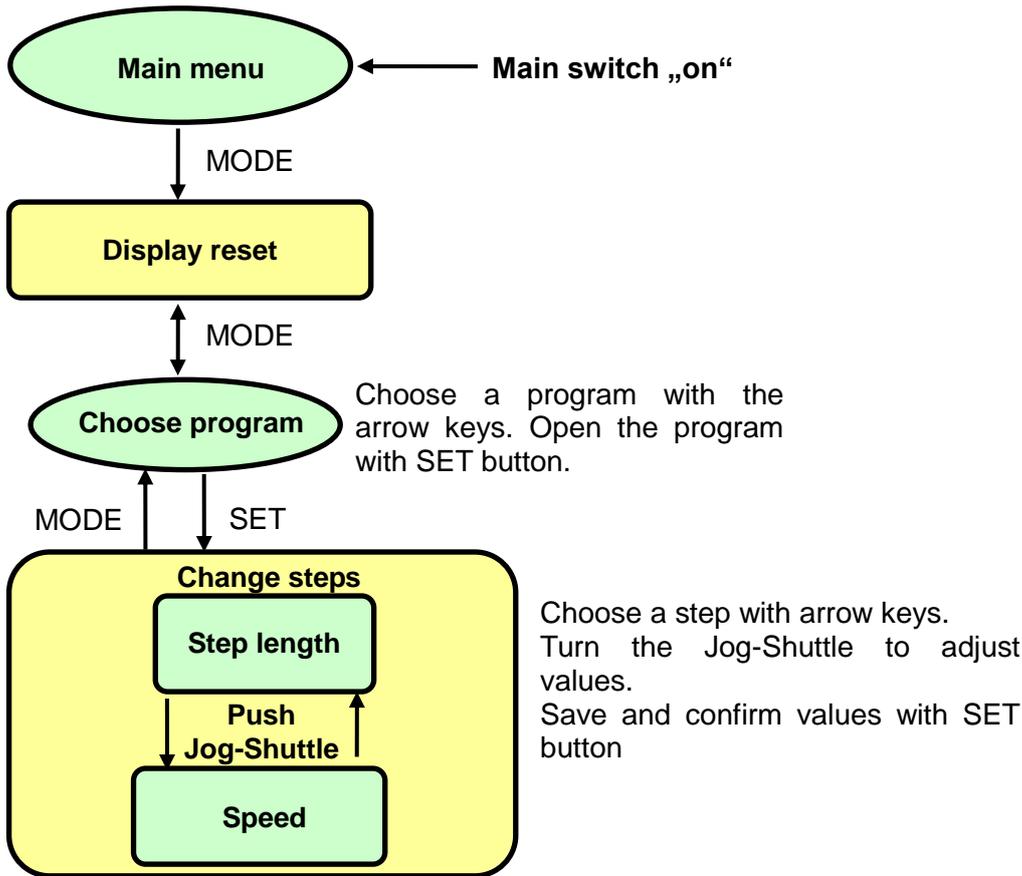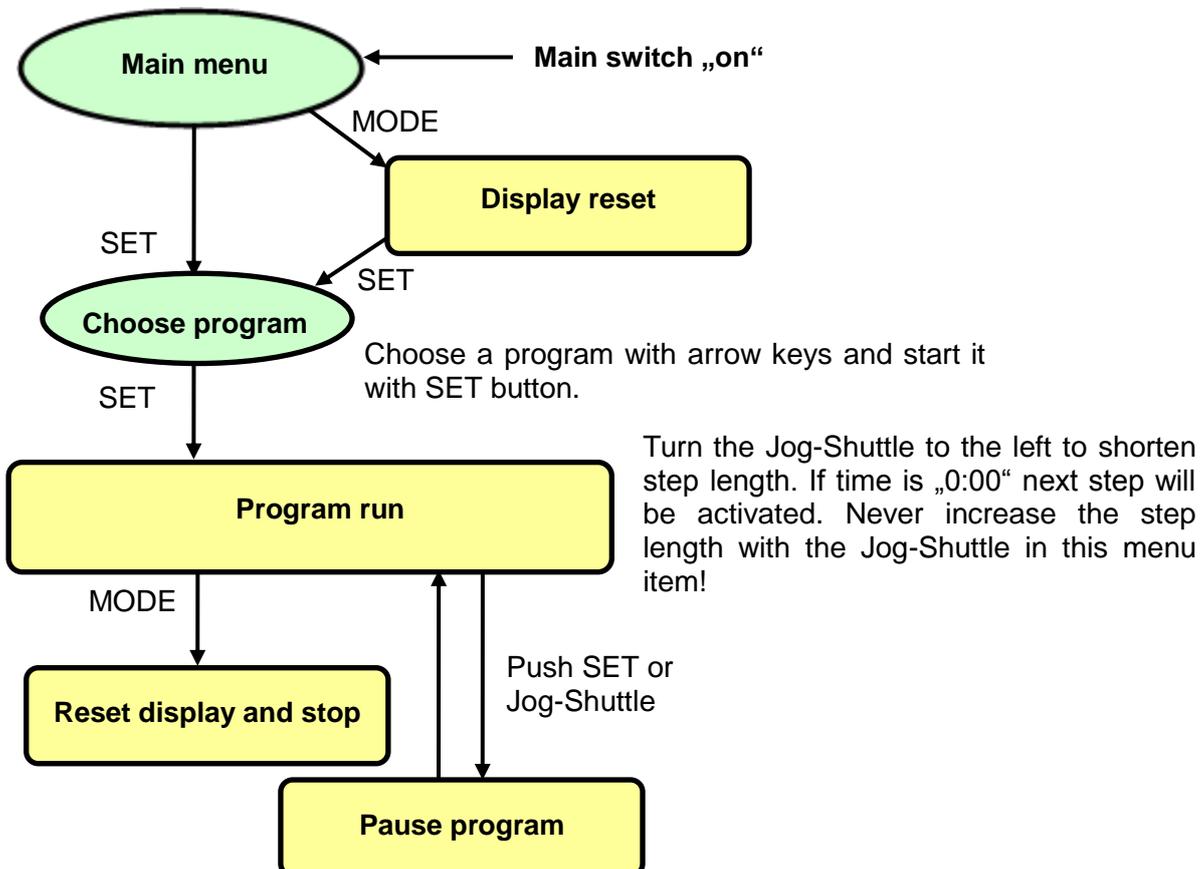- For storing values you only need to hold the Jog-Shuttle knob pushed for two seconds. Successful storing is displayed by blinking of the SV for three times.
- Any stored parameters are erasable. Turn the value with the Jog-Shuttle till "0" and then hold the knob pressed for two seconds.
- If you want to overwrite a value it is not necessary to erase the old one. Set the desired value and use the Jog-Shuttle knob for storing.

7 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device.
- Pull the power plug out of the electrical outlet.
- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- Only use original spare parts.
- Grease the device minimum after every 6 month. Check the bearings more frequent if you use the device very often and with very high load. Grease the bearings by the grease nipples and a lubrication gun.

8 Troubleshooting Guide

Situation	Confirmations and Solutions
Power Failure	<ul style="list-style-type: none"> ✓ Check the supplied voltage is present. ✓ Check the power connection and the power cord for proper connection and integrity. ✓ Check if the MAIN Power Switch is "ON". ✓ Fuse is blown.
Rolling operation without function	<ul style="list-style-type: none"> ✓ Check if the MAIN Power Switch is "ON". ✓ Check the set RPM.

If other problems arise, or one of the above not solves problems, please contact your official agent or the manufacturer.

Reparations and amendments should only made by [witeg](#) authorized people. Contraventions exclude the warranty.

9 Technical data

	BML-2	BML-6
Capacity	2 pods	6 pods
Housing (B x T x H mm)	620 x 320 x 520	1.200 x 515 x 795
Rollers (L x Ø mm)	300 x 50	800 x 70
Maximum load	40 kg	120 kg
Motor power	100 W	
Speed (rpm)	50 - 600	
Controller	Digital, programmable	
Memory	6 programs with 10 steps each	
Timer	99 h : 59 min	
Hosing material	Powder coated steel	
Weight (Kg)	32,5	80
Power supply	AC 230 V, 1N~, 50/60 Hz	

10 Certificate

Certificate

for

The quality and all features were checked by the manufacturer before the shipment.

We grant from date of purchase

two years guaranty.

This certificate excludes damages by natural disasters or incorrect usages by the costumer.

Please look on your account and complete following table:

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witeg Labortechnik GmbH