## HP-D

# Heizplatte

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# Hotplate

## Operating Manual

gilt für / written for  
HP-20D / -30D

**witeg Labortechnik GmbH**

(English manual at page 11)

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# Contents

<b>1 Introduction .....</b>	<b>13</b>
<b>2 Safety Instructions .....</b>	<b>13</b>
2.1 Description of Symbols .....	13
2.2 Instructions .....	13
<b>3 Product description .....</b>	<b>15</b>
3.1 Key features .....	15
3.2 General Survey.....	15
3.3 Operating surface .....	16
3.4 Term description .....	16
<b>4 Packing contents.....</b>	<b>16</b>
4.1 Standard score of delivery .....	16
4.2 Additional accessories .....	17
<b>5 Positioning the device and commissioning.....</b>	<b>17</b>
<b>6 Job description.....</b>	<b>18</b>
6.1 Menu structure .....	18
6.2 Set temperature.....	19
6.3 Justification.....	19
6.4 Set turn-off delay .....	19
6.5 Key lock.....	19
6.6 Storage Funktion .....	19
<b>7 Maintenance and cleaning.....</b>	<b>20</b>
<b>8 Troubleshooting Guide .....</b>	<b>20</b>
<b>9 Technical data.....</b>	<b>21</b>
<b>10 Certificate.....</b>	<b>22</b>

# 1 Introduction

Thank you for purchasing our **HP-D Hotplate**. You did a good choice buying a **Witeg** product. Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you have to follow the safety instructions of the second chapter.

Please keep this operating manual in a save place for future reference and hand it out if you consign the device to a third person.

**Witeg** reserves to modify the capsule or make amendments for the improvement of quality and efficiency without prior notice.

## 2 Safety Instructions

### 2.1 Description of Symbols

**WARNING**

Information for avoiding injuries or fatal accidents.

**ATTENTION**

Safety instruction to avoid a damage of the product.

**ATTENTION**

Information about technical parameters.

**Environment**

Information about the operating environment.

**WARNING of explosions**

Safety instruction to avoid explosions.

**ATTENTION to high temperatures**

Instructions to avoid injuries caused by high temperatures.

### 2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Connect the device only to the suitable power supply. Please have a look to the identification plate.
- Use the device only in electric circuits with ground fault circuit interrupter.
- Check the power supply cord before you stick it into the electrical outlet.
- Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.

- Do not touch the device with wet hands during the usage because this might cause an electrical shock.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.

- The hotplate could be scorching during and after the usage.
- Do not put heat-sensitive items at the hot plate.
- Warn all attendees if you leave the device unattended.
- Only use adequate sized vessels and avoid an over boil of the liquids.
- Do not clean the scorching ceramic suited hotplate.

- Reparations and amendments should only made by **witeg** authorized people. Contraventions exclude the warranty and may cause fatal injuries.

- Use this device only if you read the whole safety instructions.
- This device should only used by persons who are familiarized with safety instructions made for laboratories.
- Pay attention that no liquid reach the controller or inward of the capsule.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective work clothing during the usage.
- All circuits and electro technical components are **witeg** owned technology. In fact of this it is necessary that only **witeg** authorized personal repair the device.
- Pull out the power cable if you didn't want to use the device for a longer time.

- It is important that the device stands up on a horizontal and firm subsurface.
- Make sure that the device can not slip away.
- To avoid heat congestion or fires make sure that there is enough space for air circulations.
- Avoid direct solar radiation.
- Do not use the device near highly combustible materials.
- Do not set the device outdoors.
- It is superiorly to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- To set the device in wet or dusty places could cause overheating, short circuits or fires.

- Do not use this device in area with potentially explosive atmosphere.

## 3 Product description

### 3.1 Key features

The **Wisd HP-D Hotplate** is designed for most chemical, biological, medical and pharmaceutical drying and heating applications. The device was developed for simplest usage and a very long life.


Key features are as follows:

- Long life by simply and solid construction.
- Ceramic coated heating plate for high chemical resistance.
- Simplest operating by patented combination of MODE button and Jog-Shuttle knob.
- Fast commissioning.
- Illuminated LCD indicates temperature and time.
- Max. Temperature 380 °C, adjustable in 0,1 °C-steps.
- Turn-off delay timer (99h 59 min).
- Signal sound indicates elapsed delay time.
- Key lock against inadvertent changing of set temperature and timer values.
- Easy to connect extern PT100 temperature sensor for direct medium measurement.


### 3.2 General Survey

(HP-20D is pictured)

The functionality of the described components are the same at each type of the **HP-D-series**. Only the complexion differs from type to type.

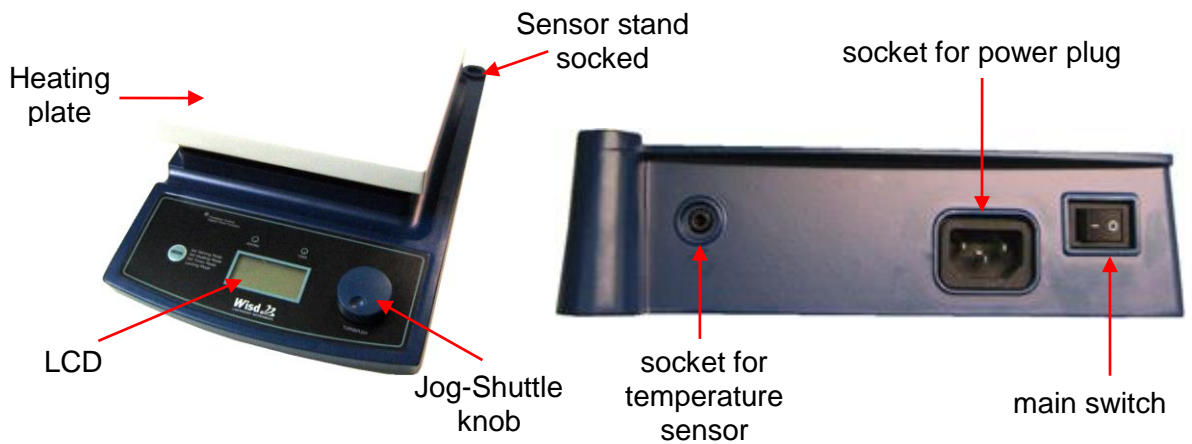### 3.3 Operating surface
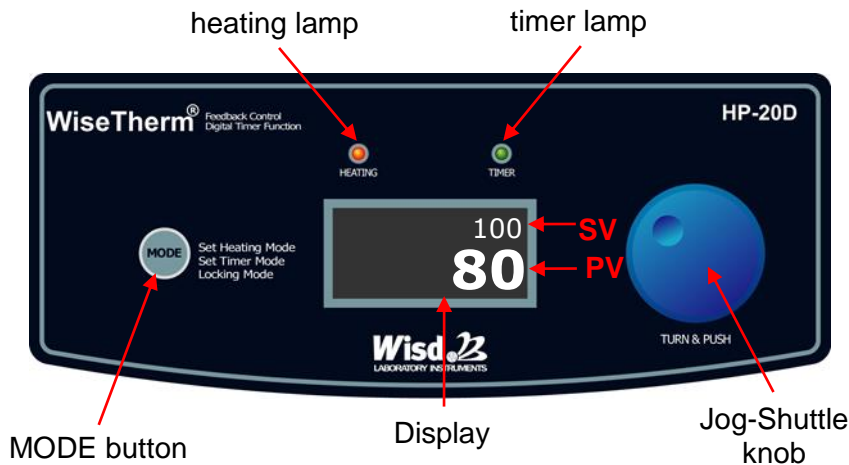
### 3.4 Term description

**SV:** set value (upper line in LCD)

**PV:** present value, measured by the device (lower line in LCD)

#### MODE button

- For keyboard shortcut and change between the menu items.

#### Jog-Shuttle knob

- Change SV-values by turning.
- Push the knob for start and stop or probate and save set values.
- Also used for keyboard shortcut.

#### Main switch

- On / off switch.

#### Timer lamp

- Is blinking in menu items for timer modulation.

#### Heating lamp

- Shines if the heater is activated.

#### Keyboard shortcut

- For reaching justify menu.

## 4 Packing contents

### 4.1 Standard score of delivery

quantity	DH.WHP03021	DH.WHP03020
	HP-20D Hotplate	
1	power supply curt	
1	operating manual	
1	display scratch cover	
1	CL220 clamp + sensor holder	
1	SS500 PT100 temperature Sensor	
1	RD200 stand rod, Ø 12 mm, length 450 mm	

quantity	DH.WHP03023	DH.WHP03022
	<b>HP-30D Hotplate</b>	
1	power supply curt	
1	operating manual	
1	<b>CL220</b> clamp + sensor holder	
1	<b>SS500</b> PT100 temperature Sensor	
1	<b>RD100</b> stand rod, Ø 10 mm, length 450 mm	


#### 4.2 Additional accessories

order number	denomination
DH.WHP503060	<b>SS500</b> temperature sensor, PT100 for HP-20D/ -30D
DH.WHP503061	<b>SS510</b> PTFE-coated temperature sensor PT100 for HP-20D/ -30D
DH.WHP503011	<b>RD100</b> stainless steel stand, Ø 12mm, length 450mm for HP-30D
DH.WHP503036	<b>RD200</b> stainless steel stand rod, Ø 10mm, length 450mm for HP-20D
DH.WHP503012	<b>CL200</b> clamp + holder for temperature sensor

## 5 Positioning the device and commissioning

1. Follow especially the safety instructions marked with following symbol:
2. Remove the package carefully.
3. If there are any unexpected transport damages, please inform your forwarder and your costumer service.
4. If you want to use the device without additional accessories jump to point 10 otherwise follow the next points.
5. Attach the stand rod to the socked.
6. Fix the clamp at the stand rod.
7. Attach the sensor holder to the clamp. Have a look at the next picture if you are unsure. Ensure that the bigger drill hole is upside.
8. Stick the sensor into the holder and connect the cable with the device's sensor socked backwards.
9. Check that each connection fits close.
10. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatised.
11. Make sure that you follow all safety instructions.
12. Connect the device with the power outlet.
13. Switch the device on.
14. Set your values and start the process (look at job description).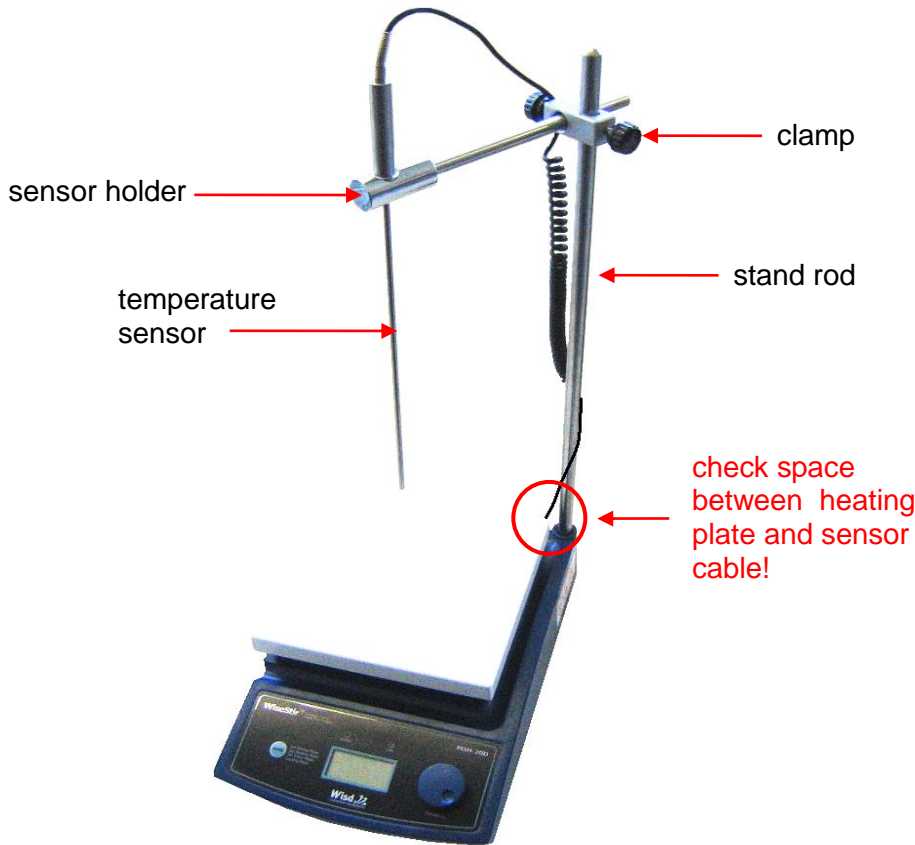


## 6 Job description

### 6.1 Menu structure

The **HP-D Hotplate** proffers you the genial combination of MODE button and Jog-Shuttle knob. That allows a simple and comfortable usage.

The menus structure is well arranged and pictured as follows:

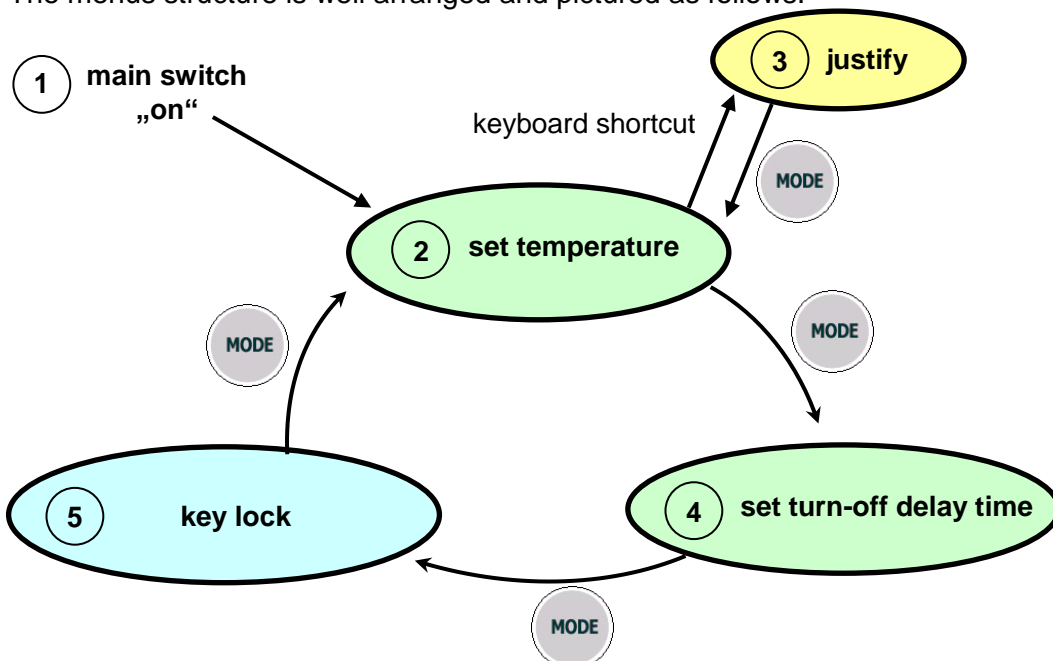## 6.2 Set temperature

- First menu item after switching the device on is “set temperature”. You can recognize this menu at “°C” behind the SV- and PV-value.
- The temperature is adjustable from 25 °C till 380 °C in 0,1 °C-steps.
- Choose your desired temperature by turning the Jog-Shuttle and confirm the set value by pushing the Jog-Shuttle.
- For changing the confirmed value again you first have to push the Jog-Shuttle knob. “HEATING” starts to blink at the LCD and the temperature is adjustable again.
- When the heating lamp is shining, the heater is activated.

## 6.3 Justification

If the temperature measured by you differs from the temperature the device is showing (PV) you have the possibility to compensate the divergence by the offset. So you can balance differences of about  $\pm 20$  °C.

Calculate the offset as follows:

**Temperature measured by you – showed Temperature (PV) = Offset**

- To reach this menu you have use the keyboard shortcut in the menu item “set temperature”.
- Set the calculated value by turning the Jog-Shuttle. The Offset will be added to the PV immediately.
- Save the offset by holding the Jog-Shuttle pushed until the values in PV and SV line starts blinking.
- Push the MODE button to leave the justification menu.

## 6.4 Set turn-off delay

The turn-off delay timer serves as automatically stop function. It turns off the heating after elapsing of the timeframe you set.

- To reach the relevant menu item you have to push the MODE button in the menu “set temperature”.
- The delay time is adjustable from one minute till over 99 hours (00:01 - 99:59) in one minute steps.
- “00:00” stands for endurance run.
- Turn the Jog-Shuttle for choosing your desired value and probate it by pushing the Jog-Shuttle. The delay time starts to run by probating the value.
- If you want to change the confirmed value, you have to push the Jog-Shuttle once again.

## 6.5 Key lock

- To avoid an accidental changing of the set values, push the MODE button in the menu item “set temperature”.
- In this menu item can you see the actually temperature in the SV-line and the remaining time of turn-off delay in the PV-line. The function of the Jog-Shuttle knob is deactivated.
- Press the MODE button once more to leave the key lock.

## 6.6 Storage Funktion

Save your precious time and store often used parameters. Stored values only need to be confirmed by pushing the Jog-Shuttle knob.

- For storing values you only need to hold the Jog-Shuttle knob pushed for two seconds. Successful storing is displayed by blinking of the SV for three times.

- Any stored parameters are erasable. Turn the value with the Jog-Shuttle till “0” and then hold the knob pressed for two seconds.
- If you want to overwrite a value it is not necessary to erase the old one. Set the desired value and use the Jog-Shuttle knob for storing.

## 7 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device and pull the power plug out of the electrical outlet.
- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- Only use original spare parts.

## 8 Troubleshooting Guide

Situation	Confirmations and Solutions
Power Failure	<ul style="list-style-type: none"> <li>✓ Check the supplied voltage is present.</li> <li>✓ Check the power connection and the power cord for proper connection and integrity.</li> <li>✓ Check if the MAIN Power Switch is “ON”.</li> <li>✓ Fuse is blown.</li> </ul>
Unit Not Heating	<ul style="list-style-type: none"> <li>✓ Check if the MAIN Power Switch is “ON”.</li> <li>✓ Check if the set temperature (SV) confirmed with the Jog-Shuttle knob (by turn the Jog-Shuttle knob cannot change the set temperature – see point 6.2).</li> </ul>
Heating, but unit not reaching temperature or temperature is fluctuated	<ul style="list-style-type: none"> <li>✓ Problem at the sensor or main controller.</li> </ul>
ERROR 1	<ul style="list-style-type: none"> <li>✓ Hotplate defect.</li> </ul>

**If other problems arise, or one of the above not solves problems, please contact your official agent or the manufacturer.**

**Reparations and amendments should only made by [witeg](http://www.witeg.de) authorized people. Contraventions exclude the warranty.**

## 9 Technical data

		HP-20D	HP-30D
<b>Dimension (mm) (BxTxH)</b>		206x307x99	286x387x99
<b>Plate size (mm)</b>		180x180	260x260
<b>Controller</b>		digital Fuzzy-controller with Jog-Shuttle knob (turn + push)	
<b>Tem. range (°C)</b>		25 - 380	
<b>Temp. control resolution (°C)</b>		± 0,5	
<b>Accuracy (°C)</b>		± 0,3	
<b>Heating power (W)</b>		600	1200
<b>Display</b>		illuminated LCD	
<b>Material</b>	<b>Plate</b>	ceramic coated Aluminium	
	<b>Housing</b>	powder coated aluminium	powder coated steel
<b>Timer</b>		turn-off delay 99h 59min	
<b>Weight (kg)</b>		2,8	3,2
<b>Others</b>		<ul style="list-style-type: none"> <li>- storage function for temperature, delay time and offset               <ul style="list-style-type: none"> <li>- key lock (Jog-Shuttle input disabled)</li> <li>- easy justification</li> </ul> </li> <li>- signal sound for elapsed delay time and sensor error</li> </ul>	
<b>Accessories</b>		<ul style="list-style-type: none"> <li>- stand for temperature sensors</li> <li>- several PT100 temperature sensors</li> </ul>	
<b>Power supply</b>		AC 230 V, 1N~, 50/60 Hz	

## 10 Certificate

# Certificate

for

**The quality and all features were checked by the manufacturer before the shipment.**

We grant from date of purchase

**two years guaranty.**

This certificate excludes damages by natural disasters or incorrect usages by the costumer.

Please look on your account and complete following table:

<b>Article</b>	<b>HP-D Hotplate</b>
<b>Type</b>	
<b>Serialno.</b>	
<b>Date</b>	

**witeg Labortechnik GmbH**

Am Bildacker 16  
D-97877 Wertheim  
TEL: +49-9342-9301-0  
FAX: +49-9342-9301-77  
Email: info@witeg.de  
**www.witeg.de**