## HP-A

# Heizplatte

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# Hotplate

Operating Manual

gilt für / written for  
HP-20A / -30A

**witeg Labortechnik GmbH**

(English manual at page 7)

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# Contents

<b>1 Introduction .....</b>	<b>8</b>
<b>2 Safety Instructions .....</b>	<b>8</b>
<b>2.1 Description of Symbols .....</b>	<b>8</b>
<b>2.2 Instructions .....</b>	<b>8</b>
<b>3 Product description .....</b>	<b>10</b>
<b>3.1 Key features.....</b>	<b>10</b>
<b>3.2 General survey .....</b>	<b>10</b>
<b>4 Package content.....</b>	<b>10</b>
<b>5 Positioning the device and commissioning.....</b>	<b>10</b>
<b>6 Job description .....</b>	<b>11</b>
<b>7 Maintenance and cleaning.....</b>	<b>11</b>
<b>8 Technical data.....</b>	<b>11</b>
<b>9 Certificate.....</b>	<b>12</b>

## 1 Introduction

Thank you for purchasing our **HP-A Hotplate**. You did a good choice buying a **Wisd** product. Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you have to follow the safety instructions of the second chapter.

Please keep this operating manual in a save place for future reference and hand it out if you consign the device to a third person.

**Witeg** reserves to modify the capsule or make amendments for the improvement of quality and efficiency without prior notice.

## 2 Safety Instructions

### 2.1 Description of Symbols

**WARNING**

Information for avoiding injuries or fatal accidents.

**ATTENTION**

Safety instruction to avoid a damage of the product.

**ATTENTION**

Information about technical parameters.

**Environment**

Information about the operating environment.

**WARNING of explosions**

Safety instruction to avoid explosions.

**ATTENTION to high temperatures**

Instructions to avoid injuries caused by high temperatures.

### 2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Connect the device only to the suitable power supply.
- Use the device only in electric circuits with ground fault circuit interrupter.
- Check the power supply cord before you stick it into the electrical outlet.
- Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.

- Do not touch the device with wet hands during the usage because this might cause an electrical shock.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.

- The hotplate could be scorching during and after the usage.
- Do not put heat-sensitive items at the hot plate.
- Warn all attendees if you leave the device unattended.
- Only use adequate sized vessels and avoid an over boil of the liquids.
- Do not clean the scorching ceramic suited hotplate.

- Reparations and amendments should only made by witeg authorized people. Contraventions exclude the warranty.

- Use this device only if you read the whole safety instructions.
- This device should only used by persons who are familiarized with safety instructions made for laboratories.
- Pay attention that no liquid reach the controller or inward of the capsule.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective work clothing during the usage.
- All circuits and electro technical components are **witeg** owned technology. In fact of this it is necessary that only witeg authorized personal repair the device.
- Pull out the power cable if you didn't want to use the device for a longer time.

- It is important that the device stands up on a horizontal and firm subsurface.
- Make sure that the device can not slip away.
- To avoid heat congestion or fires make sure that there is enough space for air circulations.
- Avoid direct solar radiation.
- Do not use the device near highly combustible materials.
- Do not set the device outdoors.
- It is superiorly to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- To set the device in wet or dusty places could cause overheating, short circuits or fires.

- Do not use this device in area with potentially explosive atmosphere.

## 3 Product description


### 3.1 Key features

The **Wisdom HP-A Hotplate** is designed for most chemical, biological, medical and pharmaceutical drying and heating applications. The device was developed for simplest usage and a very long life.

Key features are as follows:

- Simplest operating.
- Fast commissioning.
- Long life by simply and solid construction.
- Ceramic coated heating plate for high chemical resistance.
- Up to 380 °C.

### 3.2 General survey

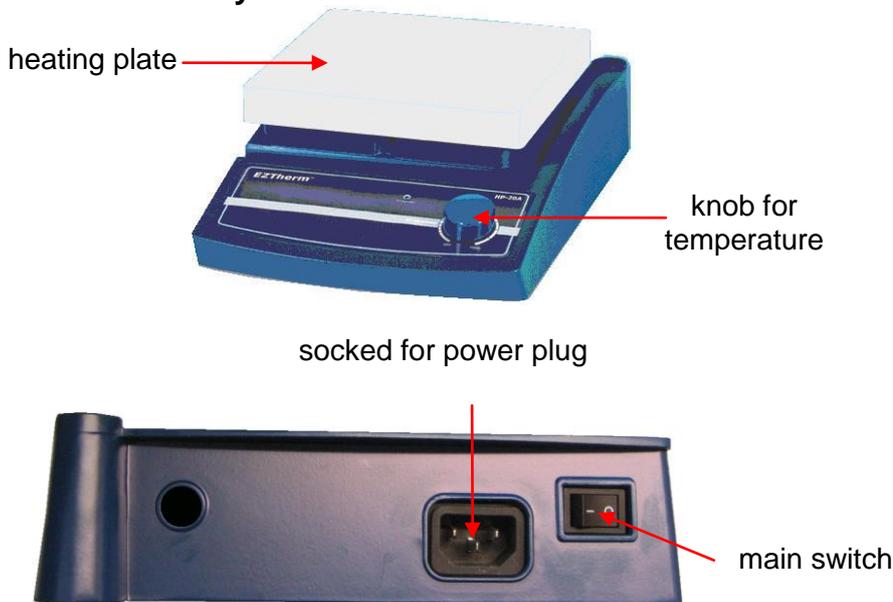## 4 Package content

quantity	denomination
1	<b>HP-A Hotplate</b>
1	power supply curt
1	operating manual

## 5 Positioning the device and commissioning

1. Follow especially the safety instructions indicated with following symbol:
2. Remove the package carefully.
3. If there are any unexpected transport damages, please inform your forwarder and your customer service.
4. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatized.
5. Make sure that you follow all safety instructions.
6. Connect the device with the power outlet.
7. Switch the device on.

## 6 Job description

Turn the knob to set the temperature. Turn it to the left stop to deactivate the heating.

## 7 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device and pull the power plug out of the electrical outlet.
- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- Only use original spare parts.

## 8 Technical data

		HP-20A	HP-30A
<b>Housing Size (mm) (B×T×H)</b>		206x307x99	286x387x99
<b>Plate size (mm)</b>		180x180	260x260
<b>Temperature range</b>		max. 380 °C	
<b>Heating power (W)</b>		600	1200
<b>Material</b>	<b>Plate</b>	ceramic coated Aluminium	
	<b>Housing</b>	powder coated aluminium	powder coated steel
<b>Weight (kg)</b>		2,8	3,3
<b>Power supply</b>		AC 230 V, 1N~, 50/60 Hz	

## 9 Certificate

# Certificate

for

The quality and all features were checked by the manufacturer before the shipment.

We grant from date of purchase

**two years guaranty.**

This certificate excludes damages by natural disasters or incorrect usages by the costumer.

Please look on your account and complete following table:

Article	HP-A Hotplate
Type	
Serialno.	
Date	

**witeg Labortechnik GmbH**

Am Bildacker 16  
D-97877 Wertheim  
TEL: +49-9342-9301-0  
FAX: +49-9342-9301-77  
Email: info@witeg.de  
[www.witeg.de](http://www.witeg.de)