JT-M6C

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Flocculation Tester

Operating Manual

witeg Labortechnik GmbH

(English manual at page 11)

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Contents

1 Introduction	12
2 Safety Instructions	12
2.1 Description of symbols.....	12
2.2 Instructions	13
3 Product description	14
3.1 Key features.....	14
3.2 General survey	15
3.4 Term description	16
4 Packing contents.....	16
5 Positioning the device and commissioning.....	16
6 Job description	17
6.1 Quick start	17
6.2 Create Program	17
6.3 Start Program	18
6.4 Storage function.....	19
7 Maintenance and cleaning.....	19
8 Troubleshooting Guide	19
9 Technical data.....	20
10 Certificate.....	21

1 Introduction

Thank you for purchasing our **JT-M6C Flocculation Tester**. You did a good choice buying a **Wisd** product.

Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you have to follow the safety instructions of the second chapter.

Please keep this operating manual in a save place for future reference and hand it out if you consign the device to a third person.

Witeg reserves to modify the capsule or make amendments for the improvement of quality and efficiency without prior notice.

2 Safety Instructions

2.1 Description of symbols

**WARNING**

Information for avoiding injuries or fatal accidents.

**ATTENTION**

Safety instruction to avoid a damage of the product.

**ATTENTION**

Information about technical parameters.

**Environment**

Information about the operating environment.

**WARNING of explosions**

Safety instruction to avoid explosions.

**ATTENTION to high temperatures**

Instructions to avoid injuries caused by high temperatures.

2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Connect the device only to the suitable power supply. Please have a look to the identification plate.
- Use the device only in electric circuits with ground fault circuit interrupter.
- Check the power supply cord before you stick it into the electrical outlet.
- Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.
- Do not touch the device with wet hands during the usage because this might cause an electrical shock.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.

- Pay attention if you open the door, hot air is coming out.
- Do not put heat-sensitive items at the device.
- Warn all attendees if you leave the device unattended.
- Only use adequate sized vessels and avoid an over boil of liquids.
- Do not clean the scorching device.

- Reparations and amendments should only made by **witeg** authorized people. Contraventions exclude the warranty.

- Use this device only if you read the whole safety instructions.
- This device should only used by persons who are familiarized with safety instructions made for laboratories.
- Pay attention that no liquid reach the controller or inward of the capsule.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective work clothing during the usage.
- All circuits and electrotechnical components are **witeg** owned technology. In fact of this it is necessary that only **witeg** authorized personal repair the device.
- Pull out the power cable if you didn't want to use the device for a longer time.

- It is important that the device stands up on a horizontal and firm subsurface to avoid vibrations.
- If you set the device in a site with vibrations, make sure that the device can not slip away.
- The device could get damaged if the vibrations are too strong.
- To avoid heat congestion or fires make sure that there is enough space for air circulations.
- Avoid direct solar radiation.
- Do not use the device near highly combustible materials.
- Do not set the device outdoors.
- It is superiorly to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- To set the device in wet or dusty places could cause overheating, short circuits or fires.

- Do not use this device in area with potentially explosive atmosphere.

- Do not put your hands in the running device.

3 Product description


3.1 Key features

The **JT-M6C Flocculation Tester** is a multi-overhead stirrer. In first line of this device was developed for Flocculation Testing in the field of wastewater treatment. The **JT-M6C Flocculation Tester** is also suitable for many other applications. The device was developed for simplest usage and a very long life.


Key features are as follows:

- Simplest control by Jog-Shuttle knob.
- Fast commissioning.
- Long life by simply and solid construction.
- Digital LCD with Back-Light.
- Central lighting for good visibility of the samples.
- Programmable: 6 programs a 10 steps.

3.2 General survey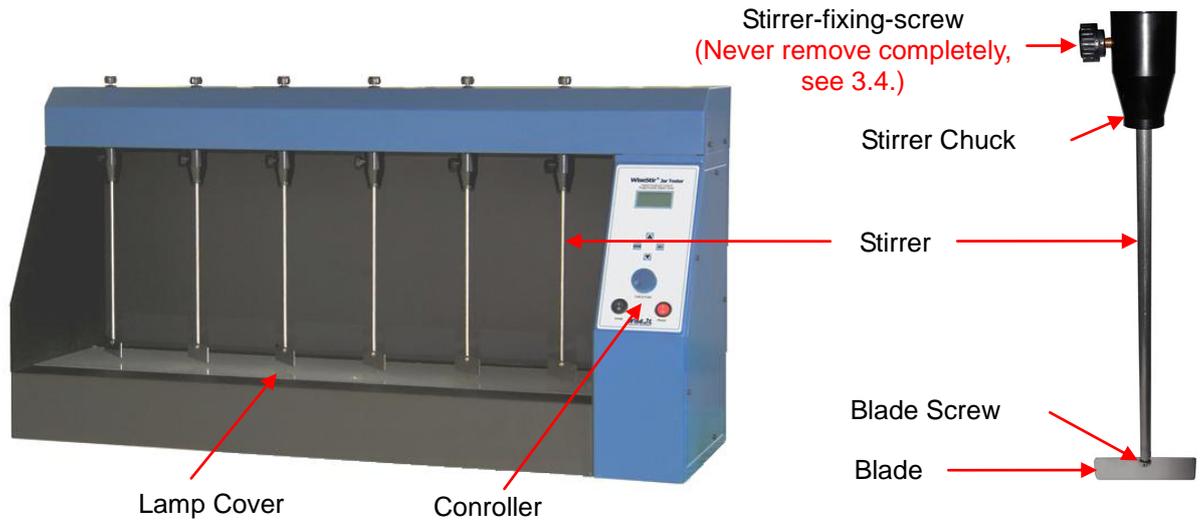

3.3 Operating surface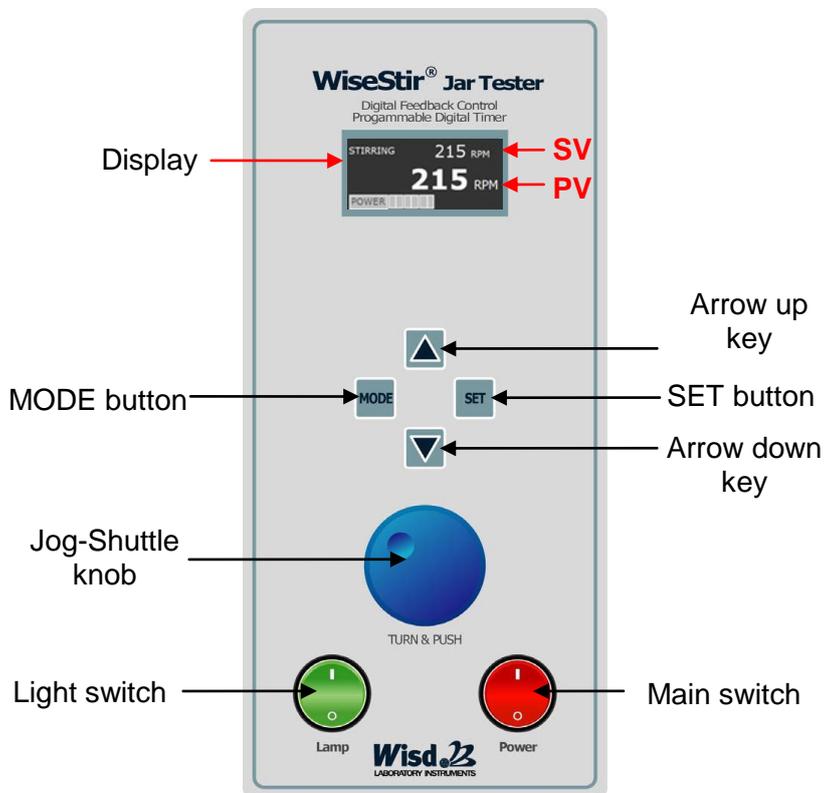

3.4 Term description

SV: set value (upper line, in key lock menu items no SV)

PV: present value, measured by the device (lower line, in key lock menu item also upper line)

Main switch

- Switches the device on and off.

Light switch

- Switches the Light on and off.

Jog-Shuttle knob

- Change SV by turning.
- Push the knob for start and stop or probate set values.

MODE button

- Display reset
- Change menu item

SET button

- Required for start, stop, storage, and menu change.

Arrow keys

- Required for selection of programs and steps.

Blade screw

- To remove the stirring rod.

Stirrer-fixing-screw

- To fix the stirrer.
- **Never remove the stirrer-fixing-screw completely out of the chuck otherwise the 4 mm steel ball will fall out.**

4 Packing contents

quantum	denomination
1	JT-M6C Flocculation Tester
1	Power supply cord
6	stirrers
1	operating manual

5 Positioning the device and commissioning


1. Follow especially the safety instructions marked with following symbol:
2. Remove the package carefully.
3. If there are any unexpected transport damages, please inform your forwarder and your costumer service.
4. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatised.
5. Make sure that you follow all safety instructions.
6. Connect the device with the power outlet.
7. Switch the device on.
8. Set your values and start the process (look at job description).

6 Job description


6.1 Quick start

Working without timer and program function.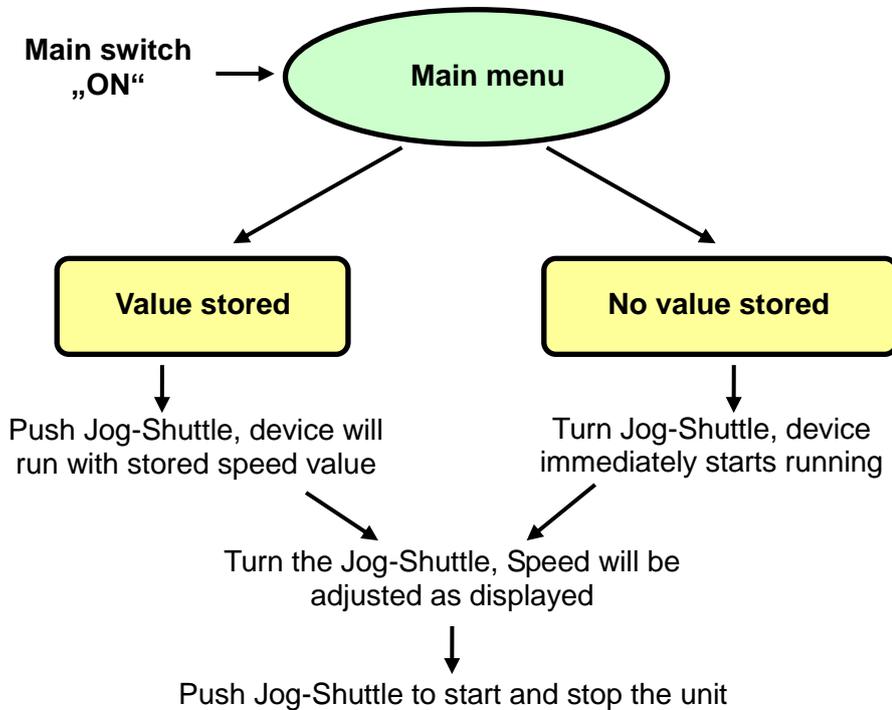


6.2 Create Program


The program function provides automatic operation with different speed levels and breaks. You can create 6 programs with 10 steps each: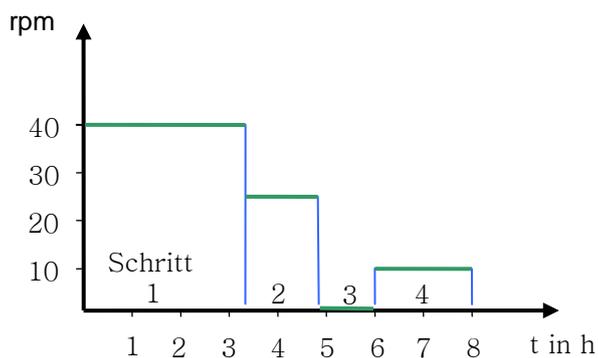

The following picture shows the menu structure. The descriptions in the diagram explain how to create a program.

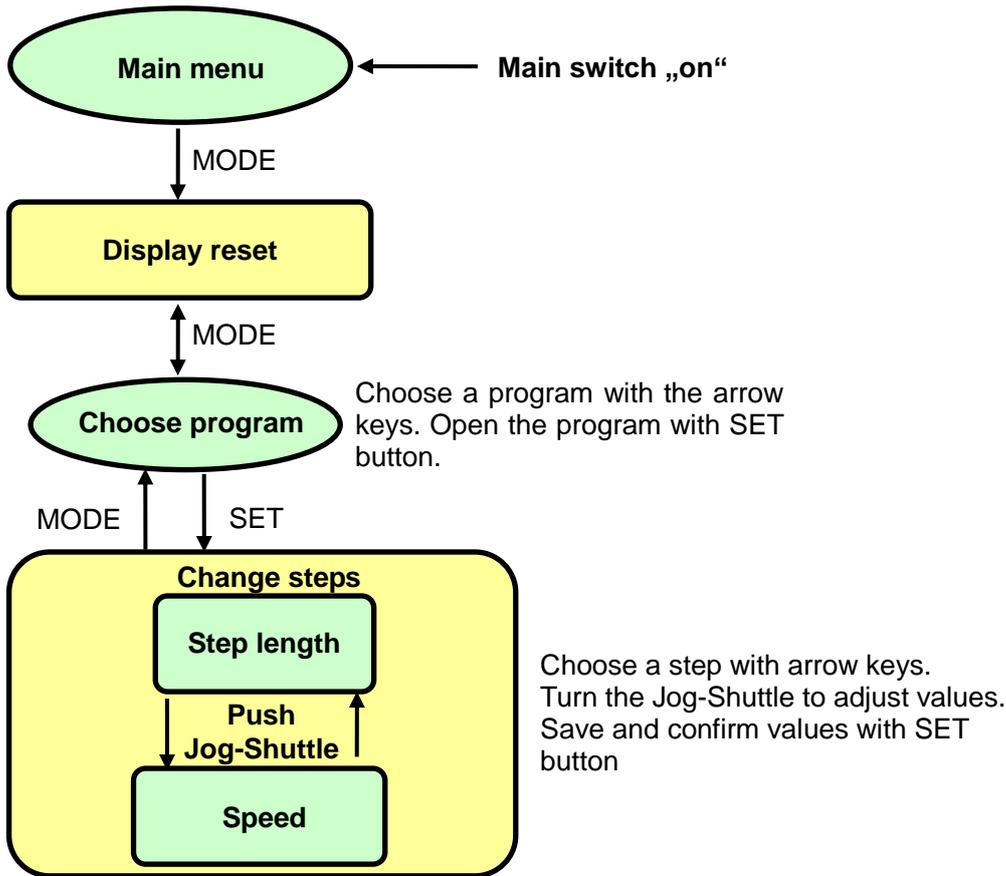6.3 Start Program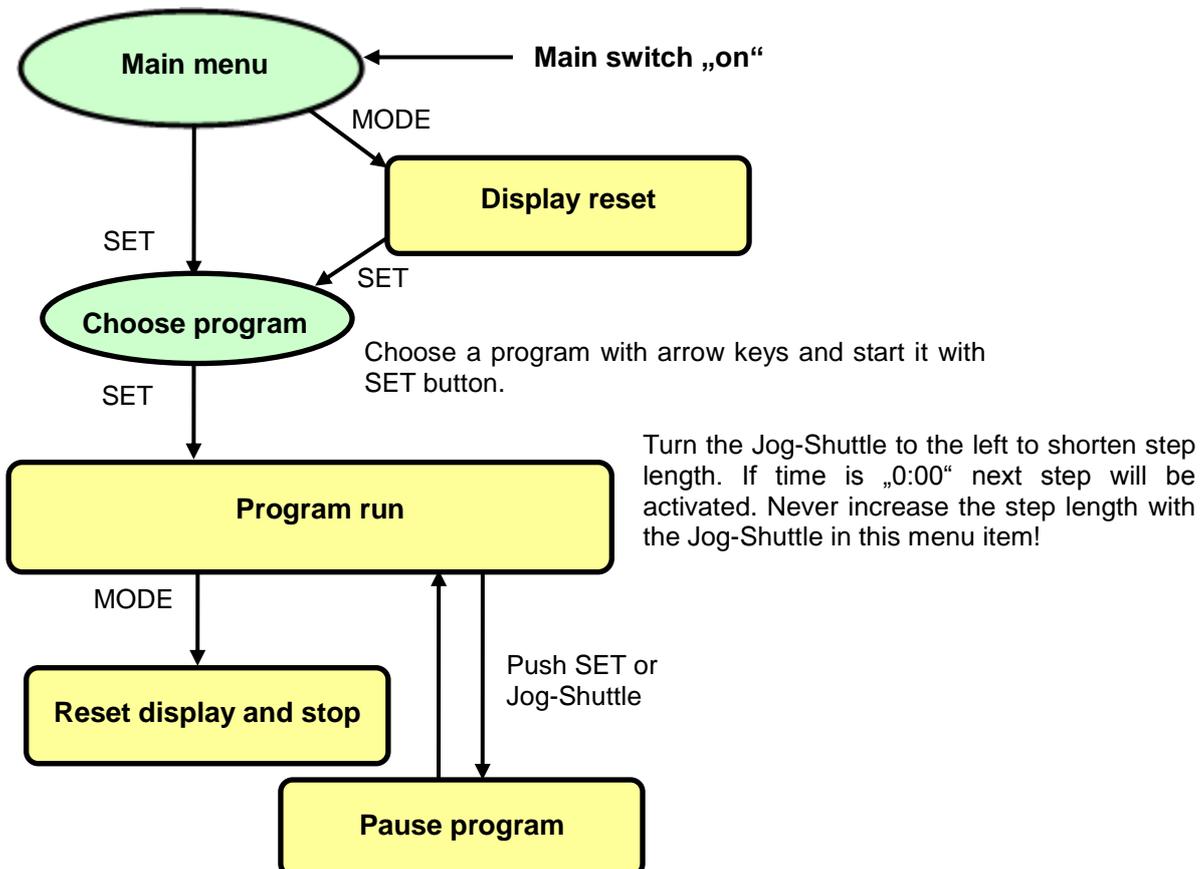

6.4 Storage function

The **JT-M6C Flocculation Tester** includes a storage function for speed values. That is practical if you want to work with the same parameters often. Then you didn't need to set them in every usage.

- For storing values you only need to hold the Jog-Shuttle knob pushed for two seconds. Successful storing is displayed by blinking of the SV for three times.
- Any stored parameters are erasable. Turn the value with the Jog-Shuttle till "0" and then hold the knob pressed for two seconds.
- If you want to overwrite a value it is not necessary to erase the old one. Set the desired value and use the Jog-Shuttle knob for storing.

7 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device.
- Pull the power plug out of the electrical outlet.
- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- Only use original spare parts.

8 Troubleshooting Guide

Situation	Confirmations and Solutions
Power Failure	<ul style="list-style-type: none"> ✓ Check the supplied voltage is present. ✓ Check the power connection and the power cord for proper connection and integrity. ✓ Check if the MAIN Power Switch is "ON". ✓ Fuse is blown.
Unit Not Stirring	<ul style="list-style-type: none"> ✓ Check if the MAIN Power Switch is "ON". ✓ Check the set RPM.

If other problems arise, or one of the above not solves problems, please contact your official agent or the manufacturer.

Repairs and amendments should only made by [witeg](#) authorized people. Contraventions exclude the warranty.


9 Technical data

	JT-M6C
Stirrers	Ø 8 mm x L 400 mm
Stirrer blade	Straight blade
Capacity	6 x 1.000 ml beakers
Stirring places	6
Housing (B x T x H mm)	1.000 x 310 x 535
Motor power	40 W
Speed (rpm)	20 - 300
Controller	Digital, programmable
Memory	6 programs with 10 steps each
Timer	99 h 59 min 59 sec
Hosing material	Powder coated steel
Weight	37 Kg
Lamps	6 x 20 W
Power supply	AC 230 V, 1N~, 50/60 Hz

10 Certificate

Certificate

for

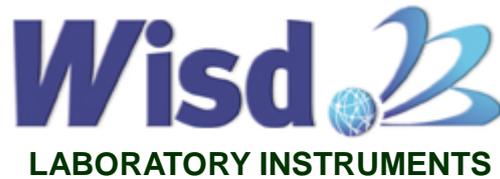The quality and all features were checked by the manufacturer before the shipment.

We grant from date of purchase

two years guaranty.

This certificate excludes damages by natural disasters or incorrect usages by the costumer.

Please look on your account and complete following table:

Article	JT-M6C Flocculation Tester
Typ	
Serialno.	
Date	

witeg Labortechnik GmbH

Am Bildacker 16
D-97877 Wertheim
TEL: +49-9342-9301-0
FAX: +49-9342-9301-77
Email: info@witeg.de
www.witeg.de