# **MS-20A**

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Email: info@witeg.de www.witeg.de

# Contents

1 Introduction	8
2 Safety Instructions	8
2.1 Description of symbols	8
2.2 Instructions	9
3 Product description	10
3.1 Key features	10
3.2 General survey	10
4 Package content	10
5 Positioning the device and commissioning	11
6 Job description	11
7 Maintenance and cleaning	11
8 Technical data	11
9 Certificate	12

## 1 Introduction

Thank you for purchasing our **MS-20A Magnetic Stirrer**. You did a good choice buying a **Wisd** product.

Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you have to follow the safety instructions of the second chapter.

Please keep this operating manual in a save place for future reference and hand it out if you consign the device to a third person.

**Witeg** reserves to modify the capsule or make amendments for the improvement of quality and efficiency without prior notice.

# 2 Safety Instructions

## 2.1 Description of symbols

#### **WARNING**

Information for avoiding injuries or fatal accidents.

#### **ATTENTION**

Safety instruction to avoid a damage of the product.

#### ATTENTION

Information about technical parameters.

## **Environment**

Information about the operating environment.

## **WARNING of explosions**

Safety instruction to avoid explosions.

## 2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Connect the device only to the suitable power supply. Please have a look to the identification plate.
- > Use the device only in electric circuits with ground fault circuit interrupter.
- Check the power supply cord before you stick it into the electrical outlet.
- > Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.
- > Do not touch the device with wet hands during the usage because this might cause an electrical shock.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.

> Do not use this device in areas with potentially explosive atmosphere.

- Use this device only if you read the whole safety instructions.
- > This device should only used by persons who are familiarized with safety instructions made for laboratories.
- > Pay attention that no liquid reaches the controller or inward of the capsule.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective work clothing during the usage.
- > Pull out the power cable if you didn't want to use the device for a longer time.

- Make sure that the device can not slip away.
- ➤ It is important that the device stands up on a horizontal, stabile and firm subsurface to avoid vibrations.
- > To avoid heat congestions or fires make sure that there is enough space for air circulations.
- Avoid direct solar radiation.
- > Do not use the device near highly combustible materials.
- Do not set the device outdoors.
- ➤ It is superiorly to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- > To set the device in wet or dusty places could cause overheating, short circuits or fires.

10

Reparations and amendments should only made by witeg authorized people. Contraventions exclude the warranty.

# 3 Product description


## 3.1 Key features

The **Wisd** MS-20A Magnetic Stirrer is designed for chemical, biological, medical and pharmaceutical stirring-applications. The device was developed for simplest usage and a very long life.

Key features are as follows:

- Simplest operating.
- > Fast commissioning.
- > Long life by simply and solid construction.
- Ceramic coated stirring plate for high chemical resistance.

## 3.2 General survey

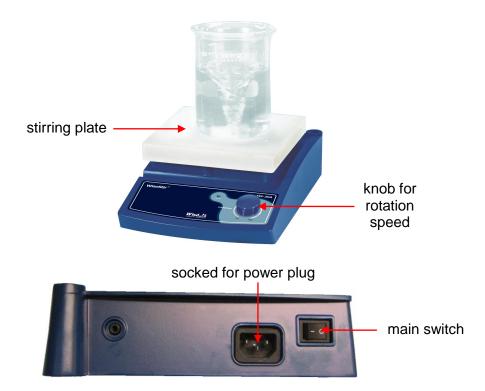# 4 Package content

quantum	denomination
1	MS-20A Magnetic Stirrer
1	power supply curt
1	operating manual

# 5 Positioning the device and commissioning

- 1. Follow especially the safety instructions marked with following symbol:
- 2. Remove the package carefully.
- If there are any unexpected transport damages, please inform your forwarder and your costumer service.
- 4. Remove perhaps soiling and package material.
- 5. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatised.
- 6. Make sure that you follow all safety instructions.
- 7. Connect the device with the power outlet.
- 8. Switch the device on and start the process (look at job description).

# 6 Job description

Turn the knob to set the speed and switch the device on or off.

# 7 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device.
- > Pull the power plug out of the electrical outlet.
- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- Only use original spare parts and accessories.

## 8 Technical data

		MS-20A
Dimension (mm) (w×d×h)		206x307x99
Plate size (mm)		180x180
Speed (U/min)		0 - 1500
Motor Rating (W)		9/4
Stirring capacity (H₂O)		max. 20 l
Matarial	Housing	ceramic coated aluminium
Material	Stirring plate	powder coated aluminium
Weight (kg)		3,1
Power supply		AC 230 V, 1N~, 50/60 Hz

## 9 Certificate

# **Certificate**

or

The quality and all features were checked by the manufacturer before the shipment.

We grant from date of purchase

two years guaranty.

This certificate excludes damages by natural disasters or incorrect usages by the costumer.

Please look on your account and complete following table:

Article	MS-20A Magnetic Stirrer.
Serialno.	
Date	

## witeg Labortechnik GmbH

Am Bildacker 16 D-97877 Wertheim TEL: +49-9342-9301-0 FAX: +49-9342-9301-77 Email: info@witeq.de

www.witeg.de