MS-MP

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Multi-Point Stirrer

Operating Manual

gilt für / written for
MS- MP4 / - MP8

witeg Labortechnik GmbH

(English manual at page 9)

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Contents

1 Introduction	10
2 Safety Instructions	10
2.1 Description of Symbols	10
2.2 Instructions	10
3 Product description	11
3.1 Key features.....	11
3.2 General survey	12
3.3 Display description	12
3.4 Term description	12
4 Packing contend	13
5 Positioning the device and commissioning.....	13
6 Job description	13
6.1 Menu structure	13
6.2 Set speed	13
6.3 Set turn-off delay.....	14
6.4 Key lock	14
6.5 Storage function.....	14
7 Maintenance and cleaning.....	14
8 Troubleshooting Guide	15
9 Technical data.....	15
10 Certificate.....	16

1 Introduction

Thank you for purchasing our **MS-MP Multi-Point Stirrer**. You did a good choice buying a **Wisd** product.

Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you have to follow the safety instructions of the second chapter.

Please keep this operating manual in a save place for future reference and hand it out if you consign the device to a third person.

Witeg reserves to modify the capsule or make amendments for the improvement of quality and efficiency without prior notice.

2 Safety Instructions

2.1 Description of Symbols

WARNING
Information for avoiding injuries or fatal accidents.

ATTENTION
Safety instruction to avoid a damage of the product.

ATTENTION
Information about technical parameters.

Environment
Information about the operating environment.

WARNING of explosions
Safety instruction to avoid explosions.

2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Connect the device only to the suitable power supply. Please have a look to the identification plate.
- Use the device only in electric circuits with ground fault circuit interrupter.
- Check the power supply cord before you stick it into the electrical outlet.
- Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.

- Do not touch the device with wet hands during the usage because this might cause an electrical shock.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.

- Use this device only if you read the whole safety instructions.
- This device should only be used by persons who are familiarized with safety instructions made for laboratories.
- Pay attention that no liquid reaches the controller or inward of the capsule.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective work clothing during the usage.
- Pull out the power cable if you didn't want to use the device for a longer time.

- Do not use this device in areas with potentially explosive atmosphere.

- Make sure that the device can not slip away.
- The device could get damaged if the vibrations are too strong.
- To avoid heat congestions or fires make sure that there is enough space for air circulations.
- Avoid direct solar radiation.
- Do not use the device near highly combustible materials.
- Do not set the device outdoors.
- It is superiorly to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- To set the device in wet or dusty places could cause overheating, short circuits or fires.

- All circuits and electrotechnical components are **witeg** owned technology. In fact of this it is necessary that only **witeg** authorized personal repair the device. Contraventions exclude the warranty.

3 Product description

3.1 Key features

The **MS-MP Multi-Point Stirrer** is designed for chemical, biological, medical and pharmaceutical stirring-applications. The device was developed for simplest usage and a very long life.


Key features are as follows:

- Long life by simply and solid construction.
- Stainless steel stirring plate.
- Jog-Shuttle Knob for easy and fast handling.
- Power bar graph, speed and timer showed on illuminated LCD.
- Storage functions for set time and speed values.
- Turn-off delay from one minute up to over 99 hours in one minute steps.
- Signal sound after end of turn-off delay.
- Key lock against inadvertent changing of set speed and timer values.

3.2 General survey

3.3 Display description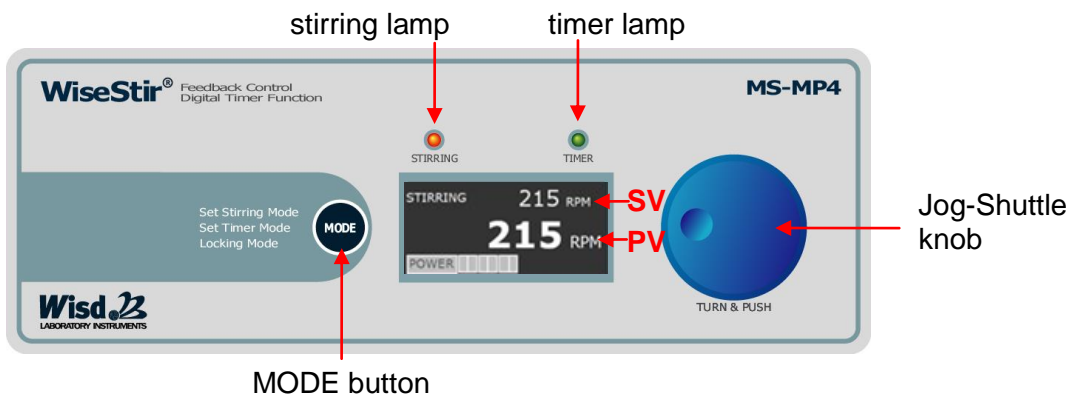

3.4 Term description

SV: set value, upper line in display, in key lock only PVs

PV: present value, lower line in display, in key lock also upper line

MODE button

- For keyboard shortcut and change between the menu items.

Jog-Shuttle knob

- Change SV-values by turning.
- Push the knob for start and stop or probate set values.

Timer lamp

- Is blinking in menu items for timer modulation.
- Shines permanent in the key lock menu if a turn-off delay time is set.

Stirring lamp

- Is blinking in menu items for speed modulation.
- Shines permanent in other menu items during the stirring process.

4 Packing content

quantity	DH.WMS03036	DH.WMS03038
1	MS-MP4	MS-MP8
1	power supply cord	
1	operating manual	
1	anti-slide pad	


5 Positioning the device and commissioning

1. Follow especially the safety instructions marked with following symbol:
2. Remove the package carefully.
3. If there are any unexpected transport damages, please inform your forwarder and your customer service.
4. Remove perhaps soiling and package material.
5. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatised.
6. Make sure that you follow all safety instructions.
7. Connect the device with the power outlet.
8. Switch the device on.
9. Set your values and start the process (look at job description).

6 Job description

6.1 Menu structure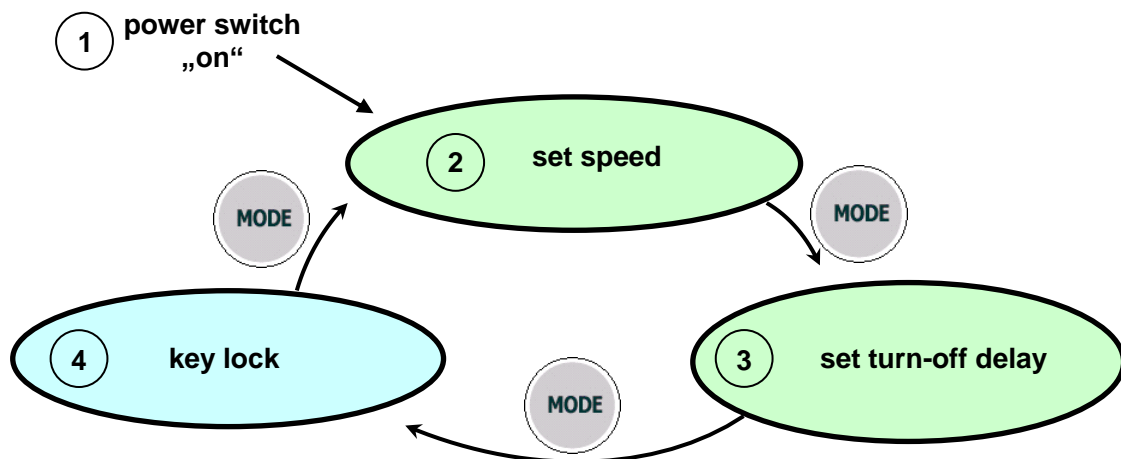

6.2 Set speed

- First menu item after switching the device on is the menu “set speed”. The stirring lamp is blinking. You can recognize the menu at „RPM“ in SV- und PV-line.
- If there is showed “O RPM” in the SV line, there is no value stored and you can start the stirring process by turning the Jog-Shuttle knob clockwise.
- If “STIRRING” is blinking on the LCD there is a speed value stored. The saved value is showed in the SV-line and you can start the stirring process by pushing the Jog-Shuttle knob.
- In both cases is the speed during the stirring process variable. To change the SV you

only have to turn the Jog-Shuttle knob clockwise to increase the speed and counter-clockwise to reduce it.

- To stop the stirring process manually, you have to push the Jog-Shuttle knob again or decrease the speed till "0" by turning the Jog-Shuttle knob counter-clockwise.
- The set speed is increasing gradually, so it could takes some time till the SV is climaxed.

6.3 Set turn-off delay

- The delay timer serves as automatically stop function. It turns off the process after elapsing the timeframe you set.
- To reach the timer menu you have to push the MODE button in the menu item "set speed". The timer lamp is blinking.
- By renewed using of the MODE button you reach the key lock without delay time setting.
- The delay time is adjustable from one minute till over 99 hours (00:01 - 99:59) in one minute steps.
- Turn the Jog-Shuttle knob to set your delay time. Push the Jog-Shuttle if you reached the desired value to confirm it. With the confirmation the timeframe is elapsing
- If you want to change the value after confirmation, you have to push the Jog shuttle once again. The PV initiates blinking and you can modify the value again.
- To confirm the value push the Jog-Shuttle knob yet again.

6.4 Key lock

- To avoid accidental changing of the set values, push the MODE button in the menu item "set delay time".
- In this menu item can you see the actually speed in the SV-line and the remaining time of delay function in the PV-line. The function of the Jog-Shuttle knob is deactivated.
- Press the MODE button once more to leave the key lock. Than you are in the menu item "set speed".

6.5 Storage function

The **MS-MP Multi-Point Stirrer** includes a storage function for speed and delay values. That is practical if you want to work with the same parameters often. Then you need not to set them in every usage.

- For storing values you only need to hold the Jog-Shuttle knob pushed for two seconds. Successful storing is displayed by blinking of the SV-value for three times.
- If you saved a speed value, "STIRRING" is blinking after the device's restart.
- If Every stored parameter is erasable. Turn the value with the Jug-Shuttle till "0" and then hold the knob pressed for two seconds.
- If you want to store a new value it is not necessary to erase the old one.

7 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device and pull the power plug out of the electrical outlet.
- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- Only use original spare parts.

8 Troubleshooting Guide

Situation	Confirmations and Solutions
Power Failure	<ul style="list-style-type: none"> ✓ Check the supplied voltage is present. ✓ Check the power connection and the power cord for proper connection and integrity. ✓ Check if the MAIN Power Switch is "ON". ✓ Fuse is blown.
Unit Not Stirring	<ul style="list-style-type: none"> ✓ Check if the MAIN Power Switch is "ON". ✓ Check the set RPM.
ERROR 2	<ul style="list-style-type: none"> ✓ Errors on the Drive.

If other problems arise, or one of the above not solves problems, please contact your official agent or the manufacturer.

Repairs and amendments should only be made by **witeg** authorized people. Contraventions exclude the warranty.

9 Technical data

		MS-MP4	MS-MP8
Number of Stirring Points		4 (2x2)	8 (4x2)
Distance between Stirring Points		100 mm	
Stirring Capacity per Point		Max. 0.5 Lit.	
Motor Rating		16/9 W	
Speed Range (rpm)		80 ~ 1,200	
Speed Control Resolution (rpm)		5	
Controller		Digital Feedback Controller with Jog Shuttle Switch (Turn + Push)	
Display		Digital LCD Display (Timer, RPM and Power)	
Timer		99hr 59min, (Continuous Function Included)	
Material	Plate	Stainless Steel (#304)	
	Body	Powder Coated Steel	
Dimension (wxdxh mm)		210x280x75	210x480x75
Weight (kg)		2.5	3.5
Others		Storage Function (RPM, Timer), Locking Mode(Jog-Shuttle Input disabled) Alarm (Error Status and Timer-end)	
Power Supply		AC 230 V, 1N~, 50/60 Hz	

10 Certificate

Certificate

for

The quality and all features were checked by the manufacturer before the shipment.

We grant from date of purchase

two years guaranty.

This certificate excludes damages by natural disasters or incorrect usages by the costumer.

Please look on your account and complete following table:

Article	MS-MP Multi-Point Stirrer
Type	
Serialno.	
Date	

witeg Labortechnik GmbH

Am Bildacker 16
D-97877 Wertheim
TEL: +49-9342-9301-0
FAX: +49-9342-9301-77
Email: info@witeg.de
www.witeg.de