Wachstumschammer STH-155/-305/-420/-800
Growth chamber STH-155/-305/-420/-800

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Please study carefully and follow step by step!

GROWTH CHAMBER STH

Content

- 1 General3**
 - 1.1 Precautions before use..... 3
 - 1.2 Product description..... 4
 - 1.3 Product equipment 5
- 2. Installation.....5**
 - 2.1 Product installation 5
 - 2.1.1 Caution of Installation 5
 - 2.1.2 Installation Process 6
 - 2.1.3 Safety Symbols on the product..... 6
 - 2.2 Installation of WiRe™ App & Web Service..... 6
 - 2.3 Product specifications..... 11
- 3. Product Usage12**
 - 3.1 Precautions for Use 12
 - 3.2 Names of Each Part 15
 - 3.3 Usage instructions 18
 - 3.3.1 Temperture setting and operation..... 18
 - 3.3.2 Humidity setting and operation..... 18
 - 3.3.3 Product SETTINGS 19
 - 3.3.4 Chart..... 27
 - 3.3.5 Offset..... 29
 - 3.3.6 Error messages 30
- 4 Products management.....31**
 - 4.1 Accessories 31
 - 4.2 Maintenance 31
 - 4.3 Troubleshooting 32
- 5 Guarantee certificate33**

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1 General

1.1 Precautions before use

Thank you for purchasing our growth chamber STH. You made a good choice buying a witeg product.

This manual describes the performance and usage instructions of the device, and the precautions in the handling. Please read this manual carefully, before using the product. The following warnings should be strictly adhered to:

(Warning signs are marked to attract the users attention, so please use the product safely after being well-acquainted with the meaning.)

	Warning Failure to follow the instructions may result in injury. Caution Failure to follow the instructions may result in product failure. Proximity Caution Keep clear of the product during use.
	Notification Do not disassemble or modify the product or use it for purpose other than its intended function.
	Caution: High Temperature High-temperature heating, risk of burns.
	Danger of Explosion Danger of explosion, if volatile or explosive chemicals are used.
	Protective conductor (ground)
	“OFF” on the power on/off switch
	“ON” on the power on/off switch
	Connection point of the neutral conductor
	Alternating current voltage

GROWTH CHAMBER STH-E

1.2 Product description

This product is an air circulation-type constant temperature and constant humidity chamber equipped with an optimal constant temperature and constant humidity controller. The air circulation is designed to control constant temperature and humidity inside the chamber by implementing a low-noise design. In addition, it can be remotely controlled by the installed Smart-Lab™ System. Through WiRe™ App & Web Service the device can be monitored anywhere, anytime.


This product was developed and manufactured considering the best possible ease for use and safety, and has the following features:

1. The actions on various menus have been simplified through 7" touch screen Smart-Lab™ controller.
2. In real time, unusual conditions are sent to the user through the WiRe™ App & Web Service support. Remote monitoring and control are possible.
3. The chamber is designed to uniformly maintain the internal environment through the forced circulation fan for the gasdynamic structure.
4. Through the built-in digital PID controller, temperature, humidity and intensity of illumination are precisely adjustable.
5. In order to maintain perfect tightness, the chamber is made of stainless steel.
6. The advanced insulator prevents overheating on the instruments surface as much as possible.
7. It is easy to observe the experimental status inside the chamber through the built-in tempered glass.
8. The graphic display has a high resolution to check data. Data backups and program updates can be made through the USB port.
9. The constant and accurate growth environment is composed by maintaining a relative humidity from 30% up to 95%.
10. The set value for temperature, illuminance and humidity can be used in repeated experiments by setting programs with a maximum number of 999 cycles.
11. CE certified product.

1.3 Product equipment

Growth chamber STH-155 /-305 /-420 /-800, 1 pc.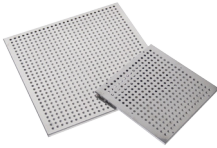

Teflon coated steel shelves, 3 pcs.

Including a copy of the operating manual

Models

- Growth chamber STH-155 : Smart-Lab™ Temperature / Humidity chamber, 155 l
- Growth chamber STH-305 : Smart-Lab™ Temperature / Humidity chamber, 305 l
- Growth chamber STH-420 : Smart-Lab™ Temperature / Humidity chamber, 420 l
- Growth chamber STH-800 : Smart-Lab™ Temperature / Humidity chamber, 800 l

2. Installation

2.1 Product installation

2.1.1 Caution of Installation

1. Please be careful when transporting the product, it is heavy.
2. Please do not lay the product on the floor or subject it to shock. This can cause damage inside the product.
3. Please install the product at a hard, flat and stable place.
4. Please do not install the product at a place where it is exposed to direct sunlight or at dangerous places.
5. To minimize the risk of short circuits, do not install the product at places where it will be subjected to moisture or organic solvents at the regulator part or inside.
6. This product was manufactured for normal operation at a rated voltage, so please check the voltage status before installing the product.
7. When installing and using the product, please use the provided power cord.


GROWTH CHAMBER STH

2.1.2 Installation Process

1. After receiving the unit and locating it in the specified location, remove the packaging.
2. Check the unit for any transport damages. If any such damage is found, please contact the deliverer of the product.
3. To prevent fire, the product should be positioned in a location where it has enough space for ventilation. The recommended distance on each side is at least 20 cm from the back wall and 100 cm from the side walls and the ceiling.
4. Place the device on a level, horizontal surface to support the devices performance without vibration.
5. Once the unit is placed at its installation location, let the unit stand for 15-30 minutes so that all the internal parts can stabilize.
6. Connect the power cord to the wall socket.
7. When installing the device, ensure the wall socket is easily accessible.
8. Do not earth to gas pipes, plastic water pipes or phone lines. This may cause electrical shocks, fire, breakdowns or the explosion of the unit.


2.1.3 Safety Symbols on the product

- Safety symbol stickers are attached to the product. For the meaning of each symbol, please refer to the chapter "Precautions Before Use".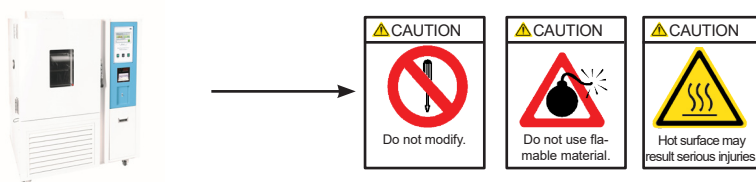


2.2 Installation of WiRe™ App & Web Service

- If you want to use the WiRe™ App & Web Service, please install it before using the product.
- The installation is progressed on the touch-screen.
- Because the wireless lan for using the internet is not installed when shipping it at the factory, please connect by finding the wireless lan stick among the components of the product.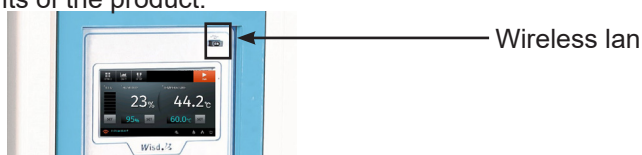

- This describes how to install WiRe™ Service.
- It only describes what must be installed to use WiRe™ Service. Please refer to chapter "3. Product Usage" p.12 for more product operation details.

- This describes the procedure on how to set the WiRe™ Service. Please follow the steps in the described order.
- Proceed to connect WiRe™ Service in the “SETTING” menu on the top left of the main screen.
- The “SETTING” menu is disabled during operation. Stop the operation by touching the “STOP” button on the top right and check if the SETTING menu is activated.

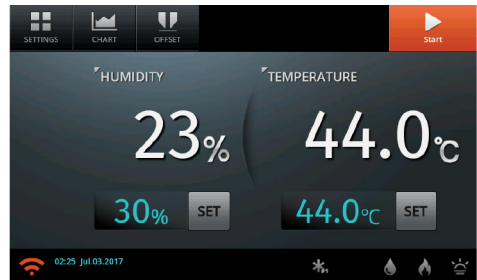- “System”, “WiRe”, “WiRe Users” are specific submenus to connect WiRe™ Service in “SETTING”. Proceed in order of description numbers as shown below.

- Proceed to set the system after pressing “System”. When you touch the return button on the top right of screen in all menu screen, the screen switches back to the previous.

I. System Settings


(1) Current Time Set

- Set the current time to enter the exact experimentation time later on. Touch the date/time box on the right.
- If you randomly change the current time supplied from the factory, please be sure to use after re-applying power to the equipment. If you have changed the time of product, its behavior may also cause problems with some functions. Therefore, if the current time is changed, be sure to use the device again after turning it off.

GROWTH CHAMBER STH


- Please use „Automatic“ only with the Check Box is unchecked and disabled.
- “Select time zone” synchronizes with the time of the country if WiRe™ Service is used at a mobile app that is connected with the device. Make sure to enter the country wherein the device is being used.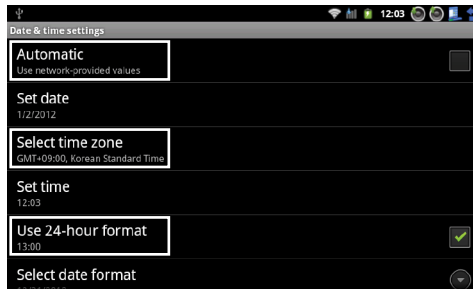

- Make sure “Use 24-hour format” is checked as shown above for a precise operation of device and never change it.
- After setting, switch back to the previous page by clicking on the blue return icon on the top right.

(2) Wifi Setup Set

- The WiRe™ Service is only possible if connected to internet. Be sure to connect to a wireless lan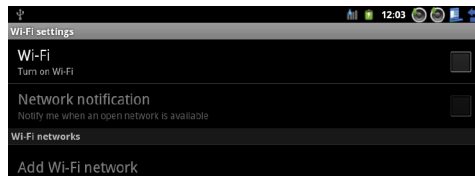

- Connect to WiFi by clicking on the check box at the Wi-Fi window, and enter the password if required. Be sure to check if the USB dongle is properly installed if the WiFi signal is not available. (Check flashing LED light)
- After connecting with a WiFi, switch back to the previous screen by clicking on the blue return icon on the top right.


WiRe Settings

- If WiRe mode button is switched to the right, WiRe™ Service is connected. This is only possible if WiFi connection is stable.


- WiRe mode is not possible if no WiFi can be connected or network is unstable. Therefore, make sure the network is stable.
- After connecting, it is possible to check serial number, model, and organization name below.

WiRe Users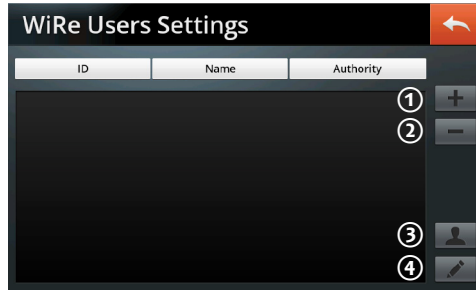


(1) Adding users

- This device can register existing user IDs, if a user is already registered on an existing Smart-Lab™ device. In such cases, a single user can check several devices from the list of registered devices that emerges when logged in with a single ID from the mobile app.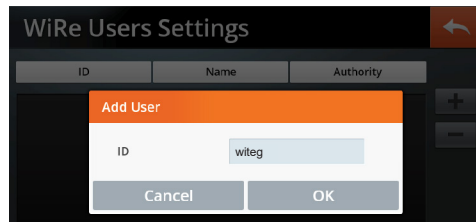

(2) Deleting users

- Users are deleted by touching the delete-button after selecting the user to delete from the user list.
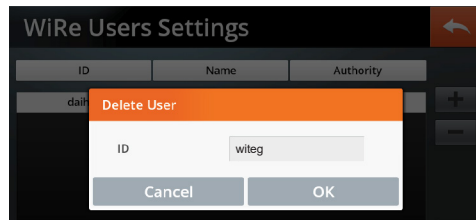


(3) Registering new users

- First-time users for Smart-Lab™ and WiRe Service™ must register user information. Enter ID, password and name.



GROWTH CHAMBER STH

- The ID must contain at least 6 and less than 20 lowercase alphabetic characters and digits. In case of a duplicate ID, less than 6 or more than 20 places or other than lower- or uppercase alphabets or numbers are entered, a pop up message will be shown alerting that this ID is not allowed.
- The password must contain at least 6 and less than 20 capital or lowercase alphabetic characters or special symbols. If not at least 6 and less than 20 capital or lowercase alphabetic characters or special symbols are entered, a pop up message not allowing registering will appear.
- Once entered, ID, password and name can not be changed.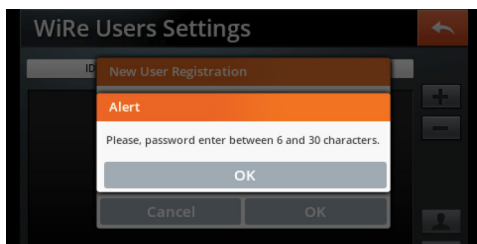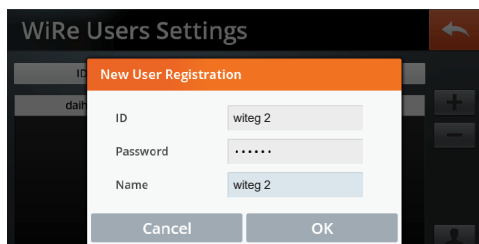

(4) Authority settings

- Set a registered user as an administrator or general user. The admin (administrator) is the only one who controls the Remote Control System. Up to 100 users (general users) can be registered and are able to monitor the status of device. (General users can not remotely control.)
- After selecting an ID from the registered user list, the authority can be set from Admin to User and vice versa.
- Since only one Admin is allowed, Admin ID other than the latest ID might be changed to a general user.

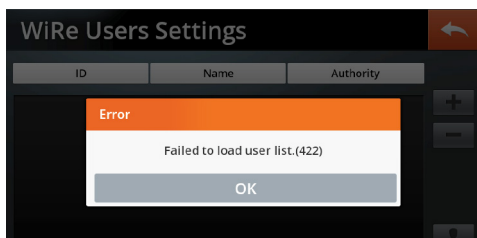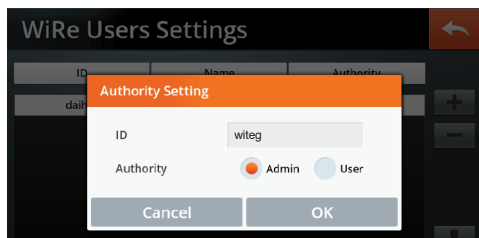- If a message that it cannot load registered user information pops up, check if WiFi connection is weak and reenter in WiRe Users menu.
- After completing the steps above, the Mobile App is connected and the WiRe™ Service is ready to use. Please refer to chapter “3. Product Usage” p. 12 for setting device functions.

2.3 Product specifications

Model		STH-155	STH-305	STH-420	STH-800
Temperature range		-20°C to 100°C			
Humidity Range		30% to 98% RH			
Compressor		7/8 HP	1.5 HP	1.5 HP	3 HP
Heater	Temperature	1.5 kW	2.0 kW	2.2 kW	2.5 kW
	Humidity	2.0 kW	3.0 kW	3.0 kW	3.0 kW
Dimensions (w x d x h)	Interior	505 x 505 x 610 mm	620 x 620 x 800 mm	700 x 700 x 820 mm	1,000 x 800 x 1,000 mm
	Exterior	1,100 x 960 x 1550 mm	1,310 x 1,080 x 1,765 mm	1,390 x 1,135 x 1,690 mm	1,570 x 1245 x 2,060 mm
Capacity		155 Lit.	305 Lit.	420 Lit.	800 Lit.
Refrigerant		CFC-free (R-404A) refrigeration system			
Temperature	Accuracy	±0.3°C			
	Stability	±0.3°C			
Humidity	Accuracy	±0.3 RH			
	Stability	±3% RH at 70%			
Heat-up Time		35 min (-20°C to 80°C)			
Cool-Down Time		45 min (+20°C to -20°C)			
Controller		Smart-Lab™ controller, 7" full touch-screen TFT LCD			
Sensor		PT100 & capacitive sensor			
Material	Internal	Stainless steel (#304)			
	External	Powder-coated steel			
Shelves		3 x perforated stainless steel shelves with adjustable position			
Safety device		High temperature and over current protector, sensor error and low level detector, leakage breaker			
Packing dimension (w x d x h)		134 x 116 x 175 cm	152 x 125 x 207 cm	160 x 130 x 202 cm	178 x 141 x 237 cm
Gross weight		353 kg	369 kg	478 kg	678 kg
Power consumption		4.7 kW	5.7 kW	6.3 kW	9 kW
		6.2 kW	7.9 kW	9.9 kW	12.7 kW
Power supply		1 Phase AC 120V, 60 Hz or 230V, 50/60Hz or 3 Phase			

GROWTH CHAMBER STH

3. Product Usage

3.1 Precautions for Use

	Please be familiar with the user manual before using this product. Operation should be performed only by qualified personnel.
	This product should be connected to a appropriate power source that it is generally used. Please use a lower power source than the regulation load permitted (AC 120, 60Hz or AC 230V, 50/60Hz).
	Place the product on a level, horizontal surface to support the weight of the product without vibration. Install the product in a place where the floor is hard and flat. Failure to do so can cause accidents by negligence.

Warnings

- Please observe the following precautions.

	This product should be operated in a room with temperature maintained between 5°C and 40°C. Long-term use in an environment that exceeds the allowable range may cause a malfunction.
	Please do not damage or distress the power cord. (Please do not hurt on the power cord, or process it, or force bend it, or pull it, and tie it.) Also, if heavy objects are placed over the power cord, or the power cord is inserted in between objects, the power cord can be damaged and may cause fire or electric shock.
	Please do not directly spray water on the product or wash it with water, to avoid short circuits or electric shock.
	Please do not touch or manipulate the electrical parts such as the power plug with wet hands electric shock.
	Please do not operate the product by manipulation of the plug to avoid electric shock or short circuits.
	Do not install the product where it can get damp or in a dusty environment. It may cause overheating or short circuit. - Pollution Degree 2 -
	Do not install the product in a location at an altitude of over 2,000 meters (6,562 feet).
	Please do not use any kind of plastic container or the goods for dry purpose. The temperature around the heater mounted on the floor of the chamber is higher than the operating temperature and if the internal temperature is rapidly risen due to the malfunction of user, there may be a fire risk.

	The internal temperature rises up to 100°C, so please be careful when opening or closing door.
	The inner chamber is hot during an operation, do not touch the inner chamber without proper hand protection during operation as burns may occur.
	Do not operate this product near a heater or an air-condition.
	Clean the bottom of inner chamber and install the shelves properly before use.
	Do not get close to the chamber with your face, hands or body. It may cause an injury or burn by high temperature.
	Only operate the unit after checking the safety it.
	The temperature of the regulator SAFETY should be set up 20 ~ 30°C higher than the operating temperature. If the regulator SAFETY is set to 0, the heater does not operate.
	If unexpected sound, smell or smog is generated, pull the main plug out from the wall socket and contact the technical assets of witeg Labortechnik GmbH.
	If irradiated or contaminated samples are placed in the chamber, the warranty is voided. Do not use this product to sterilize or disinfect objects or samples.
	Do not use outdoors avoid overheating or electrical faults. The use in places where the sun continuously shines and where it can get rained on is prohibited. It may cause overheating or electrical fault of the product.
	Please do not use flammable or volatile products in the vicinity of the product to minimize the risk of fire. Please do not use the flammable spray such as lacquers and paint, or volatile, flammable chemicals and similar products in the proximity of the product. This may cause ignition due to a spark on the switch.
	Please avoid the disassembly or modification of this product by a non-authorized person to avoid serious electrical problems and to retain the warranty of witeg Labortechnik GmbH.
	This product must be connected to a properly grounded power supply.

GROWTH CHAMBER STH

Please regularly check whether dust or dirt are on the power plug and completely connect the plug. If the plug is covered with dust, or the connection is not complete, it may cause an electrical shock or fire.

Please separate the power plug from the outlet for safety when the product is unused for long periods of time to avoid overheating or ignition due to dust.

When disconnecting the power plug, it must be grasped with the hands and then pulled. Grasping only the cord and then disconnecting it may cause overheating or ignition because internal lines may be disconnected.

Insert the plug securely into an appropriate wall socket. Loose connections result in excessive heat generation to the device and may cause arcing at the connection.

If unusual noise or signs of abnormal behavior are found during use, please immediately contact the supplier or witeg Labortechnik GmbH after turning off the power supply and disconnecting the power line.

When moving the product, please move it after disconnecting the power plug from the electrical outlet, so that the power cord and the product are not damaged in order to avoid electrical shocks or ignition due to damage on the cord.

Please maintain the clean condition of this product by removing foreign matters by gently wiping with neutral detergent and a dry cloth. Please check whether foreign matter is present or not, before and after using this product.

The fluctuation of main power voltage should be lower than $\pm 10\%$ of the nominal voltage.

The maximum relative humidity is 80% at 31°C and 50% at 40°C .


If the product was transported or stored in high humidity condition, please check the packaging condition before use. If the carton or wooden box is waterlogged, please contact the deliverer or our technical service engineer. Do not operate the unit before checking with an engineer, otherwise it may cause electrical shocks or a fire.

The sound level of the buzzer is maximum 60dB in 1 meter distance when an error is detected.

	The device should be set at a well ventilated place. Make sure the holes on the side or rear surface of the body are not blocked by a wall or an object. Do not operate the device at a poor ventilated place or use closed. It may cause fire or damage on the device as the heat is not to release or the ventilation is insufficient.
	This Symbol indicates that this product is to be collected separately. The following applies only to users in European countries: This product is designated for separate collection at an appropriate collection point. Do not dispose with the household waste. This product is required to comply with the European Union's Waste Electrical & Electronic Equipment(WEEE) Directive 2002/96/EC.

- All of the circuits, electrical or mechanical parts and configurations used in this product are the technical assets of witeg Labortechnik GmbH. Only technicians of the technical support team at witeg Labortechnik GmbH or technicians approved through the education from witeg Labortechnik GmbH should be allowed to repair this product.

3.2 Names of Each Part

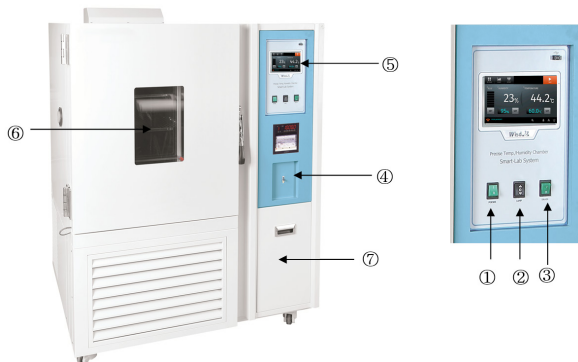No.	Names	Description
①	Power button	Power ON / OFF
②	Light button	Internal observation Lamp ON / OFF
③	Drain button	Drain ON/OFF
④	Circuit breaker	Leakage breaker
⑤	Controller	Full touch-screen TFT LCD
⑥	Window	Double glazing for internal observation
⑦	Water tank	Water tank for the humidity function


GROWTH CHAMBER STH


No.	Names	Desciption
①	Set temperature	Control and checks temperature
②	Current temperature	Indicates current temperature
③	Set humidity	Control and checks humidity
④	Current humidity	Indicates current humidity
⑤	Start/ Stop	Start and stop the operation
⑥	Program Mode control	Set programs and controls operations
⑦	Off set	Set compensating temperature values
⑧	Chart	Monitor operational status in graph
⑨	Settings	Set various functions on the device
⑩	WiFi connectivity	Indicates internet connectivity
⑪	Date	Indicates current date, time, operation hours
⑫	Operation Information	Indicates operation information
⑬	Operation ON/OFF icon	Indicates ON/OFF oft the operation with icons

	Humidity control ON		Humidity control OFF		In operation		Stopping
	Comp 1 ON		Comp 1 OFF		Heater ON		Heater OFF

Usage Instructions of water tank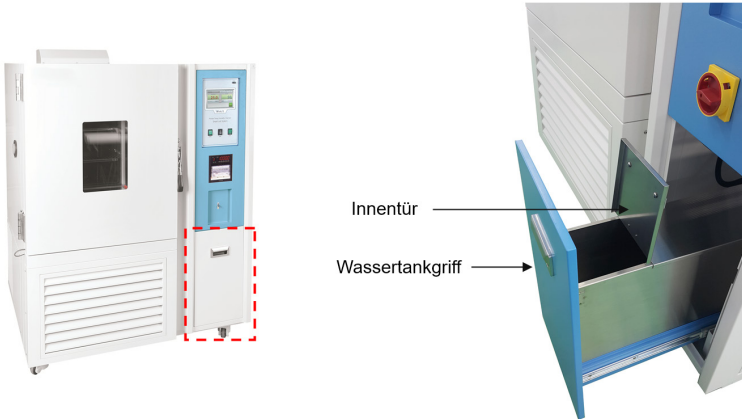

1. Please check "water tank" which locates in front of the unit
2. Pull the handle of water tank, it will be opened.
3. Please open the inner door and then fill the water until proper level.
4. Please close the inner door and then push the water tank into the unit.
5. Water tank exists to humidifier. After operation of unit, when you turn on the drain button, humidifier's water will drain to the water tank.
6. Please turn off the drain button before starting of operation. If you don't, it may cause a problem in temperature-related performance.
7. If you want to drain water from the water tank, please open the right side door and then open the manual drain valve.

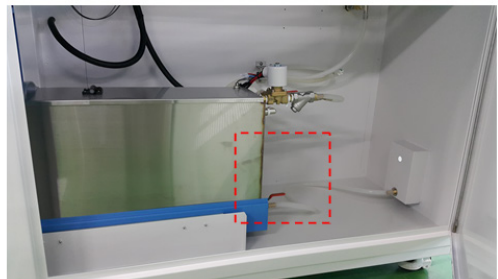
GROWTH CHAMBER STH

3.3 Usage instructions

- If turning ON the power of this product, please turn ON the main power switch after turning ON the leakage breaker located at the bottom of the rear panel.
- On the contrary, if turning OFF the power of this product, please turn OFF the leakage breaker after turning OFF the main power switch.
- Prior to perform the operation of the product, the SAFETY Knob should be set up about 20-30°C higher than the operation temperature.
- Please operate after at least one minute when you turn the power ON again.
- Please check that the internal shelf is properly installed before the first operation.
- Please set the drain button to OFF of Before the first operation.

3.3.1 Temperture setting and operation


- You can set the temperature by entering the number directly on the temperature setup screen by pressing the “SET” button on the right where the set temperature is shown.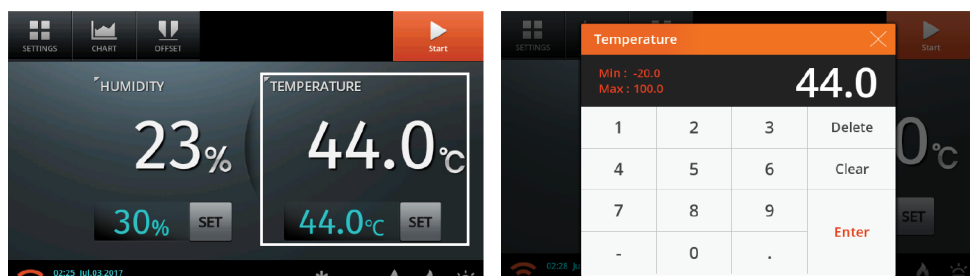

- The STH-E models can set the temperature at a 0.1°C increments and can set between -20°C to 100°C. Press “ENTER” after entering the desired temperature. It returns to the main screen when the set temperature is applied.

3.3.2 Humidity setting and operation

- You can set the Humidity by entering the number directly on the humidity setup screen by pressing the “SET” button on the right where the set temperature is shown.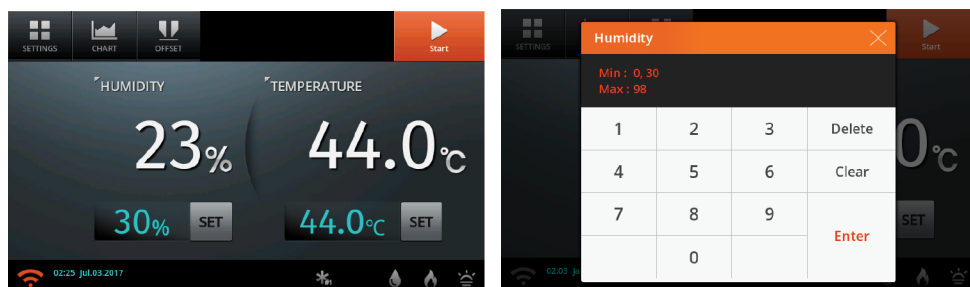

- You can set the humidity between 30% to 98%. Press “ENTER” after entering the desired humidity. It returns to the main screen when the humidity is applied.
- The humidity is not displayed at a set temperature of 15°C or lower. The water in the humidification tank is taken from the water tank.

3.3.3 Product SETTINGS



All functions on this device can be set in the SETTINGS menu. When you touch the return button on the top right of the screen in all menus, it switches back to the previous screen.

- Touch the “STOP” button on the top right to stop the operation and check if SETTINGS menu is activated since SETTING menu is disabled during operation.

- It is described in the order of the operation. Proceed with the following steps in order.
- In order to use the product's WiRe™ Service, please refer to "2.2 Installation of WiRe™ App & Web Service" on p. 6.

① Operation Settings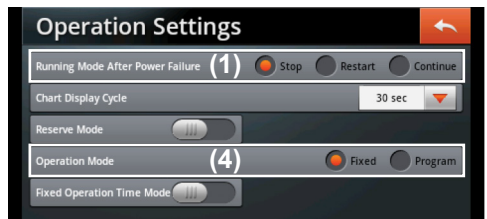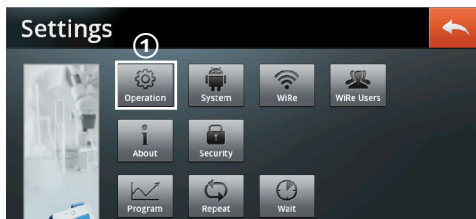

(1) Running Mode After Power Failure

- Here you can set the mode the devices operates in when the power is abruptly cut off and turned back up due to power supply issues such as power outage.
 - **Stop mode:** Is on hold at the value prior to the power outage and maintains the operation.
 - **Restart mode:** It operates depending on the (4) operation mode.
 - Fixed mode:** When the power turns back, it automatically operates at the temperature it previously operated on and the operation time on starts from 0 again.

GROWTH CHAMBER STH

Program mode: It restarts from the first segment of the previous pattern it operated on.



- **Continue mode:** It operates depending on (4) Operation Mode.

Fixed mode: When reauthorizing the power, it automatically operates at a temperature set that it previously operated on and the operation time on main screen continues its counting.

Program mode: It restarts to count from the remaining time of the first segment that stopped and restarts it.



(2) Chart display cycle

- The user can set the temperature cycle shown on the chart. Select from 30 seconds, 1 minute and 2 minutes.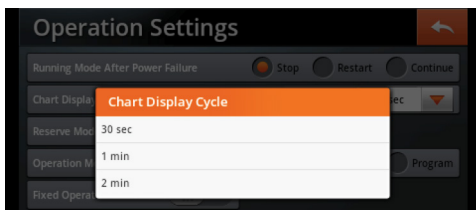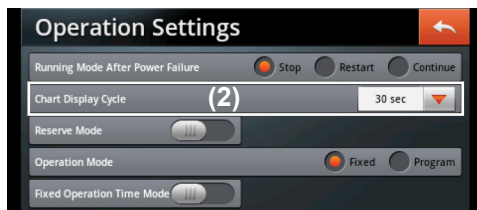

(3) Reserve mode

- Set date and time you want to operate the device and it will start at the reserved time. Minimum time required to use this mode is 5 minutes.
- It automatically turns off after being applied one time.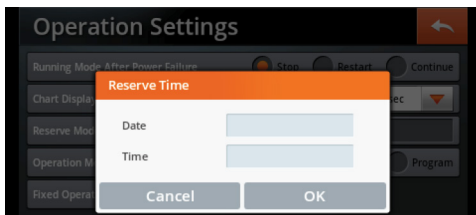
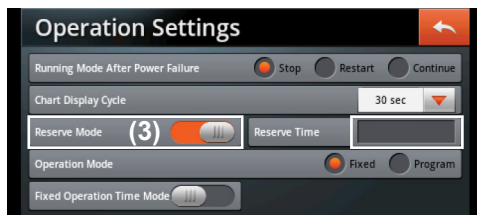


(4) Operation mode

- There are two possibilities: Fixed mode that operates within the fixed temperature range and the Program Mode that operates with various temperatures in sequence by patterns.

Fixed mode


- If this is selected, fixed operation time window is activated.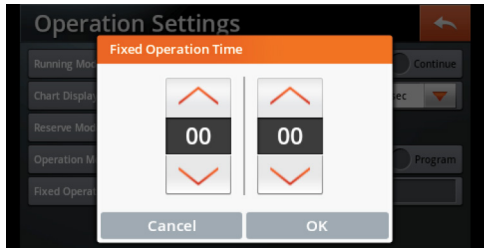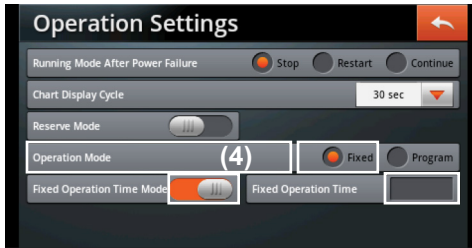

- Turn the fixed operation time mode ON and press fixed operation time to set the timer value. The device operates for the time set when you start it from the main screen.
- If fixed operation time mode is turned OFF, the set temperature keeps operating.
- the limits of the operation time is 3 minutes to 23 hours and 59 minutes.

Program Mode (in order)

- If the program is selected, the program mode window below is activated.
- In the program mode one of two different methods to reach the target temperature can be selected.
- The following describes how to set the product operation. Please refer to (4) **Program** for how to set program mode in detail.

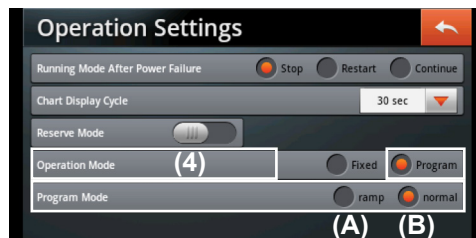A) Ramp: Set time to reach a target temperature.

Per each target temperature, the user can set a segment to reach the target temperature and a segment to maintain the target temperature. If the maintaining time for target temperature is not needed, do not make second segment for the target temperature.

B) Normal: Operates at the maximum performance to reach the target temperature.

GROWTH CHAMBER STH

② System settings

(1) Buzzer


- Select ON/OFF for the error alarm.

(2) Touch sound

- Select ON/OFF for the touch button sound.


(3) Data initialization

- Initializes operation settings and chart data stored.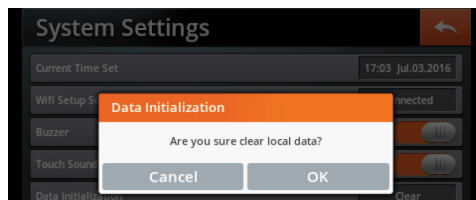


③ About

Check information on the device as serial number, model, temperature range and version of touch app.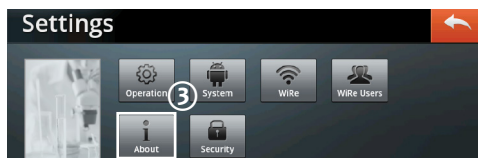

④ Security settings

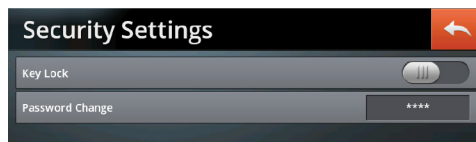
(1) Key lock

- Requests a password to access some key menu settings. Key lock is activated by sliding the button to the right.


(2) Password change

- Allows to change the password. Enter the 4 digit password stored in „Current Password“ (factory setting “0000”) and type in “New Password” and “Retype Password” to finalize the process.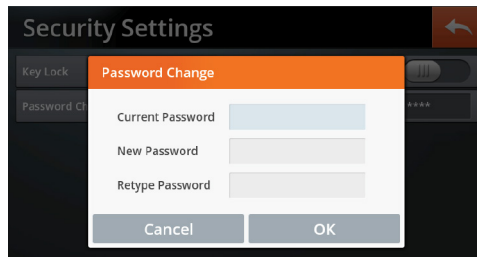


⑤ Programm (In order mode)

- You can select and store on operation program as pattern and segments.
- You can select and use one of several programs stored when using the product.
- A program is a collection of patterns and patterns are a collection of segments.

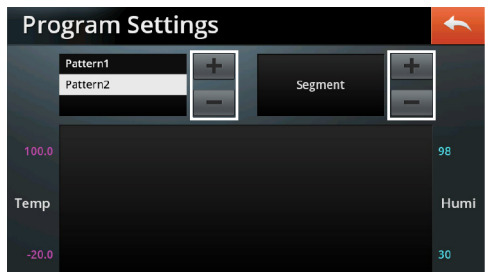


- Up to 120 patterns can be stored and a maximum of 100 segments can be stored in 1 Pattern.
- A pattern is created by using "+" button. In case you want to delete, press the number of Pattern you want to delete, verify the pattern number selected in white color and delete using the "-" button. If a pattern is deleted, the segments stored within the pattern also is deleted.
- If you want to set a segment within pattern, select the desired pattern number, create the segment by using the "+" button next to segment. To delete, press the corresponding segment number, verify the selection in white color and delete using the "-" button.
- Patterns and segments created can be immediately verified on the screen in the graph. You can enter or change values by clicking the graph.
- Enter the value of a segment after creating patterns and segments. You can verify temperature, humidity, illumination and time set for the selected segment when you select the segment on the segment window and touch.



GROWTH CHAMBER STH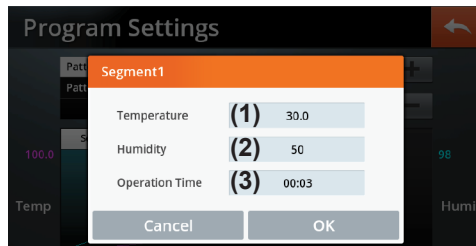

- Reset the temperature by clicking on the (1) temperature box of the segment. The Humidity can be reset by clicking on the (2) humidity box.
- Reset the timer by clicking on the (3) operation time box.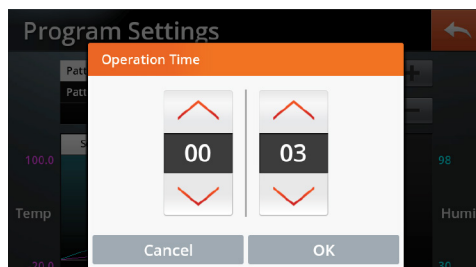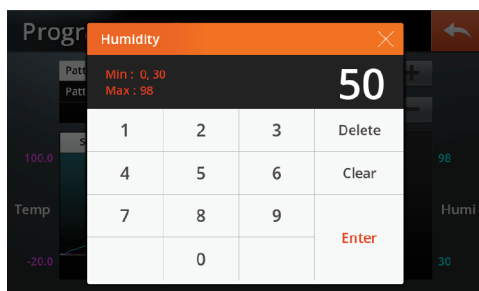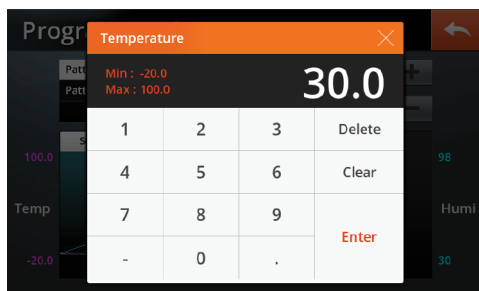

- Press OK after entering the value set to immediately verify the changes on the graph.
- The display of the segment graph depends on the operation method selected at the program mode of the operation mode.

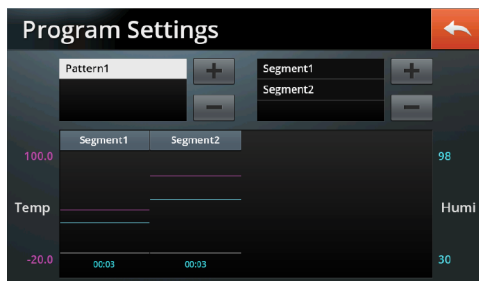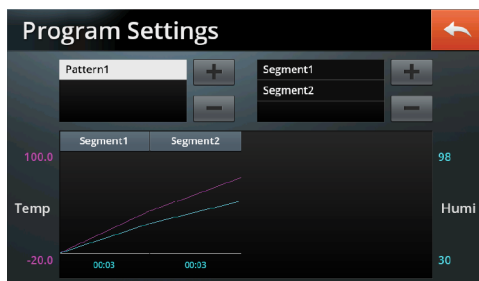


- When setting the program mode, the (1) Program menu is activated and the operation is stopped. The (2) STEP- and (3) HOLD-buttons are created in the operation of the program mode.

- (1) Program button: Enter rapidly into the Program setting screen.
- (2) Step button: Skip to next Segment during operation.
- (3) Hold button: Temporarily stops the operation.

⑥ Repeat

- Set the repeat of the operation within the selected pattern by entering the repetition after selecting the pattern from the program settings screen.
- If a pattern is created from the program settings, you can set and store the repeat mode regardless of the operation mode. However, when operating in the program mode, you must turn the repeat mode ON.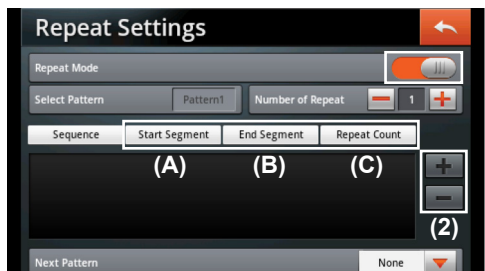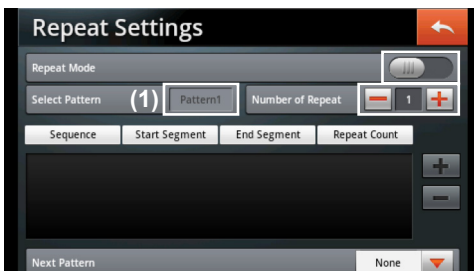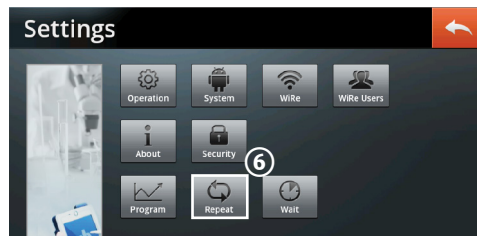

- (1) Set the number of repetition after verifying the pattern number. You can set to repeat it up to 100 times or to ∞ infinity by pressing "-" button at 1.


GROWTH CHAMBER STH

- (2) You can edit the order of the operation after setting the number of repeats. Select start and end segment by pressing the "+" button from the sequence list and set the number of repetitions between the sequences selected.



(A) Start Segment: Select first Segment for repeat operation

(B) End Segment: Select final Segment for repeat operation

(C) Segment count: Set the number of repeats from Start Segment till End Segment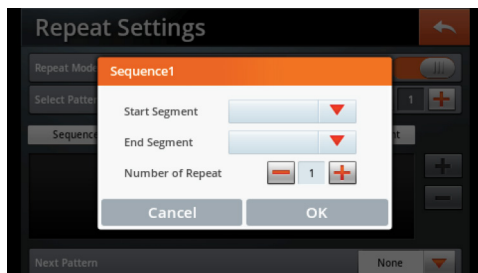




- Enter a maximum sequence of 10.
- The (3) next pattern starts to operate if the repetition of the selected pattern is completed.
- Next pattern can be selected from the stored patterns other than the pattern selected for the repeat operation. It does not allow the operation in continuation of other patterns after the repeat operation if set to none.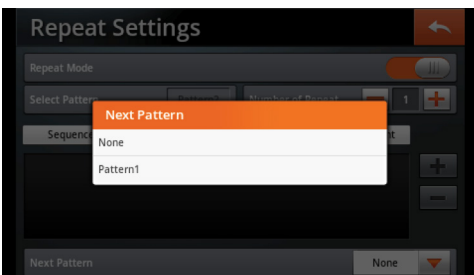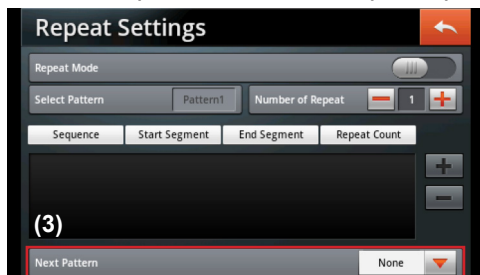


⑦ Wait

- In the program mode, if a Segment of the selected pattern has not reached the set value within time, wait time allows to wait until temperature or humidity reaches the temperature zone (set error of the target temperature/humidity $\pm$ value) without moving into the next segment.
- If the set temperature is not reached within the remaining time in ramp mode, the wait time starts.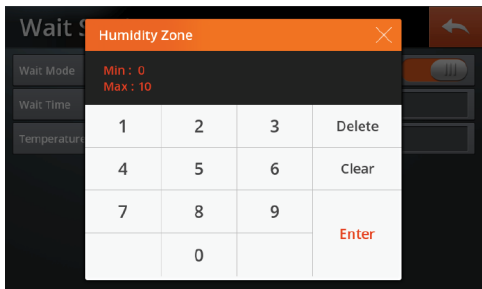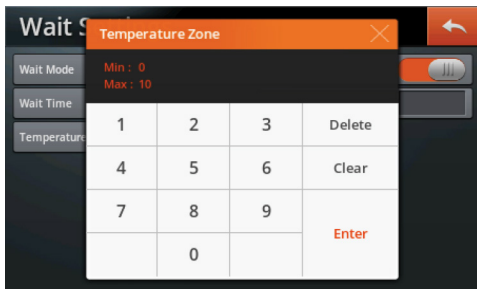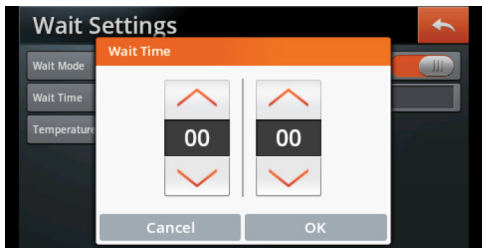


- Wait time can be set to 23 hours and 59 minutes and temperature zone can be set to $\pm 10^{\circ}\text{C}$.

3.3.4 Chart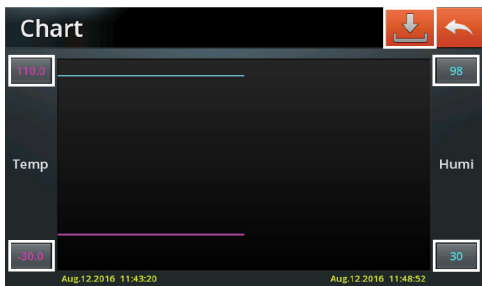


- Temperature data is stored in the internal memory every 10 seconds. A maximum of 6 months of data are kept. Please use the data backup at least once every 6 months.
- The temperature graph is displayed in purple. Temperature scope shown on graph can be changed using the set button on the low left. Maximum and minimum values can be changed within the limits shown on a pop up window.


GROWTH CHAMBER STH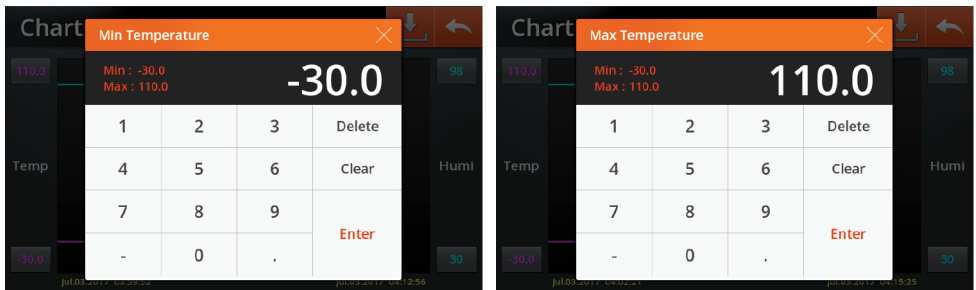


- The humidity graph is displayed in blue. Humidity scope shown on graph can be changed using the set button on the low left. Maximum and minimum values can be changed within limits shown on a pop up window.

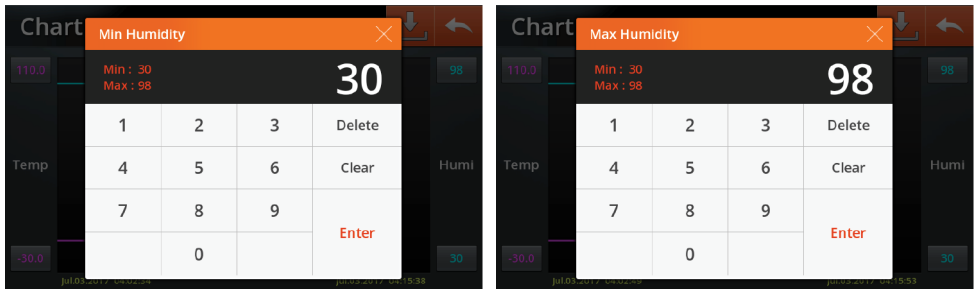- After inserting a USB flash drive into the port on top of controller, stored on the USB by pressing on the store icon on top right.
- An error message appears if the flash drive is not recognized. Press “OK”, remove and reinsert the USB, then press the store icon.


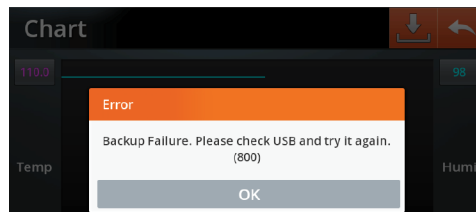


- You can see the detailed information such as date, time, temperature, error, etc. if you press a point on the graph. If you drag right and left or press two different points at the same time, detailed information of the location emerges on a detail display.
- Check the information on graph by moving at a small increment using the arrows on the detail display. The area showing detailed information is displayed in a straight line on the overall graph in yellow vertical lines.

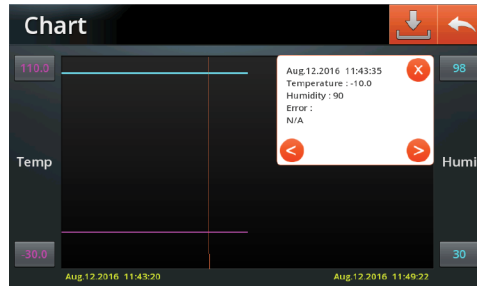3.3.5 Offset

In case a user must test the device by matching temperature, the humidity gauge can be used as a standard to test the temperature. The humidity gauge is shown on the controller. You can set the value of offset and increase or decrease the temperature displayed on the screen.

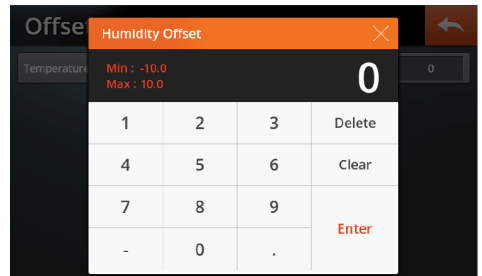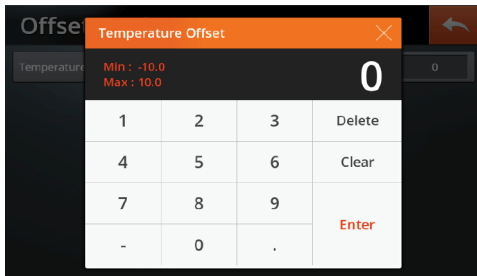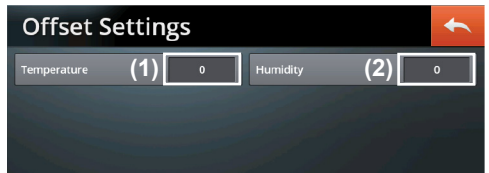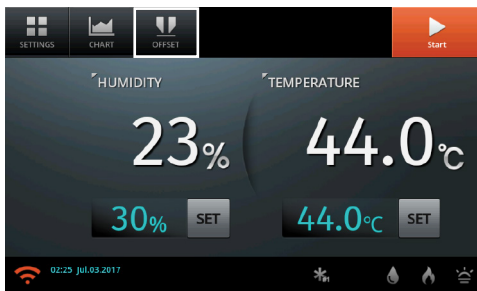- The limits of the offset values are -10.0°C to +10.0°C, and current temperature after setting the offset value displays the value including the offset.
- Once the offset value is set, it is stored and reflected in the current temperature even if you press ON/OFF button.

GROWTH CHAMBER STH

3.3.6 Error messages

- If an error happens, an error notice will be showed in the display. Click on it to see what happened in detail.

ERROR Message	Causation	Solution
Temp. Error	Display shows an unreasonable temperature value through an interference of the temperature sensor	Call the A/S Center.
Humi. Error	Display shows an unreasonable humidity value through an interference of the humidity sensor	Call the A/S Center.
Out of Control Error	Present temperature is over 15 degree (Limit can be adjusted in the User Config. manu)	Check the set value and the re-frigeration system
Water Error	Notifies when there is not enough water in the water tank.	Fill the water tank up to the mark.
Door Open	This error is generated when the door is open for more than 1 minute.	Check if the door is closed and if the problem persists when the door is closed. If so, please contact the A / S Center.
Revolution Error	Communication error problem of equipment and touch board.	Check the connection of the touch board.

* If malfunctions in the touch screen occur, please press the reset button by inserting an appropriate tool (such as a paper clip) into the reset hole, which is located on the right side of the touch-screens frame as shown below. The touch screen is rerun by pressing the reset button.

4 Products management

4.1 Accessories

Ordering information

Image	Order number	Information
 Validation (IQ, OQ) service	DH.WTH05100	Temperature & humidity recorder, "WTH5100"
	DH.WTH05105	Chart Paper, "WTH5105", 33 Days
	DH.STH31155	for "STH-155", "STHV1155"
	DH.STH31305	for "STH-305", "STHV1305"
	DH.STH31420	for "STH-420", "STHV1420"
	DH.STH31800	for "STH-800", "STHV1800"

4.2 Maintenance

This temperature and humidity chamber was manufactured with the consideration of the various situations that may occur during the operation. However, for a long-term use with stable performance, please observe the following precautions:

1. Make sure to use a rated power source of AC 120V, 60Hz or AC 230V, 50/60Hz, single phase.
2. Operate in a dry indoor location if possible and maintain the temperature between 5°C and 40°C.
3. If not used for a long period, unplug and dry it before packaging for storage.
4. If replacement parts are required due to the failures in use, only use the original parts supplied by witeg Labortechnik GmbH.
5. Failures occurred from using its outer limits cannot be repaired.
6. Authorized power sources must be used.
7. Do not touch power plugs and electrical devices with wet hands.
8. Do not put conductive or flammable materials into the vents on the outside of product for this may cause fire or electric shocks.
9. Wipe with soft towels after unplugging and moisten a towel with a lower-temperature freezing solvent to remove dirt.

GROWTH CHAMBER STH

4.3 Troubleshooting

Situation	Confirmations and Solutions
Power does not turn ON.	• Make sure that Power Cord is plugged well. • Check that there are no problems in the power supplied. • Make sure that the circuit breaker is in the ON position. • Check the short circuit status of the fuse.
Power failure	• Check the power cord is plugged into wall socket. • Check the supplied voltage is proper and regulated.
Heating, but temperature fluctuates	• Check the currently set temperature. • Make sure that the power switch is ON. • Check that the SAFETY knob is not set to "0". If the SAFETY knob is in the "0" position, set it 20°C - 30°C higher than the service temperature.
Heats but unit does not reach the temperature	• Check if the set temperature is higher than ambient temperature (recommended minimal set temperature: 10°C higher than ambient temperature). • Problem in the sensor or main controller.
Fuse is blown OR circuit breaker is flickered.	• Check the power source of wall socket and see what other loads are on the wall circuit.

If you have any problem other than the above, please call the technical service of the manufacturer or an official agent.

5 Guarantee certificate

CERTIFICATE

of
quality tested by manufacturer
and free warranty up to 2 years
for laboratory instrument

limited warranty

1. This instrument has gone through our testing process about all functions before shipment.
2. This instrument is warranted for 2 years from purchasing date.
3. A breakdown caused by customer's improper use or natural disaster is excepted from this warranty.
4. This certification is only for international customers.

Item	Smart-Lab™ Temperature Humidity Chamber
Model No.	STH
Serial No.	
Date	
Customer	

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