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Overhead Stirrer

OS-26 Digital Pro

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■ SAFETY PRECAUTIONS

◆ Safety symbol used in the manual

Danger  The exclamation point within an equilateral triangle is intended to alert the user to the presence of important operating and maintenance instructions in the literature accompanying the appliance.	Dangerous voltage  The lightning flash with arrowhead symbol, within an equilateral triangle, is intended to alert the user to the presence of uninsulated "dangerous voltage"
Explosion hazard  If volatile and explosive chemical materials are used, the explosion may occur.	High temperature  Be careful not to use at high temperature.
Notice  If the user arbitrarily dismantles or transforms the Product, the fatal damage of the Product may be induced.	Approach Prohibition  Prohibit the user from approaching to this area.
warning  If the Product is mishandled while ignoring this symbol, the danger by which human beings are dead or severely injured may occur.	Contact Prohibition  Prohibit the contact at the moment that leads to the accident.
Caution  If the Product is mishandled while ignoring this symbol, the possibility in which human beings are injured or Property damage occurs exists.	 Observe the safety rules.
 Be careful to maintain the horizontality of the Product and not to fall to the ground.	 Earth
 Keep the Product out of the direct rays of the sun.	 OFF position of power switch  ON position of power switch  Connection point of the neutral wire of installation  AC voltage

WARNING:

TO PREVENT FIRE OR SHOCK HAZARD, DO NOT EXPOSE THIS APPLIANCE TO RAIN OR MOISTURE.

CAUTION
RISK OF ELECTRIC SHOCK
DO NOT OPEN

CAUTION: TO REDUCE THE RISK OF ELECTRIC SHOCK, DO NOT REMOVE COVER. NO USER SERVICEABLE PARTS INSIDE. REFER SERVICING TO QUALIFIED SERVICE PERSONNEL.

● Precautions at transportation and installation

- Please never tilt or lay the Product during transportation.
- Be careful not to drop when transporting the Product.
- This Product was manufactured to move normally at the rated voltage. Confirm the voltage state before the installation of Product.

- Install the Product in a place where the floor is hard and flat.
- Keep the set free from moisture, water and dust. It may cause overheating or short circuit. (Pollution Degree 2)
- Do not install the Product in a location at an altitude of over 2,000 meters. (6,562 feet).

- Avoid high temperatures: Install the Product out of the direct rays of the sun.
- When the Product is moved, separate the power plug from a wall outlet, and then move not to damage the Product. The cord damage becomes the root cause for an electric shock or an ignition.

- If the Product was moved or stored in high humidity state, identify the packaging state of the Product before use.
In case that the carton box or wooden box is soaked in water, consult with an engineer about the matter. Do not operate the equipment before the safety is confirmed. This wetting state may be the cause of an electric shock or a fire.
- When installing and using the Product, please certainly use the power cord Provided with the Products.

Precaution in use

- Please be familiar with the user manual before using this Product, and operation should be performed only by qualified personnel.
- This Product must be used within the permitted standard load resistance.
- This Product should be operated in a room with temperature maintain between 5°C and 40°C. Long-term use in an environment that exceeds the allowable range may cause a malfunction.

- Do not cut or Process the power cord or do not bend it by force or pull or tie up it. In addition, if heavy things are placed on the power cord, it may cause damage and it may become the root cause of a fire or an electric shock.

- Please do not touch or manipulate the electrical parts such as the power plug with wet hands electric shock.
- Please do not operate the Product by manipulation of the plug to avoid electric shock or short circuits.

- Do not operate the Product in the outdoors. It is not allowed to use it at the place that is exposed to the sun or at the place that is gotten in the rain. This leads the Product to overheating or electric malfunction.

- Do not use the Product on a high place or at high temperature. Do not use the Product near the heater or the wind of air conditioner.
The accident by the overheating of the equipment can occur.

- Do not operate the equipment at the place of an inflammable danger or a fire danger. This triggers the severe accident such as an explosion or a fire.

- Please do not use the flammable spray such as lacquers and paint, or volatile, flammable chemicals and its similar Product, in the Proximity of the Product. They may cause the ignition due to the spark of switch.
- If you detect any noise, which is unlike any other time, or it exhibits wrong movement symptom, immediately turn off the power of the Product and separate the power cable from the body.
- Please regularly check whether dust or dirt are on the power plug and completely connect the plug. If the plug is covered with dust, or the connection is not complete, it will cause an electrical shock or fire.
- Please separate the power plug from the outlet for safety when the Product is unused for long periods of time to avoid overheating or ignition due to dust.
- Please handle the power cord carefully. Hold the plug when unplugging the cord
- The fluctuation of main power voltage should be lower than +/-10% of the nominal voltage.
- The maximum relative humidity is 80% at 31°C and 50% at 40°C.
- Please maintain the clean condition of this Product by removing the foreign matter while gently wiping with the neutral detergent and a dry cloth after checking whether the foreign matter present or not, before and after using this Product.

■ INTRODUCTION

Thank you for purchasing the OS-26 Digital Pro Overhead Stirrer.

witeg Labortechnik GmbH is one of the leading manufacturers and suppliers of glassware, liquid handling articles and electrical laboratory equipment.

The company was founded in 1960 for the development of glassware and today covers almost all laboratory requirements. The products fulfil the highest requirements in chemistry, pharmacy and medicine - and are sold in over 100 countries worldwide.

Our company philosophy prioritises quality for the benefit of our customers.

This product is manufactured in sub-miniature size to allow effective use in the customer's laboratory and guarantees low noise and vibration levels to ensure a comfortable laboratory environment. In addition, it has been equipped with electrical and instrumental safety precautions to ensure safe use. We thank you!

TABLE OF CONTENTS

① Before Using	28
② Features	28
③ Product Installation	29
• Composition identification	29
• How to install	30
• Product specification	32
④ Check before operation	33
• Product Structure	33
• Control Panel & Setup	34
⑤ Operation	36
• Stirring, Favorite and Timer operation	36
• Program Set and Operation	38
• Real-Time Monitoring Function	39
⑥ Troubleshooting	40
• Error identification	40
⑦ Product management	42
• Regular inspection	42
• Cleaning and storage	43
⑧ Accessory	43

① BEFORE USING

Pay attention to the following before using this unit:

- **Moving the set**

To prevent short circuits or damaged wires in the connection cords, always unplug the power cord safely.

- **Before turning the power switch on**

Check once again that all connections are Proper and that there are not Problems with the connection cords.

- **About operating manual**

Read and understand all the instructions in the user manual before using the Products.

In any case, we don't bear responsibility for the damage of Product which occurs by no observance for the requirements of user manual. The contents of user manual may be changed or improved. Use only for the intended purpose which is described in the user manual. After reading, store this manual along with the warranty in a safe place.

- **Note that illustrations in this manual may differ from the actual set for explanation purposes.**




② FEATURES

The OS-26 Digital Pro Overhead Stirrer provides Superior function for customer's convenient and Proper experimental results. This Product was developed and manufactured considering the best possible ease for use and safety, and has the following features.

- Ideal for Stirring Material of Viscosities Ranging from Light to Mid-Viscosity Fluids with High-precision measurement of torque in real time.
- Precise constant speed control is achieved through the adoption of the "**Digital PID Feedback Controller**", and the "**Smooth Start function**" prevents solution overflow and equipment damage due to sudden movements.
- Digital Feedback Control: Steady and constant Stirring Speed against Viscosity Change of a sample during Reaction.
- Supports "Reverse Motion" to create a larger vortex during operation and makes it easier to wet samples such as fine powder. In addition, the program operation function can be used to automatically switch from forward operation to reverse operation.
- High Fidelity Full Touch Control 3.5" TFT LCD Display with Jog Dial – The wide display enables users a convenient checking on current range of set values for the operation of the product.
- Provides 9 Program Settings and operation functions. Each program can create a total of 11 steps, and the operation time, speed, and forward/reverse operation method can be determined for each sept, allowing precise and convenient experiments to be completed.
- The timer can be set up to 99 hours and 59 minutes, providing a wide range of uses.
- Real-Time Monitoring Function - Monitor the stirring operation information in real time using the USB port.
- Chrome-Plated Anti-Rust Chuck and Precise Impeller holding part with Stainless Ste.
- Over-Load Protection and Over-Heat Protection Functions for Motor.
- Anti-Trembling function realize High Stability of Stirring.

③ PRODUCT INSTALLATION


Composition identification

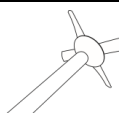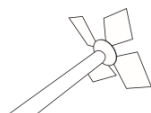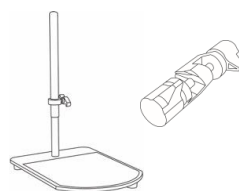Basic Components		Qty
1	 "OS-16/26/41/61" Digital Pro Overhead Stirrer Unit (Support Rod is included)	1
3	 Impeller, Propeller, 3 x bladed, Ø8 rod, Ø50 mm blade In OS-16/26	
4	 Impeller, Propeller, 4 x bladed, Ø8 rod, Ø70 mm blade In OS-41	1
5	 Impeller, Diamond, 4 x bladed, Ø8 rod, Ø90 mm blade In OS-61	
6	 Telescopic Stand & Clamp In OS-* Set package	1
6	 Operating Manual (Drill Chuck Key & Spanner are included)	1


How to install – Telescopic Stand

1 Components

Before installation, check the components below. The nut is mounted on the rod.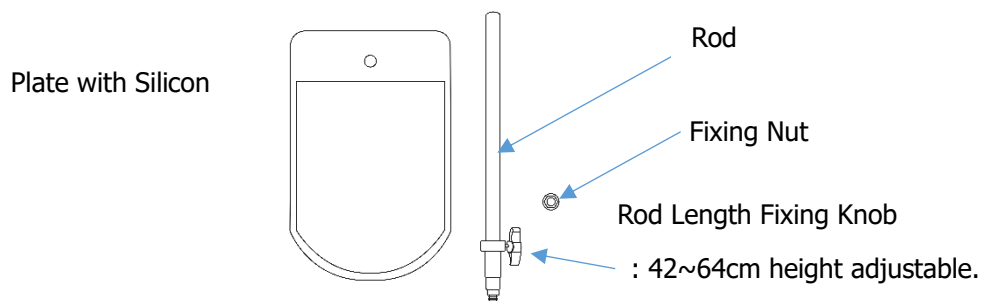





2 Installation - Disassembly method is in reverse order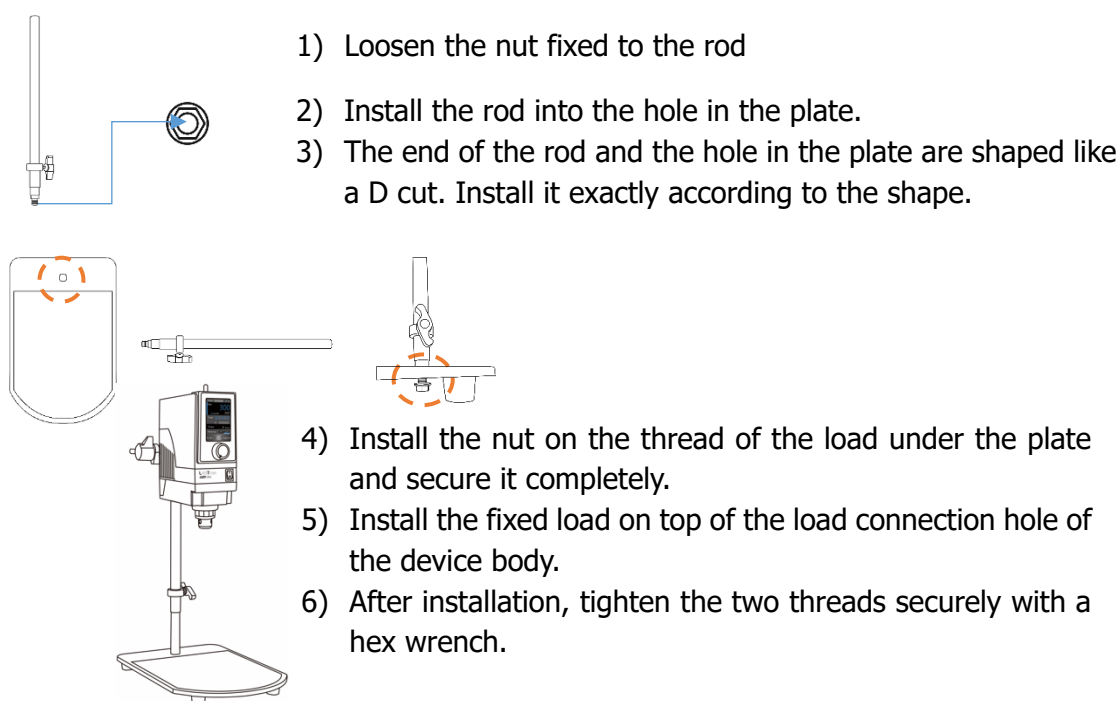

3 Caution

- The rod must be completely fixed to the plate. Otherwise, there may be a safety issue when the mounted device is in operation.
- When adjusting the length of the rod during use, be sure to stop the device, hold the bottom of the device safely with your left hand, loosen the rod length fixing knob with your right hand, and carefully adjust the length of the rod.
- After adjusting the rod length, be sure to tighten the rod length fixing knob firmly to secure it.

How to install – Standard Plate Stand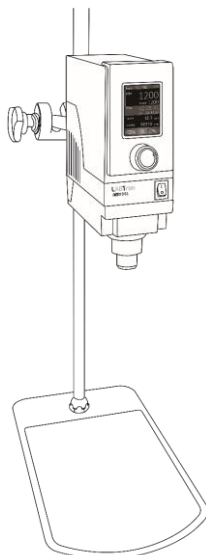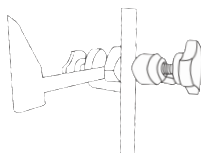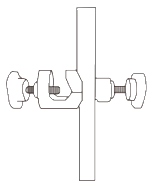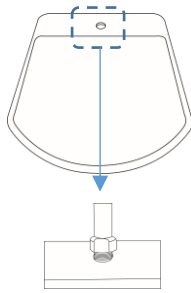


1. Place stand in to the appropriate location.
2. Connect Rod into stand and then fasten it by spanner.
3. Connect Clamp into Rod by turning it to clockwise direction.
4. After installing the support rod with the nut in the threaded hole on the back of the Product by turning it clockwise and then tighten the nut clockwise with the included spanner.
5. After connect support Rod into Clamp and screw it by turning it to clockwise direction.
6. Insert Impeller into the bottom of Drill Chuck. Impeller Length: up to 3~4cm from top of unit, if not the Impeller will be shaking when the unit is stirring.
7. After that, fasten Drill Chuck Key.
8. Complete assembly.

Product Specification

Model	OS-16 Digital Pro	OS-26 Digital Pro	OS-41 Digital Pro	OS-61 Digital Pro
Stirring Capacity	Max. 15 Lit. (H ₂ O)	Max. 25 Lit. (H ₂ O)	Max. 40 Lit. (H ₂ O)	Max. 60 Lit. (H ₂ O)
Maximum Viscosity	50,000 mPas	70,000 mPas	100,000 mPas	150,000 mPas
Maximum Torque	85 Ncm	130 Ncm	195 Ncm	293 Ncm
Gear Rate	1.5:1	2.0:1	3.0:1	4.5:1
Speed Range	50 to 3,000rpm	30 to 2,000rpm	20 to 1,200rpm	15 to 800rpm
Motor Rating (Input / Output)	200 / 160 W			
Timer	99hr 59min			
Display	High Fidelity Touch Color 3.5inch LCD Display			
Controller	Digital PID Feedback Control High Performance MCU			
Max. Impeller Shaft Diameter	10 mm			
Program Operation	9 Program, 11 Steps			
Working Direction	Forward & Reverse (Can change automatically by program operation)			
Others – Special function	Torque Measuring Function in real time Real Time Monitoring by USB port			
Permissible Temp. & Humidity	0 ~ 40°C, 80%			
Safety Device	Overload & Heat Protector for Motor Error Code Display			
Material	ABS and Aluminum			
Extension Arm Size	Ø12.7 X L250mm			
Power Consumption	280W			
Dimension (w x d x h) & Net Weight	80 X 185 X 260mm, 4.0Kg /10.5Kg			
Packing Size (w x d x h) & Gross Weight	310 X 365 X 135mm, 4.7Kg /290 X 800 X 180mm, 12.3Kg			
Power Supply & Cord Plug	1 Phase, AC 120V, 60Hz or AC 230V, 50/60Hz, with Cord/Plug			

④ Check before operation


Product structure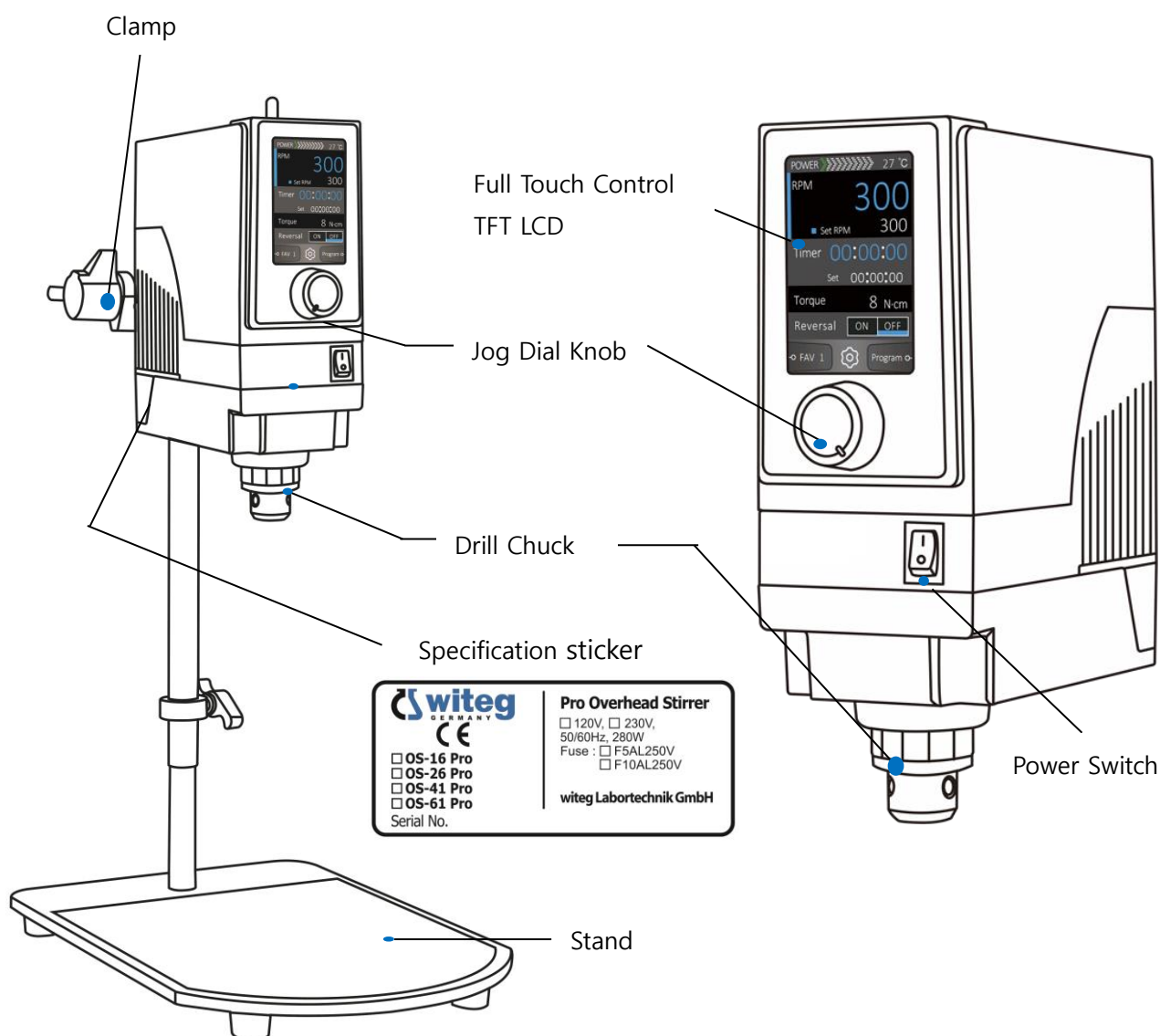

<OS-26 Digital Pro Overhead Stirrer>

- Full Touch Control TFT LCD: Viscosity and Torque measurement, RPM and Timer control.
- Jog Dial Knob: Set Target Viscosity, Stirring RPM and Timer value.
- Power Switch: Turn on / off the power.
- Specification sticker: Recording the specification and serial number of Product.

Control panel & Setup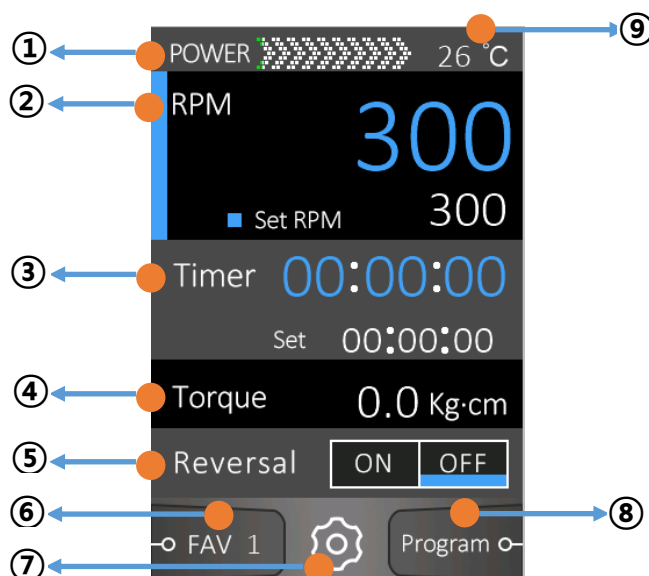

① **Motor Power Bar:** Meaning of displayed colors

- Green: Recommended of use.
- Yellow: Intermittent use, requires user's attention.
- Red: Non-recommended use (operation may be stopped for safety when used for a long time).

② **Displays the currently set speed and operating speed.**

- When you touch the RPM area, a blue bar on the left side appears at the RPM position and you can set the RPM.
-

③ **Displays the timer set by the user and the elapsed timer time.**

- When you touch the timer area, a blue bar on the left side appears at the timer location and you can set the timer.


④ **Real-time measured Torque value**

⑤ **Reverse Operation button:** Press the button to mix in the reverse direction. This can be done while in motion and helps ensure rapid wetting.

⑥ **Favorite Setting:** This is a favorite function and shows that favorite number 1 is currently being used.

⑦ **System Setting:** Set the main functions of the device.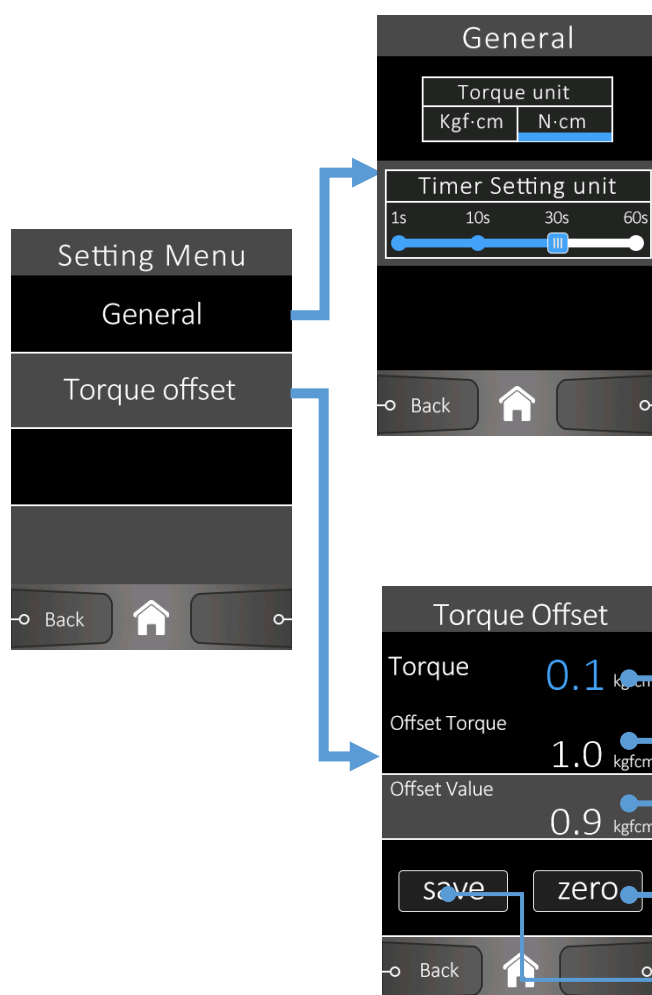

General Setting:

- Change the Torque unit.
 - a. Default: kgfcm
 - b. Main torque unit changed.
 - c. Torque offset value is initialized when the unit is changed if it is set.
- Change the Timer Setting unit.
 - a. Default: 30s
 - b. When setting the timer, you can set it to change in units of 1 second / 10 seconds / 30 seconds / 60 seconds.

Torque offset:

- When you want to arbitrarily offset the torque depending on the experimental conditions.
 - a. Current Torque
 - b. Offset applied torque value.
 - c. Set value of Offset
 - d. Offset Value is set automatically (current value is "0")
 - e. It is saved even when the power is turned off. The same effect is achieved by pushing and holding the Jog Dial Knob.


⑧ **Program Operation Function** – Provides 9 programs settings and operation functions. Each program can set a total of 11 points and the number of repetitions for that program can also be set.

⑨ **Overhead Stirrer temperature display.**

⑤ Operation

- Confirm whether the power cable is normally connected or not.
- Confirm whether or not the Impeller into Overhead Stirrer.
- Confirm whether or not the setting of sample.

Stirring and Timer Operation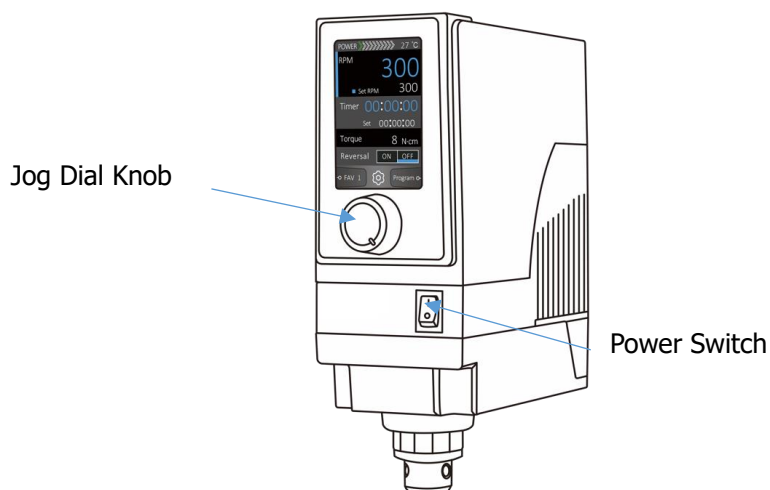

1. When the **Power Switch** is turned on, it is in the stirring RPM mode and the stirring operation is stopped at 0 RPM.

2. How to set Stirring and operation.


- Turn clockwise to increase the set RPM value and increase the speed.
- Turn counterclockwise to decrease the speed.
- If you push the Jog Dial Knob for more than 2 seconds, the current setting value is saved, and you can turn on the power again and stir with the previous setting value by simply pushing the Jog Dial Knob.

3. How to set Favorite, Timer and operation

- If you push the Jog Dial Knob for more than 2 seconds, the current setting value will be saved in the current "Favorite Setting" number. You can turn the power back on and simply push the Jog Dial Knob to perform a stirring operation with the previous setting value.
- You can also press the "Favorite Setting" button to load other settings and operate them immediately.

4. Favorite Setting: This is a favorite function and shows that favorite number 1 is currently being used.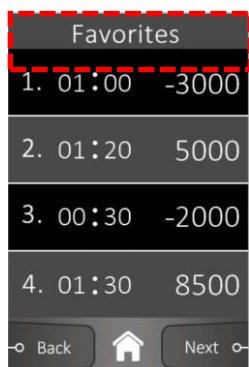


- You can store up to 8 frequently used operation times, speeds, and forward/reverse operation methods and conveniently perform experiments.
- If the reverse direction is set, the speed value is displayed as a negative value.
- Press the Next button to display the next favorites, and press the center home button to return to the main screen.

5. Timer settings and Operation


- To set and operate the timer, click the timer area on the LCD. A blue bar on the left side will appear in the timer area and you can set and operate it.
- Turn the Jog Dial Knob left and right to set the timer value and then push to start the timer.
- If you push the Jog Dial Knob for more than 2 seconds, the current setting value will be saved in the current "Favorite Setting" number. You can turn the power back on and simply push the Jog Dial Knob to operate the timer with the previous setting value.

6. Program Settings and Operation: Provides 9 programs settings and operation functions. Each program can set a total of 11 points and the number of repetitions for that program can also be set.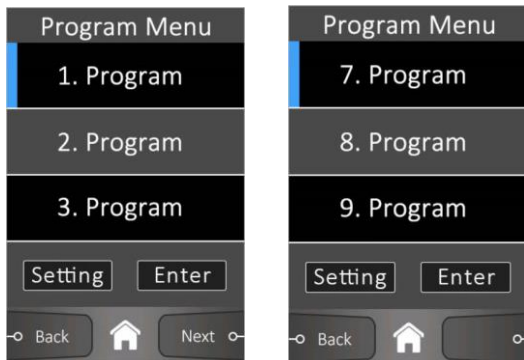

Program Menu Setting:


- If you want to set one of the 9 programs that can be set, press the "Setting" button. You can set the value of each point setting inside the program and operate it right away.
- If you want to immediately run a program that has already been saved among the 9 programs, press the "Enter" button.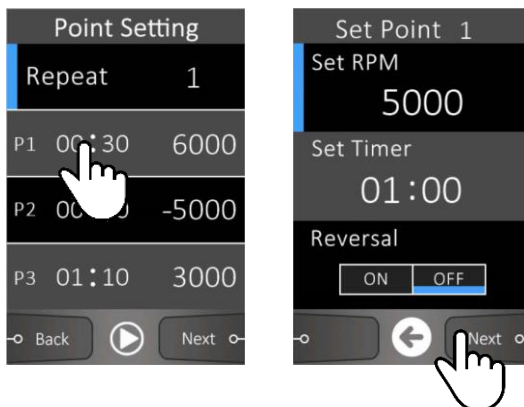

Set Point Setting:


- You can set the speed and timer operation direction starting from point 1 for a total of 11 settable points.
- Click the P1 box to set it and after completion, press the "Next" button to set the next point 2.
- If you no longer want to set the next point, press the center arrow key to return to the point setting screen.
- Click "Repeat" to set the number of repetitions of the set point settings.

➤ When setting the program mode time and repetition count, a warning screen will appear if the total time exceeds 99 hours, 59 minutes, and 59 seconds. Please adjust the settings.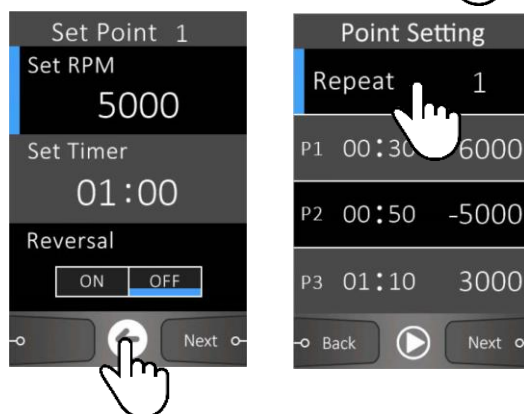

Program Operation:

- When you select a program from the program menu and press the "Enter" button or click the action icon at the bottom after completing the point settings in the program, an action screen appears with a graph depicting the program you set.
- Click the "Start" button to start the operation.
- You can go to the previous screen by pressing the Back button or go directly back to the main screen by pressing the "Home" icon.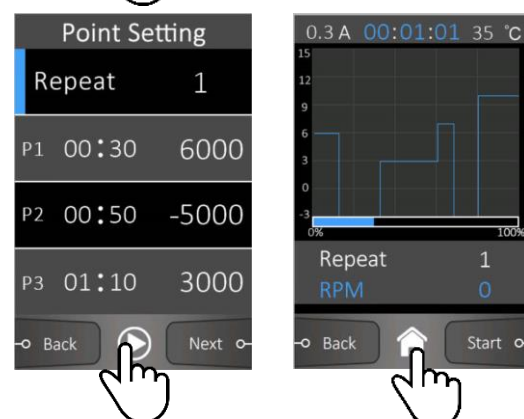


Real-Time Monitoring Function (USB Port Connection)

Digital stirrer can communicate serially with PC using USB cable (B type). Agitation status can be monitored in real time using the Provided Program, and information about agitation and equipment can be easily stored in a database.

■ Installation and usage for operation

- A. Install serial communication Program on PC: Search for "Serial Debug Assistant" app in 'MicroSoft Store' and install it.
- B. PC and Stirrer connection: Turn on the stirrer and connect the stirrer to the PC with the 'USB B to A Type' cable. Connect type B to the Stirrer and type A to PC.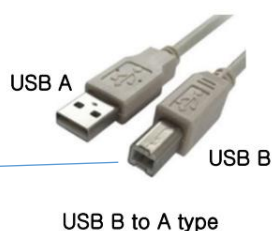

* Caution:

It is recommended not to use a USB hub, but to connect directly to the PC.

C. How to use "Serial Debug Assistant" app

- A. 'Select 'USB Serial Device' in 'Serial Port' setting.
- B. Select Baud Rate as "115200" and then click the 'Open serial port' button to connect to the stirrer and start data logging to retrieve data in real time.
- C. You can pause or restart data logging by checking 'Pause receiving display'.
- D. Click 'Save data' to save the data in a file type of ".txt" or ".csv", then convert these files to ".xlsx" and use it for data organization.
- E. After pausing data logging, you can copy the data to be saved by dragging it with the mouse and then paste it directly into 'Excel' to save it.

⑥ Troubleshooting

Error Identifications



■ ERROR Message

ERROR occurs when the stirrer is overloaded or operates abnormally. Please take Propriate action referring to the information displayed on the LCD or the manual.

Error types include 'Power Error / Current Error', 'Overheat Error', 'Driver Error', and 'System Error'.



A. Power Error / Current Error

- Occurs when a load greater than the maximum load of the Stirrer is measured.
- Check the impeller for jamming or foreign substances in the stirring solution.
- If caused by high viscosity, it is recommended to use a lower RPM or use a larger capacity stirrer.
- After solving the cause, push the Jog Dial Knob to release it.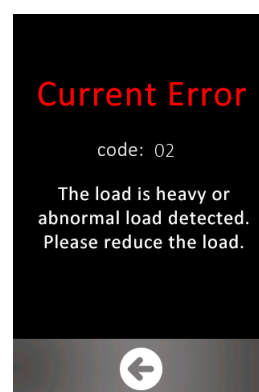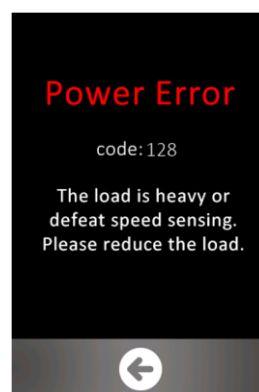


B. Overheat Error

- Occurs when the stirrer is used for a long time and overheats. (Occurs when the motor temperature on the main screen is over 60°C)
After resting (cooling) for a certain period of time recommended to use.
If the temperature is lower than 60°C after resolving the cause, the return button is created and returns to the main screen when pressed.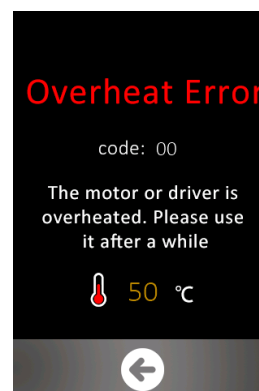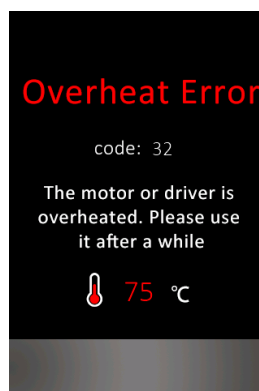


C. Motor Driver Error

- Unusable due to damage to the internal circuit of the stirrer.
- If the Problem persists even after turning the power off and on again, contact the Customer Service.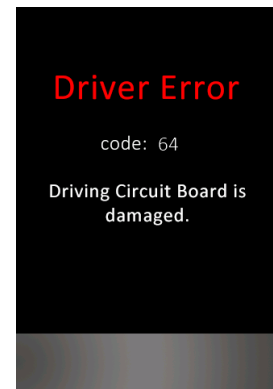

D. System Error

- Unusable due to damage to the internal circuit of the stirrer.
- If the Problem persists even after turning the power off and on again, contact the Customer Service.

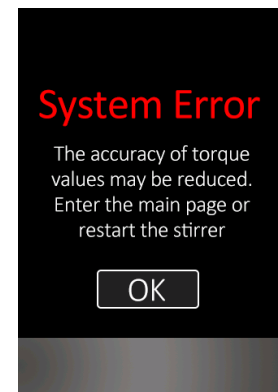■ **Guide to selecting an Overhead Stirrer model based on sample viscosity.**

Material	Viscosity	Recommend model			
Water	1	OS-16 Digital Pro	OS-26 Digital Pro	OS-41 Digital Pro	OS-61 Digital Pro
Milk	3				
Soybean oil	160				
Varnish	500				
Lacquer	1,000				
Honey	3,000				
Standard Viscosity	10,000				
Standard Viscosity	15,000				
Ketchup	30,000				
Standard Viscosity	50,000				
Mustard	70,000				
Corn Syrup	100,000				
Standard Viscosity	150,000				
Peanut Butter	250,000				

4 Product management

Regular inspection

Classifications	A period of checking time		
	Day	Week	Month
Connection state of the source of electric power	✓		
Damage of power cable	✓		
Cleaning state of power cable	✓		
Assembly state of accessories	✓		
Cleaning state of regulator	✓		
Shape preservation of regulating foot		✓	
Extent of noise		✓	
Normal control of movement mode		✓	
Normal control of regulating knob		✓	
Separate and clean the Impeller		✓	

Cleaning and storage

● Cleaning of the Product

Maintain the cleaning state of the Product by regular cleaning. The regular maintenance of the Product extends the use period of the Product and minimizes the contamination of specimen.

6. Confirm whether or not any foreign matters exist, and if yes, immediately wipe off them.
7. Do not use abrasive or detergent based on solvent.
8. If a chemical substance stains the Product, wear gloves and a mask which are composed of chemical resistance material and wipe off them with a piece of dry cloth.
9. When cleaning the Product, you must separate the power cord from the body.
10. For the part from which the contaminant is not easily erased, wipe off only the part with the towel wetting with the solvent that has low freezing point.

● Disposal of the Product



- When the warranty period of the Product is terminated, it should be disposed in accordance with the management regulation of wastes of the corresponding territory.

● Storage of the Product






3. If the Product is not used for long time, separate the power socket from the body, wrap and store it.
4. Store them in the same place not to lose the components.

5 Accessory

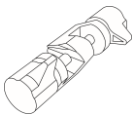Component	Information
	Standard Plate Stand, 25x37cm Plate with Rod and Silicon Pad
	Clamp, up to Ø23mm Rod

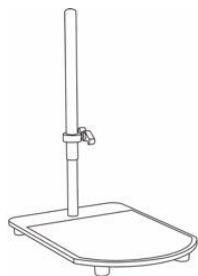Telescopic Stand for Overhead Stirrer, Homogenizer.

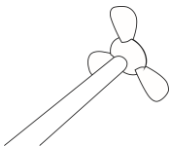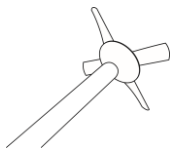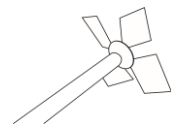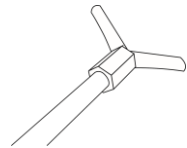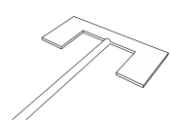	Impeller, Propeller, 3 x bladed, Ø8 rod, Ø50 mm blade
	Impeller, Propeller, 3 x bladed, Ø8 rod, Ø70 mm blade
	Impeller, Propeller, 3 x bladed, Ø10 rod, Ø60 mm blade
	Impeller, Propeller, 4 x bladed, Ø8 rod, Ø70 mm blade
	Impeller, Propeller, 4 x bladed, Ø8 rod, Ø90 mm blade
	Impeller, Propeller, 4 x bladed, Ø10 rod, Ø90 mm blade
	Impeller, Diamond, 4 x bladed, Ø8 rod, Ø90 mm blade
	Impeller, Diamond, 4 x bladed, Ø8 rod, Ø110 mm blade
	Impeller, Diamond, 4 x bladed, Ø10 rod, Ø90 mm blade
	Impeller, Diamond, 4 x bladed, Ø10 rod, Ø110 mm blade
	Impeller, Centrifugal, 2 x bladed, Ø8 rod, Ø100 mm blade
	Impeller, Centrifugal, 2 x bladed, Ø8 rod, Ø90 mm blade Teflon-coated
	Impeller, Centrifugal, 2 x bladed, Ø8 rod, Ø120 mm blade Teflon-coated
	Impeller, Anchor, 2 x bladed, Ø8 rod, Ø70 mm blade
	Impeller, Anchor, 2 x bladed, Ø8 rod, Ø90 mm blade
	Impeller, Anchor, 2 x bladed, Ø10 rod, Ø90 mm blade
	Impeller, Anchor, 2 x bladed, Ø10 rod, Ø140 mm blade

■ Free warranty service

If any malfunction due to the flaw of the Product occurs within 2 years from the purchasing date, the Product is free repaired or replaced it with the same kind Product after inquiring our head office.

■ Onerous service

- In case after a lapse of warranty service period of 2 years. In case the Product is arbitrarily remodeled or repaired.
- In case the malfunction by natural disaster. (fire, damage of sea water, gas disaster, flood damage and so on).
- In case the malfunction occurs by not observing the safety caution and warning described in the user manual.
- In case the malfunction occurs by the handling carelessness of the user.

witeg Labortechnik, GmbH

Am Bildacker 16 – 97877 Wertheim - Germany
Tel : +49 9342 9301-0 / E-Mail: info@witeg.de