**RK**

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# Contents

<b>1 Introduction .....</b>	<b>12</b>
<b>2 Safety Instructions .....</b>	<b>12</b>
2.1 Description of Symbols .....	12
2.2 Instructions .....	12
<b>3 Product description .....</b>	<b>14</b>
3.1 Key features.....	14
3.2 General survey .....	14
3.3 Display description .....	15
3.4 Term description .....	15
<b>4 Packing contents.....</b>	<b>15</b>
4.1 Standard scope of delivery.....	15
4.2 Additional accessories .....	15
<b>5 Positioning the device and commissioning.....</b>	<b>16</b>
<b>6 Job description .....</b>	<b>16</b>
6.1 Menu structure .....	16
6.2 Set speed .....	16
6.3 Set turn-off delay.....	17
6.4 Create programs .....	17
6.5 Key lock .....	17
6.6 Storage function.....	18
<b>7 Maintenance and cleaning.....</b>	<b>18</b>
<b>8 Troubleshooting Guide .....</b>	<b>18</b>
<b>9 Technical data.....</b>	<b>19</b>
<b>10 Certificate.....</b>	<b>20</b>

# 1 Introduction

Thank you for purchasing our **RK-D Programmable Rocker**. You did a good choice buying a **Wisd** product.

Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you have to follow the safety instructions of the second chapter.

Please keep this operating manual in a save place for future reference and hand it out if you consign the device to a third person.

**Witeg** reserves to modify the capsule or make amendments for the improvement of quality and efficiency without prior notice.

## 2 Safety Instructions

### 2.1 Description of Symbols

**WARNING**

Information for avoiding injuries or fatal accidents.

**ATTENTION**

Safety instruction to avoid a damage of the product.

**ATTENTION**

Information about technical parameters.

**Environment**

Information about the operating environment.

**WARNING of explosions**

Safety instruction to avoid explosions.

**Risk of injury**

Avoid injuries of your hands by following these instructions.

### 2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Connect the device only to the suitable power supply. Please have a look to the identification plate.
- Use the device only in electric circuits with ground fault circuit interrupter.
- Check the power supply cord before you stick it into the electrical outlet.
- Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.
- Do not touch the device with wet hands during the usage because this might cause an electrical shock.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.

- Make sure that the device can not slip away.
- To avoid heat congestions or fires make sure that there is enough space for air circulations.
- Avoid direct solar radiation.
- Do not use the device near highly combustible materials.
- Do not set the device outdoors.
- It is superiorly to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- To set the device in wet or dusty places could cause overheating, short circuits or fires.

- Use this device only if you read the whole safety instructions.
- This device should only used by persons who are familiarized with safety instructions made for laboratories.
- Pay attention that no liquid reaches the controller or inward of the capsule.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective work clothing during the usage.
- Pull out the power cable if you didn't want to use the device for a longer time.

- Reparations and amendments should only made by **witeg** authorized people. Contraventions exclude the warranty and may cause fatal injuries.

- Do not use this device in areas with potentially explosive atmosphere.

- Do not put your hands under the shaking plate while the shaking process runs.
- Pull out the power supply cord before you clean the device.

### 3 Product description


#### 3.1 Key features

The **RK-D Programmable Rocker** is designed for chemical, biological, medical and pharmaceutical shaking-applications. The device was developed for simplest usage and a very long life.

Key features are as follows:


- Simplest control by Jog-Shuttle knob.
- Fast commissioning.
- Long life by simply and solid construction.
- Quietly and powerful brushless DC motor for very long life.
- Over temperature and over current protection.
- Maximal tilt angle amounts 9° at standard scope delivery. 6° up on your request.
- 5-50 swings per minute.
- Second platform is mountable.
- Time, power and speed showed on illuminated LCD.
- Turn-off delay from one minute up to over 99 hours in one minute steps adjustable.
- Six customized programs with ten steps each.
- Signal sound after finished program or elapsed delay time.
- Storage functions for adjusted time and speed values.
- Key lock against inadvertent changing of set speed and timer values
- Smooth start and change of shaking speed.

#### 3.2 General survey

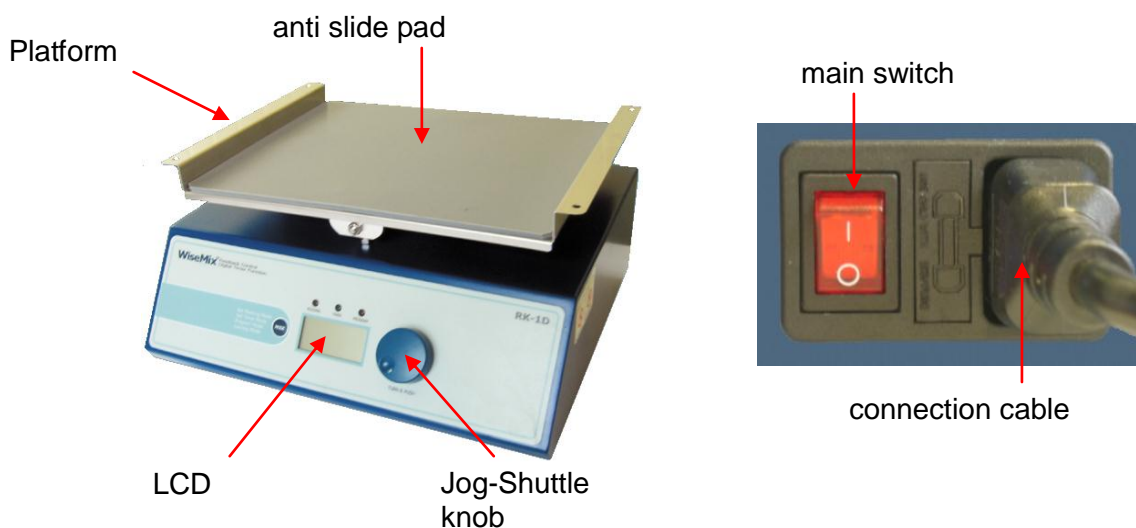### 3.3 Display description
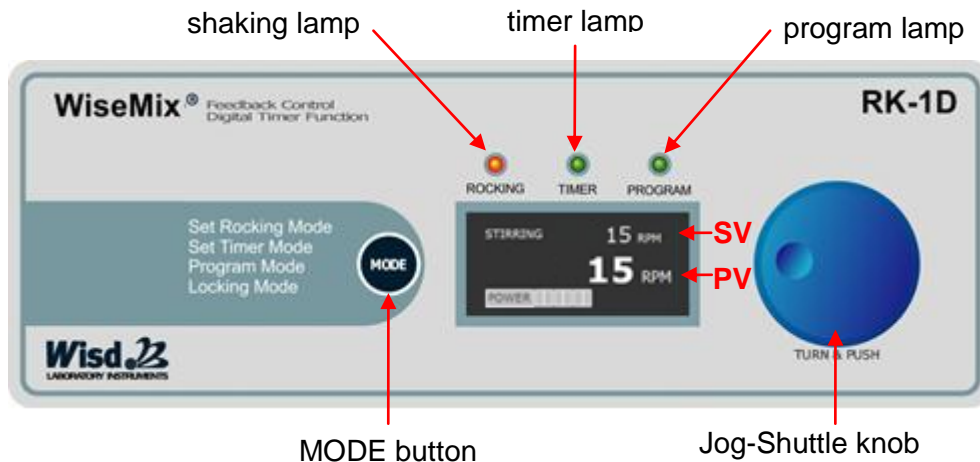
### 3.4 Term description

**SV:** set value

**PV:** present value, measured by the device

#### Main switch

- On / off switch.

#### Jog-Shuttle knob

- Change SV by turning.
- Push the knob for start and stop or probate set values.

#### MODE button

- For changing between the different menu items and storing program steps.

#### Timer lamp

- Is blinking in menu items for timer modulation
- Shines continuous in the key lock menu item if a delay time is running.

#### Shaking lamp

- Is blinking in menu items for speed modulation
- Shines continuous in the key lock menu item if the motor is running.

#### Program lamp

- Is blinking in menu items for program modulation.
- Shines continuous in the key lock menu item if a program is running.

## 4 Packing contents

### 4.1 Standard scope of delivery

quantity	DH.WRK01010	DH.WRK01020
1	<b>RK-1D</b>	<b>RK-2D</b>
1	power supply curt	
1	operating manual	
1	anti-slide pad	

### 4.2 Additional accessories


order number	denomination
DH.WRK05011	second platform for RK-1D, 290 mm x 200 mm, 4 pillars included
DH.WRK05021	second platform for RK-2D, 300 mm x 300 mm, 4 pillars included

## 5 Positioning the device and commissioning

1. Follow especially the safety instructions marked with following symbol: 
2. Remove the package carefully.
3. If there are any unexpected transport damages, please inform your forwarder and your costumer service.
4. Remove perhaps soiling and package material.
5. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatised.
6. Make sure that you follow all safety instructions.
7. Connect the device with the power outlet.
8. Switch the device on.
9. Set your values and start the process (look at point 6).

## 6 Job description

### 6.1 Menu structure

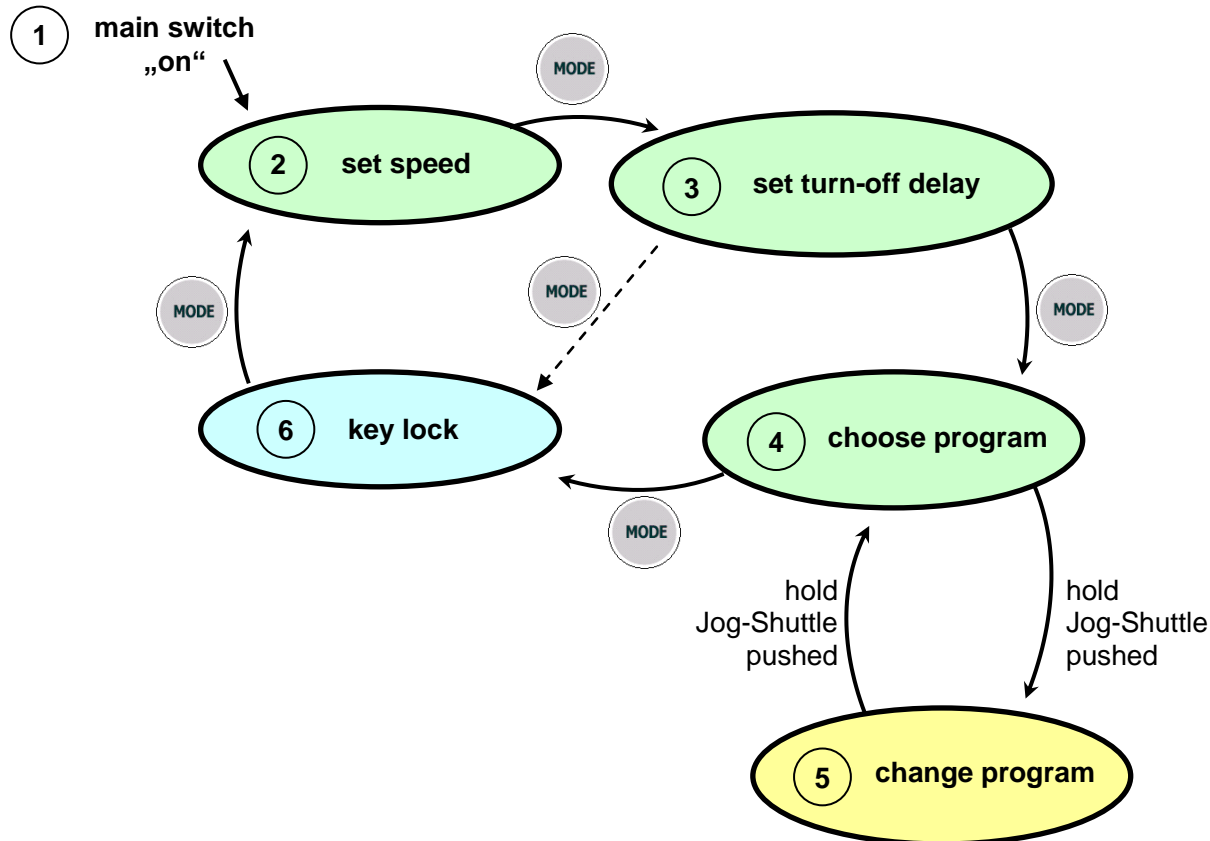### 6.2 Set speed

- First menu item after switching the device on is the menu “set speed”. The shaking lamp is blinking. You can recognize this menu at „RPM“ in SV- und PV-line.
- If there is showed “O RPM” in the SV line, there is no value stored and you can start the stirring process by turning the Jog-Shuttle knob clockwise.
- If “STIRRING” is blinking on the LCD there is a speed value stored. The saved value is showed in the SV-line and you can start the stirring process by pushing the Jog-Shuttle

knob.

- In both cases is the speed during the shaking process variable. To change the SV you only have to turn the Jog-Shuttle knob clockwise to increase the speed and counter-clockwise to reduce it.
- To stop the stirring process manually, you have to push the Jog-Shuttle knob again or decrease the speed till "0" by turning the Jog-Shuttle knob counter-clockwise.

### 6.3 Set turn-off delay

- The delay timer serves as automatically stop function. It turns off the process after the timeframe you set is elapsed.
- To reach the timer menu you have to push the MODE button in the menu item "set speed". The timer lamp is blinking.
- The delay time is adjustable from one minute till over 99 hours (00:01 - 99:59) in one minute steps.
- Turn the Jog-Shuttle knob to set your delay time. Push the Jog-Shuttle if you reached the desired value to confirm it. With the confirmation the timeframe starts elapsing.
- If you want to change the value after confirmation, you have to push the Jog-Shuttle once again. The PV initiates blinking and you can modify the value again.
- To confirm the value push the Jog-Shuttle knob yet again.

### 6.4 Create programs

- If there is no delay time running and the platform stands still, you reach the menu item "choose program" by pushing the MODE button in the menu "set turn-off delay".
- You can recognize this menu at "Pro" is showed on LCD and at the blinking program lamp.
- Turn the Jog-Shuttle knob to choose the program you want to use.
- Hold the Jog-Shuttle pushed for two seconds to reach the menu "change program" for modulation of the chosen program. "StP0" is showed on LCD now and means the first step of the program.
- First you have to set the speed by turning the Jog-Shuttle. Push the MODE button to probate and save the set value.
- Now you have to set the step duration by turning the Jog-Shuttle. Push the MODE button to probate and save the set value and jump to the next step.
- In every step you have to set speed first and then the step duration afterwards.
- It is not possible to set durations of "0", the program will only run to the previous step and stop there. But setting the speed "0" is possible.
- If you want to leave the program modulation you have to hold the Jog-Shuttle pushed for two seconds again.
- Start and stop programs by short pushing of the Jog-Shuttle.
- There is a signal sound after the program's end.

### 6.5 Key lock

- To avoid accidental changing of the set values, push the MODE button in the menu item "set turn-off delay" or in the menu "choose program" (have a look at the menu structure if you are unsure).
- In this menu item can you see the actually speed in the SV-line and the remaining time of delay function in the PV-line. The function of the Jog-Shuttle knob is deactivated.

- Press the MODE button once more to leave the key lock. Then you are in the menu item “set speed”.

## 6.6 Storage function

The **RK-D Programmable Rocker** includes a storage function for speed and delay values. That is practical if you want to work with the same parameters often. Then you need not to set them in every usage.

- For storing values you only need to hold the Jog-Shuttle knob pushed for two seconds. Successful storing is displayed by blinking of the SV for three times.
- Every stored parameter is erasable. Turn the value with the Jug-Shuttle till “0” and then hold the knob pressed for two seconds.
- If you want to store a new value it is not necessary to erase the old one.

## 7 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device and pull the power plug out of the electrical outlet.
- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- Only use original spare parts.

## 8 Troubleshooting Guide

Situation	Confirmations and Solutions
Power Failure	<ul style="list-style-type: none"> <li>✓ Check the supplied voltage is present.</li> <li>✓ Check the power connection and the power cord for proper connection and integrity.</li> <li>✓ Check if the MAIN Power Switch is “ON”.</li> <li>✓ Fuse is blown.</li> </ul>
Unit Not Stirring	<ul style="list-style-type: none"> <li>✓ Check if the MAIN Power Switch is “ON”.</li> <li>✓ Check the set RPM.</li> </ul>
ERROR 1	<ul style="list-style-type: none"> <li>✓ Errors on the Drive.</li> </ul>

**If other problems arise, or one of the above not solves problems, please contact your official agent or the manufacturer.**

**Reparations and amendments should only made by [witeg](#) authorized people. Contraventions exclude the warranty.**

## 9 Technical data

		RK-1D	RK-2D
<b>Dimension (wxdxh)</b>	<b>Platform (mm)</b>	290x200	300x300
	<b>External (mm)</b>	290x240x150	300x330x150
<b>Tilt angle (degree)</b>		±6 / ±9 (±9 as Factory Default)	
<b>Motor</b>		35 W geared DC motor	
<b>Speed range (rpm)</b>		5 - 50	
<b>Controller</b>		programmable controller with Jog-Shuttle knob (turn + push)	
<b>Program function</b>		6 programs with 10 steps each	
<b>Display</b>		digital LCD, timer, rpm and power bar-graph	
<b>Timer</b>		99hr 59min (programmable)	
<b>Safety device</b>		over load protection for motor	
<b>Others</b>		<ul style="list-style-type: none"> <li>- smooth start and change of stirring speed for stable running</li> <li>- storage function (rpm and timer),</li> <li>- key lock (Jog-Shuttle Input disabled)</li> <li>- alarm ( sensor error and timer-end )</li> </ul>	
<b>Power Supply</b>		AC 230 V, 1N~, 50/60 Hz	

## 10 Certificate

# Certificate

for

**The quality and all features were checked by the manufacturer before the shipment.**

We grant from date of purchase

**two years guaranty.**

This certificate excludes damages by natural disasters or incorrect usages by the costumer.

Please look on your account and complete following table:

Article	<b>RK Programmable Rocker</b>
Typ	
Serialno.	
Date	

**witeg Labortechnik GmbH**

Am Bildacker 16  
D-97877 Wertheim  
TEL: +49-9342-9301-0  
FAX: +49-9342-9301-77  
Email: info@witeg.de  
[www.witeg.de](http://www.witeg.de)