SWOF

Smart Ofen
Bedienungsanleitung
Smart Oven
Operating Manual
gilt für / written for
SWOF -50/ -105/ -155/ -W105/ -W155

witeg Labortechnik GmbH

(English manual at page 19)

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Contents

1 Introduction	20
2 Safety Instructions	20
2.1 Description of symbols.....	20
2.2 Instructions	20
3 Packing contents.....	22
4 Product description	22
4.1 Key features.....	22
4.2 General survey	23
4.3 Operating surface.....	23
4.4 Term description	24
5 Positioning the device and commissioning.....	24
6 Installation of WiRe™ App & Web Service	24
7 Product Usage.....	28
7.1 SETUP	28
7.2 Temperature Setting.....	31
7.3 Offset	32
7.4 Graph	32
8 Maintenance and cleaning.....	33
9 Troubleshooting Guide	34
10 Technical data.....	35
11 Certificate.....	36

1 Introduction

Thank you for purchasing our **SWOF Smart Oven**. You did a good choice buying a **Wisdom** product.

Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you have to follow the safety instructions of the second chapter.

Please keep this operating manual in a safe place for future reference and hand it out if you consign the device to a third person.

Witeg reserves to modify the capsule or make amendments for the improvement of quality and efficiency without prior notice.

2 Safety Instructions

2.1 Description of symbols

**WARNING**

Information for avoiding injuries or fatal accidents.

**ATTENTION**

Safety instruction to avoid a damage of the product.

**ATTENTION**

Information about technical parameters.

**Environment**

Information about the operating environment.

**WARNING of explosions**

Safety instruction to avoid explosions.

**ATTENTION to high temperatures**

Instructions to avoid injuries caused by high temperatures.

2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Connect the device only to the suitable power supply. Please have a look to the identification plate.

- Use the device only in electric circuits with ground fault circuit interrupter.
- Check the power supply cord before you stick it into the electrical outlet.
- Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.
- Do not touch the device with wet hands during the usage because this might cause an electrical shock.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.

- Pay attention if you open the door, hot air is coming out.
- Do not put heat-sensitive items at the device.
- Warn all attendees if you leave the device unattended.
- Do not clean the scorching device.

- Reparations and amendments should only made by **witeg** authorized people. Contraventions exclude the warranty.

- Use this device only if you read the whole safety instructions.
- This device should only used by persons who are familiarized with safety instructions made for laboratories.
- Pay attention that no liquid reach the controller or inward of the capsule.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective work clothing during the usage.
- All circuits and electrotechnical components are **witeg** owned technology. In fact of this it is necessary that only **witeg** authorized personal repair the device.
- Pull out the power cable if you didn't want to use the device for a longer time.

- It is important that the device stands up on a horizontal and firm subsurface to avoid vibrations.
- If you set the device in a site with vibrations, make sure that the device can not slip away.
- The device could get damaged if the vibrations are to strong.
- To avoid heat congestion or fires make sure that there is enough space for air circulations.
- Avoid direct solar radiation.
- Do not use the device near highly combustible materials.
- Do not set the device outdoors.

- It is superiorly to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- To set the device in wet or dusty places could cause overheating, short circuits or fires.

- Do not use this device in area with potentially explosive atmosphere.

3 Packing contents

quantum	denomination
1	SWOF Smart Oven
2	stainless Steel Shelves
1	operating manual
1	connection cable

4 Product description

4.1 Key features


The **Wisdom SWON Smart Oven** are outfitted with high performance heating mechanism that is optimized respectively for capacity of chamber, power of heating element and air circulation type (Gravity Convection or Forced Convection). The temperature displayed on the controller is accurately consistent with the actual temperature in the very center of the chamber or the average temperature in the entire area of the chamber, which maintains precisely within fluctuation specification. In addition, the temperature in the entire area of the chamber shows a very high stability keeping within a very slight variation specification in terms of temperature uniformity. Simultaneously, through the optimization of Fuzzy-PID controller, the 'heat-up time' to reach the Set Temperature after operating and the 'recovery time' to get back to the set temperature after opening and closing of the door are very short. The newly developed high performance heating mechanism has achieved high degree of temperature accuracy and uniformity in the chamber. This is achieved by the heating elements installed on three sides - that is, the left, right and bottom parts of a chamber - which directly heat the stainless-steel chamber. This either enhances its material thermal conductivity to maximum (in case of the Incubator) or the heated air from the pre-heating zone in air-jacket type that is built in the three sides outside a chamber circulates first.

Key features are as follows:

- Suitable for Drying, Baking, Conditioning, Curing, Pre-Heating and Aging.
- The Best Temp. Uniformity & Accuracy by High Performance Heating Mechanism : 3 Side-Heating.
- Optimized Air Flow by Forced Convection Mechanism.
- Ergonomic Design 4" Full Touch Screen TFT LCD (Smart-Lab™ Controller).
- WiRe™ App & Web Service.
- Automatic Data Recording.
- Self-Diagnostic Function.
- Compact New Body Design.
- Variable Program Setting for Multiple Patterns with Multiple Segments.
- Internet Connectivity with WiFi.
- Controller SW can be upgraded through USB port.

- The Recorded Data could be transferred to PC by USB memory.
- Corrosion Resistant 304 Stainless Steel Chamber.
- Over Temperature & Over Current Protection, Sensor Error Detection.

4.2 General survey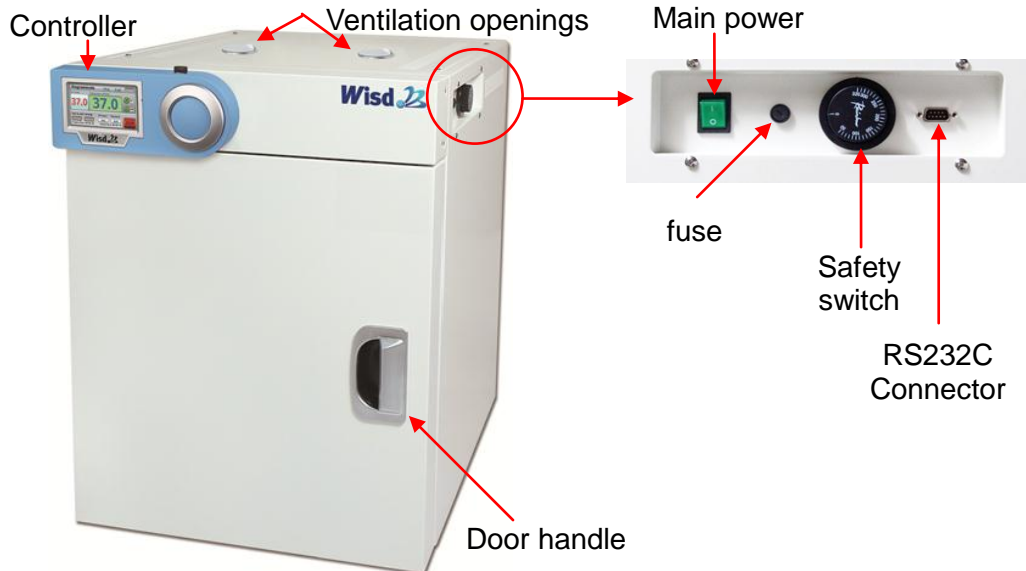

Main power

- On / off switch.

Safety switch

- The safety switch turns the heating off if your set maximal temperature is overshoot.
- Set the maximal temperature rough 30 – 40°C over the desired maximal temperature, otherwise the device maybe not reaches your set value.


Fuse

- Protect unit from High Current

RS232C Connector

- Monitoring and Controlling with PC by connecting RS232C cable.

4.3 Operating surface

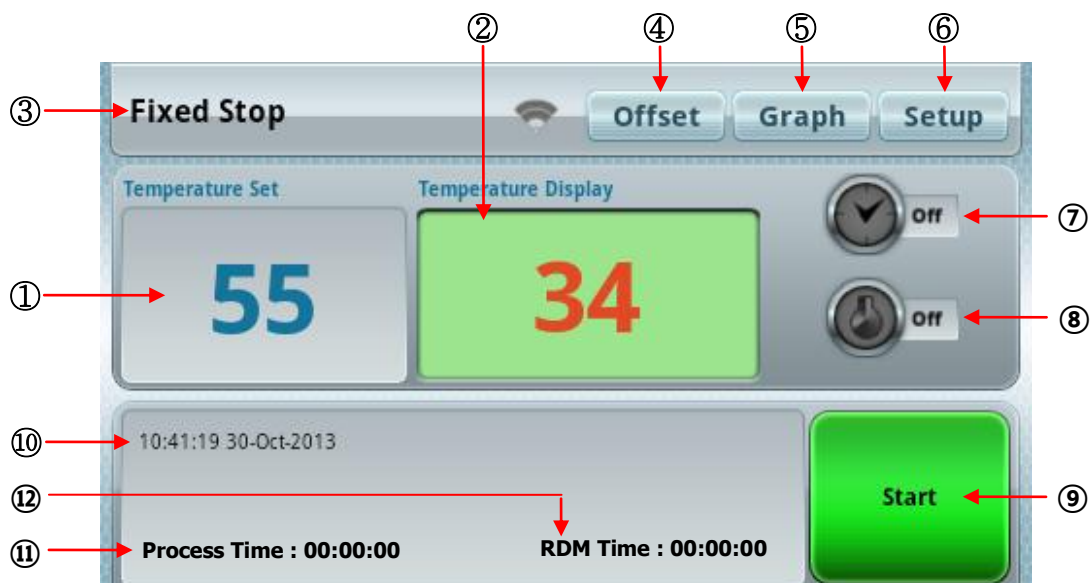4.4 Term description

NO	Names	Description
①	SET TEMP	Adjust and check the setting temperature.
②	CURR TEMP	Display the current temperature.
③	Operation Mode	Show the state of the operation mode.
④	OFF SET	Set the value of the calibration temperature.
⑤	GRAPH	Monitor the operating status by graphs.
⑥	SET UP	Set up and adjust various functions of the instrument.
⑦	Timer LED	Show the state of the timer operation through LED Lamp.
⑧	Power Failure Mode	Show the operating mode in the case of that the power is reapplied.
⑨	START/ STOP	Adjust the start and the stop of operation.
⑩	Data	Display the current date and time.
⑪	Process Time	Show the elapsed time.
⑫	RDM Time	Show the time progressed.

5 Positioning the device and commissioning

1. Follow especially the safety instructions marked with following symbol:
2. Remove the package carefully.
3. If there are any unexpected transport damages, please inform your forwarder and your costumer service.
4. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatised.
5. Make sure that you follow all safety instructions.
6. Connect the device with the power outlet.
7. Switch the device on.
8. Set your values and start the process (look at job description).


6 Installation of WiRe™ App & Web Service

- After installing this product in a static position, the user who uses WiRe™ App & Web Service, please use this product after installing WiRe Service before using the product installed.
- This installation is progressed on the regulator of front panel of the product.
- Because the Wireless LAN for using the internet is not installed in the product when shipping it at the factory, please connect by finding the wireless LAN among the components of the product.



- This description is for the installation method for users who want to use WiRe™ App & Web Service. Only matters that should be installed when use WiRe Service are written, and the detailed description on the product is listed in detail on "Instruction Manual".
- All sorts of functions of this instrument can be set through "SETUP" menu on the Touch Screen on the regulator in the front panel.

- The initial page of the description screen is the same as follows, and to press “BACK” button of the left top when the set-up is completed is switched back to the main screen.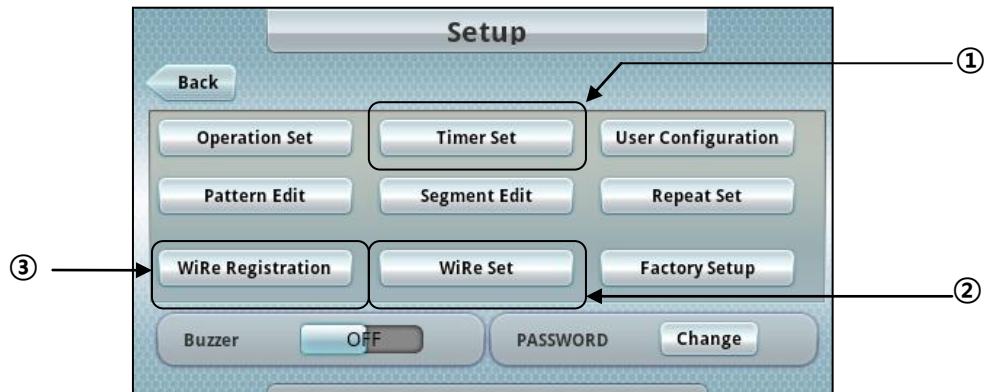


- This description is listed in order to be set when install WiRe™ App & Web Service. Please progress in order of this description.
- First, the item that you must set is Current Time. Please fill in the correct Current Time, before that the product operates through Time Set Menu.

① **Timer Set:** Set up firstly the Current Time to enter user's correct test time.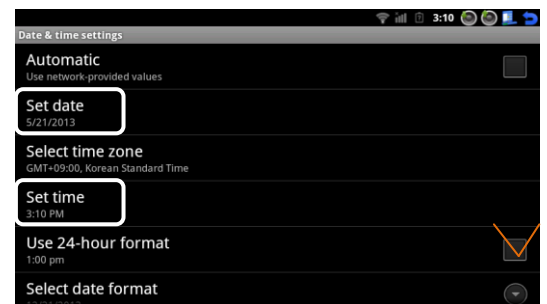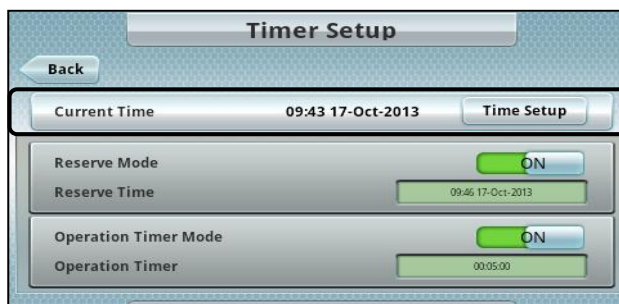

- Enter and save by entering the correct date and time by clicking "Set date" and "Set time", and then checking "Use 24-hour format" for the correct operation of the instrument.
- Please Don't check in “Automatic”, if you done it may cause unable set date and time.
- Pressing the blue returning icon at the top right of the screen return you to the previous page.
- Once you have entered the correct current date and time, connect WiRe operation.

② **WiRe SET:** Connect the WiRe™ App & Web Service, and set up its function.

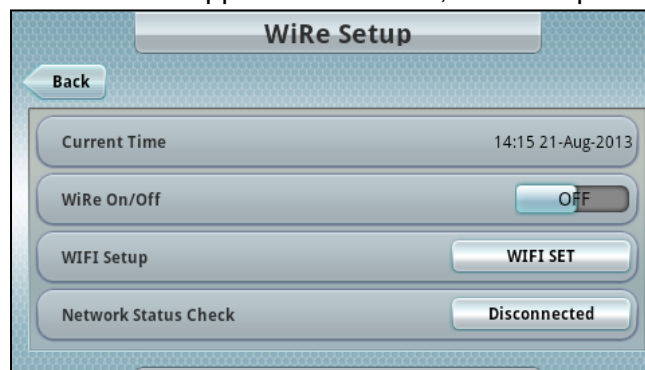- Current Time: Display the current date and time.
- Wire On/Off: Set up the connection and the disconnection to WiRe™ App & Web Service.
- WIFI Setup: Check the status of WiFi.
- Network Status Check: Change the status of network.
- Network Status Check: Displays the status connected to the network. When connecting

Wi-Fi, if the connection with the server of own company is successful, it becomes "Connect" status, and if the connection of the network or the server connection with the own company has been disconnected, displays "Disconnected".

- Sets up WiRe to On, and checks the connection status of WIFI in "WIFI SET". If the connection is completed, WiRe™App & Web Service is registered, after checking the connection with network.
- ③ **WiRe Registration:** Set up and store the record of WiRe™App & Web Service.

Factory Password: 1871

To press the green space of Touch Screen, can fill in the information.

WiRe Service Open: Store the information about the products by filling the Serial Number, Model Name, Nation, and Customer.

- Serial No and Model name have been registered when shipping the product from the factory.
- Nation information is the information for transferring the status of the equipment to near A/S Center when the problem of the product is occurred. Please enter the correct information of the corresponding country and region that are using.
- It is the item who the customer enters the used place that is using the product Please fill in the affiliated organization such as Company name and Institution name, etc. It is intended to transfer the used place of the product to A/S Center when the error is occurred in the product.
- Because the information that enters in WiRe Service Open menu is necessary to

accurately judges the product in question based on Serial No and Model name if the problem of the product is occurred, and to quickly treat by transferring the judgment information based on the information of the region and the place used to near A/S Center, you must enter the correct information.

Administrator Sign: Store the user's information about E-mail, Password, Name, and Phone.

- E-mail is the information used by ID. Please enter the mail that user mainly uses so that it is not forgotten when using.
- Password can be entered by including the numbers and characters more than four digits or less than 12 digits.
- The Create Check Box in the right of the input window is the item that should be checked when you register the Administrator ID, and Please clear the Check on the Check Box when will control other equipment with the Administration ID.
- If you want to register to different equipments with one Administrator ID, one Administrator can control different equipments, and if you want to control each the equipment by the other Administrator, Administrator ID is differently set up and each the equipment can be controlled by the Check on the Create Check Box.
- The information for Administrator can not be corrected, and when you want to modify it, it must be changed through the equipment management system of **witeg Labortechnik GmbH**.
- The Administrator generally conducts the Remote Control System as a manager.
- The Administrator manages a large number of Users, and changes the User's Active and Inactive, and can limit the use.
- If a large number of Users are set to Active and the operations are carried out at the same time, the management of the Administrator is important because an error can be occurred.

Users Sign: Many users can be stored by adding user information.

WiRe Alias: Store by filling in the information of a user who uses the WiRe™ App & Web Service. Customer can recognize different information by this from all registered units.


- Alias Sign can be registered and modified by only administrator who is registered to the Administrator Sign.
- When registering and modifying it, the Password filling in the Administrator Sign is entered and then it can be changed.

7 Product Usage

- Set the SAFETY knob to 30~40°C higher than operation temperature.
- Connect the power cord to the proper power socket and Turn on the Main Switch.
- For secure operation, the major function of controller is protected by password. Default password is "0000" and we recommend change it in "Set Up".

7.1 SETUP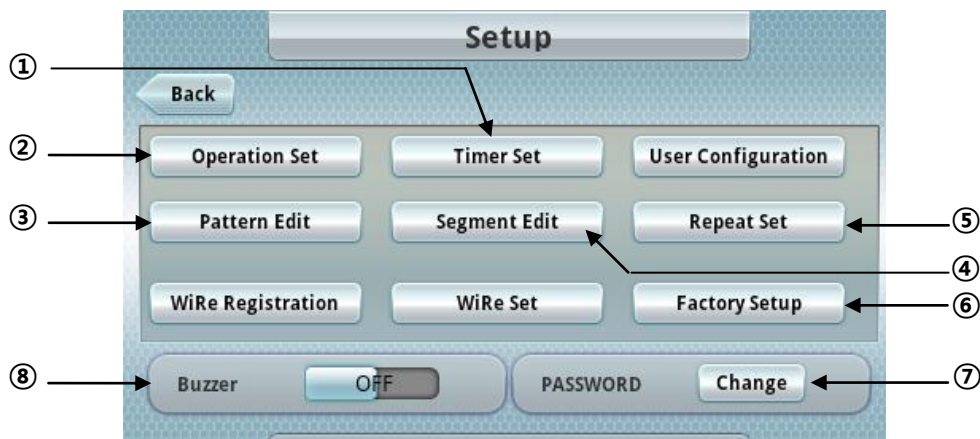

- Each button in the SET UP Mode on the left side of the Controller sets and adjusts various functions of this instrument.

① **Timer Set:** Set up the operation timer and the reservation timer.


- Current Time – Set up the current Data.

- Reserve Mode – Set up ON/OFF of the operation of reservation timer.
- Reserve Time – When Reserve Mode is set to ON, the reservation timer value can be set by pressing the green space.
- Operation Timer Mode – Set up ON/OFF of the operation timer.
- Operation Timer – When Operation Timer Mode is set to ON, the operation timer value can be set by pressing the green space.

② **Operation Set:** Select the mode of operation, and sets up the function.

- Operation Mode: It is possible to select the operation mode among both the Fixed Mode and Program Mode.
- Repeat: ON/OFF of the product are possible in only the Program Mode, and before operating, the repetition number of Pattern and Segment can be set up.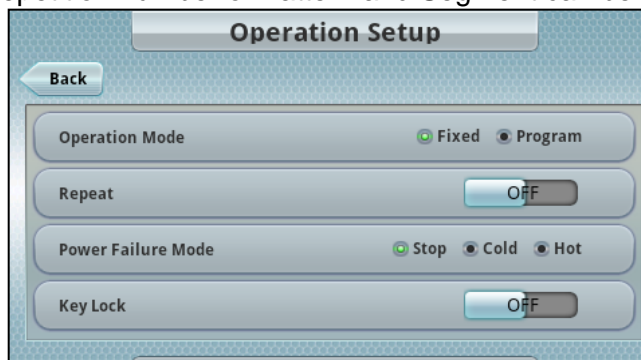

The Operation Mode can be changed by pressing “Program Stop” or “Fixed Stop” in the top of the initial screen, and its change and setup are possible in “Operation set” of “Setup”.

- Operation Mode can be selected among both the Fixed Mode that operates after setting the temperature, and the Program Mode that can operate the various stages by using patterns and segments.
- In the case that selects the Fixed Mode, the mode is activated when pressing “Start” button in the main control screen after setting “Temperature set”.
- In the case that selects the Program Mode, the pattern finally selected is indicated, and you can activate the mode by pressing the “Start” button if you want that the mode is activated.


- In the information window, the first field displays the Pattern information, and the next field displays the Segment information window.
- If you want to stop the function for the moment, press the right “Hold” button in the information window.
- If you want to allow just to go beyond to next segment, press the “Step” button.

Power Failure Mode: Set up the operating mode in the case of that the power is reapplied due to the power supply problem such as a power outage.


- **Stop:** If the power is reapplied, it becomes the standby state in the setting value before operating.
 - **Cold:** In the case of the Fixed Mode that power is reapplied, it re-operates by the setting value of the previous temperature, and if you set the timer, the timer is counted again from the beginning. In the case of the Program Mode, the timer is counted from the remaining time of the segment of the pattern that has been stopped while it has been operated previously, and it is re-operated
 - **Hot:** When power is reapplied, it re-operates by the setting value of the previous temperature in the case of the Fixed Mode, and if you set the timer, the timer counts the remaining time, and. In the case of the Program Mode, the first segment setting of the pattern that operates previously is activated again
 - **Key Lock:** ON/OFF of the Display Touch function can be set up.
- ③ **Pattern Edit:** Edit the pattern of the program operation.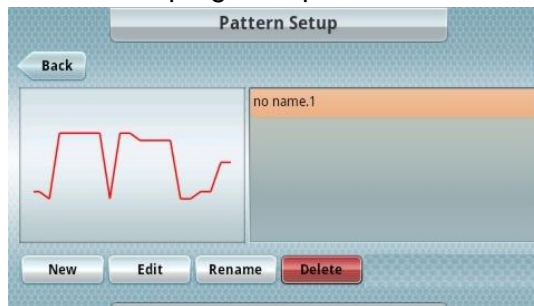


- **New:** Prepare the new Pattern.
- **Edit:** Modify the segment entered in the saved Pattern.
- **Rename:** Modify the Name of the saved Pattern.
- **Delete:** Delete by selecting the saved Pattern.

④ **Segment Edit:** Edit the Segment of Pattern.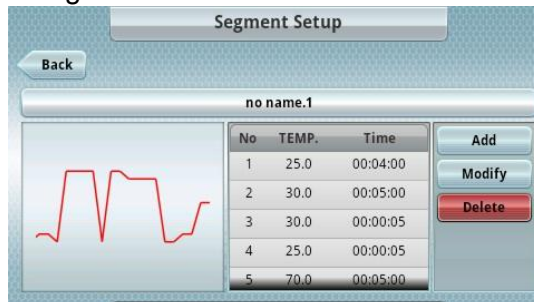

- **Add:** Prepare the new Segment by selecting the Pattern that you want.
- **Modify:** Modify the existing Segment.
- **Delete:** Select the desired Pattern and delete the Segment registered in that Pattern.

- ⑤ **Repeat Set:** When Program Mode is activated, set up the repeated numbers of the Pattern and Segment.

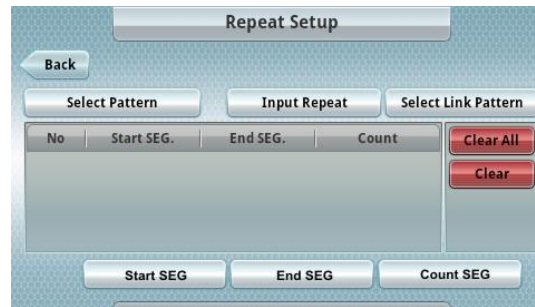
- **Select Pattern:** Select the Pattern that you want to repeat.
 - **Input Repeat:** Set up the repetition number of the Pattern.
 - **Select Link Pattern:** Select the Pattern that wants the Link.
 - **Clear:** Delete the setting repetition number, and delete all repetition number by selecting "Clear All".
 - **Segment:** Each Segment in the bottom is the beginning Segment and the ending Segment when repeatedly operate within the selected Pattern, and select the beginning and the ending by these buttons.
 - **Count:** Also set up the repetition number of the beginning and ending segment that are selected at the corresponding number.
- ① **Factory SET:** Set up the shipment information of factory, and safe operation of the product is protected by a password, and detail menu is disabled. If you need change thing, please call technical service of manufacturer or an official agent.
- ② **PASSWORD:** Menu that change/sets up a user's password.

- Enter the existing four-digit password (default value is "0000") in "Current Password", and finally save by entering a new password that you want to assign in "New Password" and "Check Password".
- ③ **BUZZER:** Allows you to stop various warning sounds. The alarm sound will ring until the user turns OFF the alarm.

7.2 Temperature Setting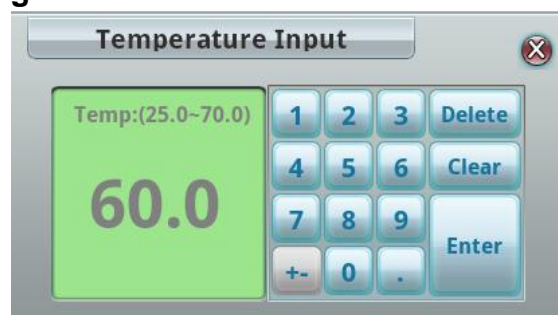

- If pressing "Temperature Set" window in the initial screen of the product, you can specify the value of the setting temperature.
- You can setup the temperature by directly entering using the number buttons, and the range of the setting temperature is adjustable in units of 1°C, and it is possible to setup up to 250°C.

7.3 Offset


- If the experiments should be performed by matching a thermometer used as a reference with the thermometer displayed on this regulator, this function is available that can reduce and add the constant values by setting the calibration value.



- If pressing "Offset" in the top of the initial screen, the correction value that you want to setup in the Offset Input window can be entered.
- The range of the compensation -10.0°C ~ +10.0°C is displayed on the top of the display window, it is possible to enter the correction value in units of 0.1°C using the number buttons.
- The display of the current temperature after setting of the compensation value is indicated as the temperature including the compensation value.
- If setting the compensation value, its value will be saved, and at ONN/OFF of the product, the current temperature that the compensation value is reflected will be displayed.

7.4 Graph

- The touch screen TFT LCD built-in in this product equips with the information function as well as the convenience for use, so the state data of operating temperature is stored in the controller for a long period.
- The Graphic display with high resolution can check the temperature record graph on the screen, and verify temperature, time and abnormal operation with simple touch, and the recorded data can be backed up to USB memory and checked on your PC.
- Data is displayed on the X-axis, and the temperatures are indicated on the Y-axis.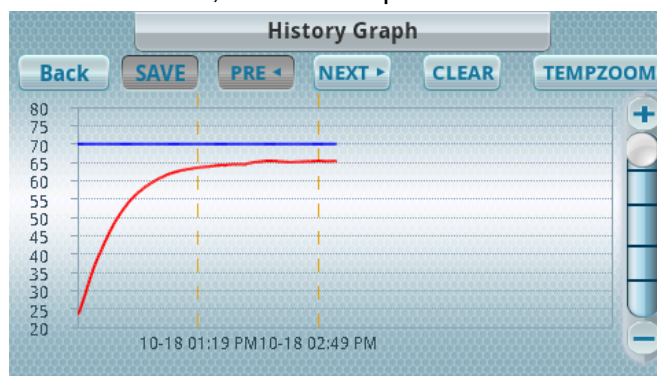

- **SAVE:** Used when store the saved data up to date in the USB memory, and at this time, the memory stick must be inserted into the port.
- **Pre/Next:** The page of graph is moved to the left (previous time) or to the right (latest hour).
- **CLEAR:** It is the button that deletes all the date stored up to now. When deleting, an alert window pops up, and if you press OK and then enter a password, all data that current equipment has will be deleted. Once the data that is deleted cannot recovered, so please note this point before deleting.
- **TEMP ZOOM:** Show by zoom in/out the desired temperature span.
- The current temperature is represented in the red graph, and the setting temperature is displayed in blue, and when you want to know more detailed status, the details window is opened if you double-click the desired field.

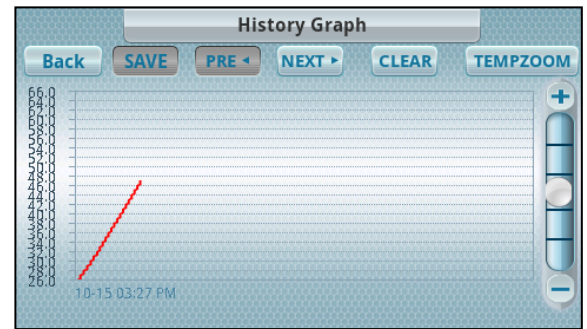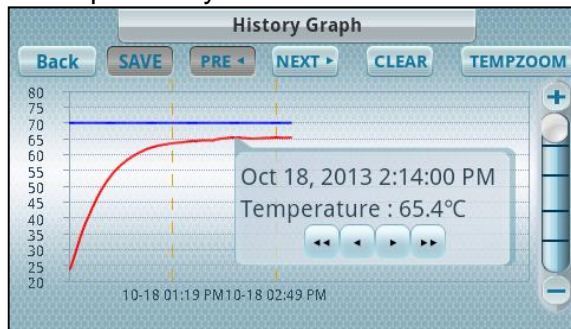8 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- The Air Duct is controlled by user's needs, and if it is unused, it should be closed the entrance and then stored in order to prevent the external contaminants entering from the entrance.
- Before cleaning the unit, disconnect the power cord from the wall socket. Otherwise, it may cause an electric shock or fire.
- To clean the unit, a neutral detergent and soft cloth is recommended. Do not use coarse cloth, strong chemicals or organic solutions.
- Do not use water, Benzene, Thinner or any alcohol for cleaning the product. It may cause discoloration, damage, an electric shock or fire.
- If you expect damage by a chemical during cleaning, call technical support team.
- Do not pour water directly into the unit. It may cause an electric shock or fire.
- During cleaning of the chamber, be careful cleaning the corners to avoid injury.

9 Troubleshooting Guide

Situation	Confirmations and Solutions
Power Failure	<ul style="list-style-type: none"> ✓ Check the supplied voltage is present. ✓ Check the power connection and the power cord for proper connection and integrity. ✓ Check if the MAIN Power Switch is "ON". ✓ Fuse is blown.
Unit Not Heating	<ul style="list-style-type: none"> ✓ Check if the MAIN Power Switch is "ON". ✓ Check the Safety switch is set higher (min. 30°C-40°C) than the set temperature (SV).
Heating, but unit not reaching temperature or temperature is fluctuated	<ul style="list-style-type: none"> ✓ Check the Set Temperature is higher than ambient temperature (recommended min set temp: 10°C higher than ambient temp). ✓ Check the Safety switch is set higher (min. 30°C-40°C) than the set temperature (SV). ✓ Check the Ventilation openings are closed. ✓ Problem in the sensor or main controller.

If other problems arise, or one of the above not solves problems, please contact **your official agent or the manufacturer.**

Reparations and amendments should only made by **witeg** authorized people. Contraventions exclude the warranty.

10 Technical data

		SWOF-50	SWOF-105	SWOF-155
			SWOF-W105	SWOF-W155
Dimension (wxdxh)	Internal (mm)	370x345x420	485x415x535	550x480x600
	External (mm)	518x647x724	653x712x882	718x777x947
Capacity		50 Lit.	105 Lit.	155 Lit.
Heater		650W	1.4kW	1.6kW
Power Consumption		674 W	1390 W	1566 W
Energy Consumption, at 100°C		117 Wh	187 Wh	204 Wh
at 150°C		230 Wh	337 Wh	398 Wh
Temp.	Range	Ambient +5 °C ~ 250 °C		
	Accuracy	±0.3°C at 100°C, ±0.5°C at 150°C		
	Uniformity	±1°C at 100°C ±1.8°C at 150°C	±1°C at 100°C ±2°C at 150°C	
Heat-up Time		25min. to 100°C 43min. to 150°C	25min. to 100°C 40min. to 150°C	
Sensor		PT100		
Control Resolution		±0.1°C		
Recovery Time (Door open 30sec)		8min. to 100°C 8min. to 150°C	6min. to 100°C 6min. to 150°C	
Controller		Smart-Lab™ Controller		
Display		4"Full Touch Screen TFT LCD		
Timer		99hr 59min, (Delay & Continuous Function Included)		
Material	Internal	Stainless Steel (#304)		
	External	Powder Coated Steel		
	Insulation	Glass Wool		
Shelves		2 Stainless Steel Shelves		
Ventilation		Stainless Steel Cap		
		Dia. 40mmx1ea	Dia. 40mmx2ea	
Circulation		Forced Air Convection Type		
Safety Device		Over Temp. and Over Current Protector, Sensor Error Detector		
Others		Storage Function (Temp., Timer), Alarm (Error Status and Timer-end)		
Net Weight (kg)		44	69	78
Packing Size (wxdxh) (mm)		622x754x807	743x819x1.068	808x884x1.133
Shipping Weight (kg)		59	79	92
Power Supply		AC 230 V, 1N~, 50/60 Hz		

11 Certificate

Certificate

for

The quality and all features were checked by the manufacturer before the shipment.

We grant from date of purchase

two years guaranty.

This certificate excludes damages by natural disasters or incorrect usages by the costumer.

Please look on your account and complete following table:

Article	SWOF Smart Oven
Typ	
Serialno.	
Date	

witeg Labortechnik GmbH

Am Bildacker 16
D-97877 Wertheim
TEL: +49-9342-9301-0
FAX: +49-9342-9301-77
Email: info@witeg.de
www.witeg.de