TITREXSOFT 2

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Titration software Operating Manual

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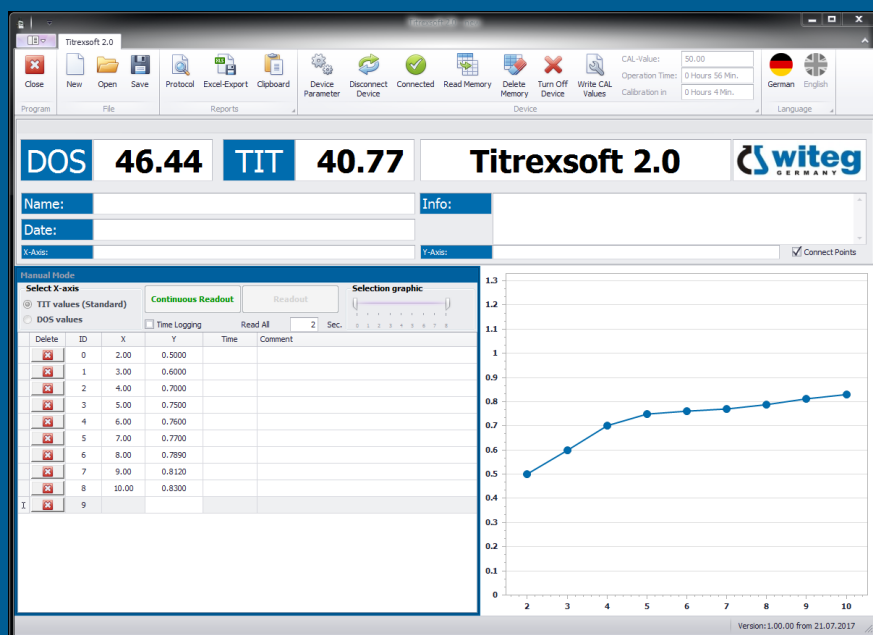
supports all manual titration applications

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Content

1. Download instructions	3
2. Installation requirements & instructions	3
3. Licensing instructions	4 - 5
4. Interface description	6 - 7
5. Project area	8
6. Report functions and saving projects	8 - 9
7. Device connection	10 - 11
7. Device features	12 - 13
9. Working modes	14 - 15
a. Manual mode	14
b. Automatic mode	14

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1. Download instructions

- Please download your Titration software at **www.witeg.de/Downloads**
- After the download is finished, extract the ZIP file into any folder on your hard drive
- To start the installation please, double-click on the extracted Setup file (EXE)

Please check the filename of the downloaded and extracted file. It should be named:

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2. Installation requirements & instructions

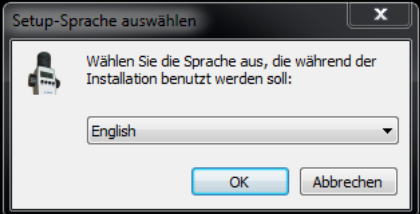
Titrexsoft requirements:

- Microsoft Windows 7 or higher
- Microsoft .NET Framework 4.0 or higher
- Screen resolution 1024x768 or higher

Requirements for connection of digital burette Titrex with Titrexsoft:

- RS232 connection cable (included in delivery) and free RS232 (COM) port on computer
- USB to RS232 adapter (on request) if no RS232 (COM) port is available

After starting the installation by double clicking on the setup file please follow the steps shown in the windows:

	<p>Choose from the following program languages:</p> <ul style="list-style-type: none">■ English■ German <p>Click on „Ok“ to get to the next window and follow the instructions shown there.</p> <p>After the installation is finished please launch the software.</p>
---	--

3. Licensing instructions

ATTENTION:

In order to use the software, you have to follow a short licensing process. You need to send some information to witeg to obtain a working license key.

The process is free of charge.

Your data will only be used for registering purposes!

License Information

To use the software you have to do the registration process.

Please fill out all relevant information under "Request Licence...". If you're using Microsoft Outlook you can send us all information with the automatically created mail. Otherwise please send a mail with all relevant information together with the Registration Key (7 digits) to license@witeg.de.

Your request will then be checked and the licence key sent to the given mail address.

This process is free of charge. The sent data will only be used for the registration process.

Request Key: 3037904

License Key: - -

Please read through the licence dialog and click on "Request license" to get to the licensing formula.

Request license key

Please fill in at least the required fields to start the request.

Your Data

Company:

Name:*

Tel.:

Email:*

serial number:

☐ No device available

Request Values

Key: 3037904

Receiver: pstockmann@witeg.de

* Required Fields

Fill out any needed information (marked with a *).

Please make sure to use a valid mail address in order to obtain your license key.

If you're using the software without a Titrex digital burette from witeg Labortechnik GmbH, please check "no device available".

Finally, click on "Start your inquiry".

For Microsoft Outlook users	For all other users
<p>After filling out any needed information, please click on „start request“. A new mail will be automatically opened which can be sent to license@witeg.de</p>	<p>After filling out any needed information please follow these instructions:</p> <ol style="list-style-type: none"> 1. If you don't have Outlook installed all information will be saved to the clipboard 2. Open a new mail in your mail software. 3. Insert copied information into the text field of your mail. <i>Make sure, that the value for "Key" (7 digits long) is available!</i> 4. As subject, please write „License request“. 5. As receiver, please take license@witeg.de 6. Send mail

After your license request, the information will be checked manually at witeg Labortechnik GmbH. As soon as possible you'll receive a mail with your license key.

PLEASE NOTE:

- The provided license key is connected to your machine and can't be consigned to other computers. If you want to move the software to another computer please install the software according to the instructions in this manual and repeat the licensing process.
- The licensing process is completely free of charge.
- Your provided information won't be used for other purposes than the licensing process.
- The license will be created manually by witeg Labortechnik GmbH within the regular working hours:

Central European time (CET) in Germany

Monday - Thursday: 7.00 AM till 4.00 PM (7.00 - 16.00 o'clock)

Friday: 7.00 AM till 1.00 PM (7.00 - 13.00 o'clock)

- If you don't receive a mail within 48 hours please contact us and tell us all needed information from the licensing process on phone or via mail at info@witeg.de.

License Information

To use the software you have to do the registration process.

Please fill out all relevant information under "Request Licence...".
If you're using Microsoft Outlook you can send us all information with the automatically created mail.
Otherwise please send a mail with all relevant information together with the Registration Key (7 digits) to license@witeg.de.

Your request will then be checked and the licence key send to the given mail address.

This process is free of charge. The sent data will only be used for the registration process.

Request Key:

License Key: - -

After receiving our license mail with your license key, please fill in the license key after starting Titrexsoft
The format of the license key must be:

XXXX - XXXX - XXXX

Please note:

You have to switch the fields manually. The cursor doesn't jump into the next field automatically.

After typing in the license key please click on "Activate license".

The software is now ready to use!

Please read the information provided on following pages to get in touch with the software. If you have any problems or suggestions for optimization feel free to contact us under license@witeg.de.

4. Interface description

Save-Button

Save current project

Open-Button

Open saved project

New-Button

Create new project

Close-Button

Close the software

Protocol-Button

Open printable protocol

Excel-Export-Button

Open Microsoft Excel with values from current project

Clipboard-Button

Copy values from current project to clipboard

DOS - TIT

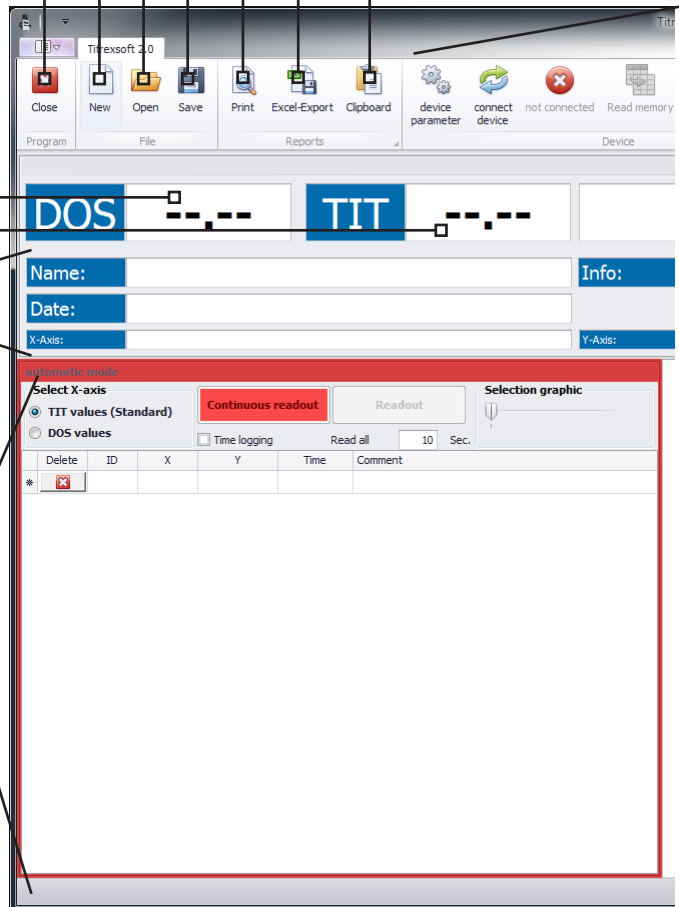
Live values from Titrex (if connected)

Information section

Editable information for current project.
The caption of x- and y-axis can be individually set

Project section

Will be explained in chapter **5. Project area**

Device section

This section will be explained in chapter **8. Device features**

Language-Buttons

Change language on the fly
German - English

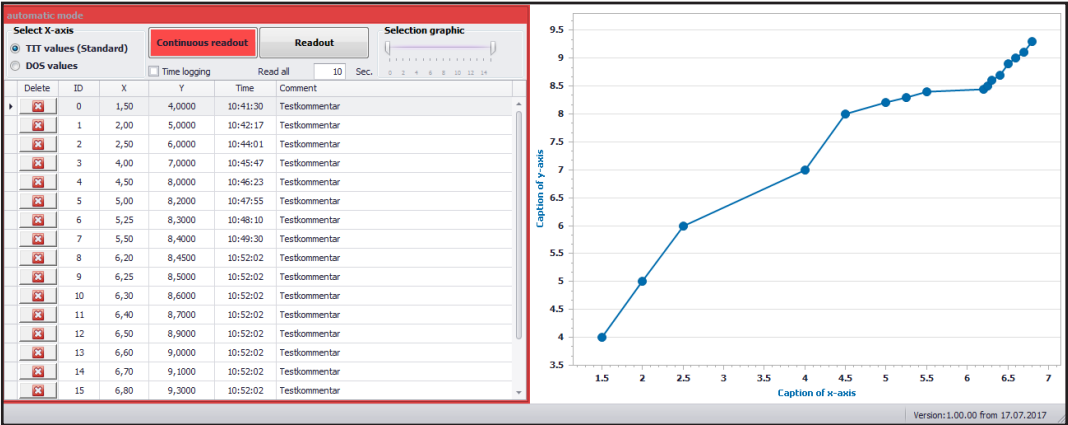
Graph section

Will be explained in chapter **5. Project area**

5. Project area

Please note: The shown values are just for explaining purposes and don't match any realistic results!

The project area is separated in two fields:


Value section	Graph section
<ul style="list-style-type: none">■ This section contains all values and information for each available measuring point■ There are two different modes: automatic and manual mode (see chapter 9. Titrexsoft modes) which can be used to add values	<ul style="list-style-type: none">■ This section is created automatically as soon as values will be added to the „value section“■ Via „Selection graphic“ (in the “value section”) the view can be changed

To get used to Titrexsoft there is an example project which can be opened via “Open” in the menu. The file is called “Beispiel.tit” and shows the same values as in the screenshot above.

6. Report functions & saving projects

Titrexsoft has three options to process all results shown in the “value section” and “graph section”:

Protocol function	Excel-Export function	Clipboard function
-------------------	-----------------------	--------------------

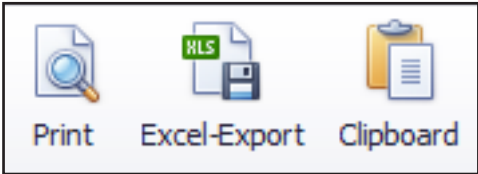Protocol function:

The built-in protocol function gives you a detailed report of all information from the “information section”, “value section” and the full graph of the “graph section” which can be printed directly or saved as PDF, RTF or XLSX file on the hard drive.

Excel-Export function:

The Excel-Export function exports all information from the “value section” directly into Microsoft Excel for further processing.

Please note: The graph from the “graph section” won’t be exported with the Excel-Export function!

Clipboard function:

The Clipboard function stores all information from the “value section” to the clipboard. The stored values are tab-separated and can be used in different programs for further processing.

Please note: The graph from the “graph section” won’t be exported!

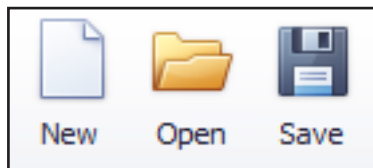
Saving projects:

Titrexsoft gives you the possibility to save your projects.

To save a project, click on the “Save” button in the menu and choose a location on any drive you want. Assign a clear project or file name.


Titrexsoft uses its own file extension *.TIT. This file extension can be directly linked to Titrexsoft and opened by a double-click*.

To open saved projects, double-click on the previously saved file in your file explorer (if linked via Microsoft Windows) or press on “Open” and choose the file within Titrexsoft.

** To link the file extension, double-click on an existing *.TIT file. A Microsoft Windows prompt will ask you which program to use to open the files in future. Please choose the option where a program can be selected manually. Titrexsoft is no standard Microsoft Windows software and must be linked manually by looking for the Titrexsoft2.EXE. The standard installation path is C:\Titrexsoft2 on your harddrive. Please make sure, that after you have selected the Titrexsoft2.EXE the option for “open these extensions always with the selected program” is checked to link it permanently.*

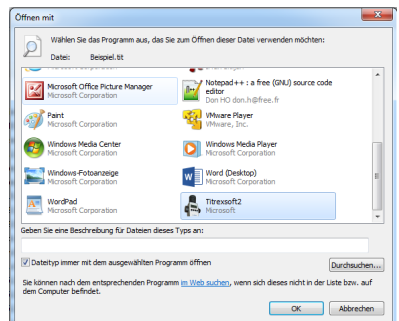More information about linking a file extension to a specific program can be found in the Internet. witeg Labortechnik GmbH isn't responsible for any loss of data or operating system crashes.

7. Device connection

To use Titrexsoft together with the digital burette Titrex, you need to follow these steps:

1. Connect the device with the supplied RS232 connection cable.
2. Connect the RS232 cable with your digital burette Titrex (*image #1*) and computer (an optional USB to RS232 adapter can be ordered separately - article number "5 497 920").
3. Turn on the digital burette Titrex by pressing "MEM" or "DOS".
4. In Titrexsoft choose from the menu "device parameter" (*image #2*) and pick the COM port from the digital burette Titrex (*image #3*); all other settings shouldn't be changed.

Please note: The drop down menu only contains available COM ports. If your COM port isn't shown you have to check the connection!

If you're using a RS232 to USB adapter please make sure, that Microsoft Windows has installed the drivers correctly. In general, the drivers will be downloaded automatically via Windows Updates if not available. If the driver installation fails please, contact your local IT department or responsible person for further assistance.

5. Click on "Save" to save settings (*image #3*).
6. Click on "Connect device" to establish the connection to your digital burette Titrex. If the connection was successful it will be shown in the menu with state "connected" (*image #4*).

image #1

image #2

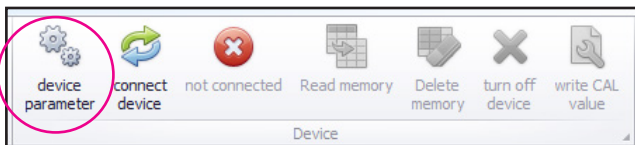
image #3


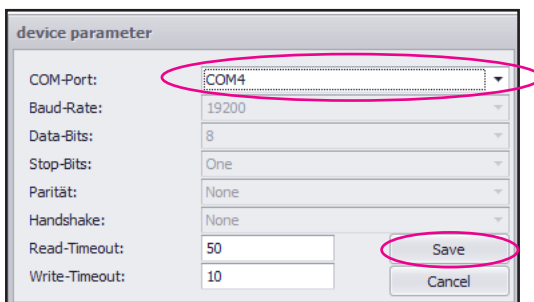
image #4

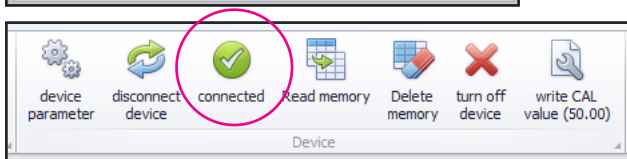To ensure that the device is connected properly with the computer and Titrexsoft, you can change the shown values from the “MEM” and “DOS” mode. The displayed values on the digital burette Titrex should match the shown values in Titrexsoft:

DOS

10.00

TIT

05.00

Titrexsoft

DOS mode	TIT mode
<ul style="list-style-type: none"> Indicated by a flashing dot between the first and second digit Showing 10.00ml in the display of the Titrex Showing 10.00ml in the software 	<ul style="list-style-type: none"> Indicated by a flashing dot between the third and fourth digit Showing 05.00ml in the display of the Titrex Showing 05.00ml in the software

Example values

Trouble shooting:


1	No COM port available after plugging in the digital burette Titrex <ul style="list-style-type: none"> Is the cable plugged in correctly? If a RS232 to USB adapter is used: Are the drivers correctly installed? Is the digital burette Titrex activated?
2	Connection not possible <ul style="list-style-type: none"> Repeat steps at chapter 7

If you have tested all possible actions to resolve a connection problem without success, please contact your IT department for further assistance!

8. Device features

After a successful connection of the digital burette Titrex with Titrexsoft, you gain access to different features:

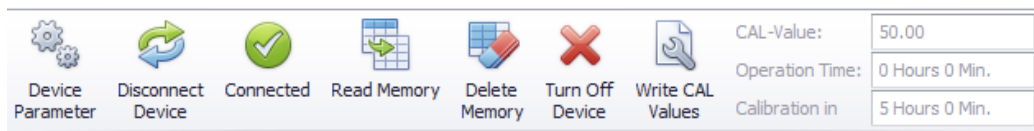- Read memory
- Delete memory
- Turn off device
- Write CAL values
- View working hours of digital burette Titrex since last reset
- View remaining working hours until next calibration

Read memory:

By selecting "Read memory", stored values will be directly exported from the digital burette Titrex into Microsoft Excel (will be opened automatically).

The storing process is explained in the Titrex operation manual provided with the scope of supply of the digital burette Titrex. If you don't have the manual please contact us.

Delete memory:

By selecting "Delete memory", all stored values in the digital burette Titrex will be deleted.

Please note: Deleted values can't be restored! You get prompted two times before the values are deleted permanently.

Turn off device:

By selecting "Turn off device" the digital burette Titrex will be turned off. This helps enhancing battery life-time.

By turning off the device, no data will be deleted from the internal memory.

Write CAL values:

By selecting "write CAL values", two options can be changed:

1. **CAL value:** changes the standard CAL value of 50.00.
Please note: The CAL value should only be changed in case of usage with liquids of varying density and viscosity or for a re-calibration. The meaning of the calibration value is explained in the Titrex operation manual provided with the scope of supply of the digital burette Titrex. If you don't have the manual please contact us. To change the CAL value for different liquids please use existing exchange values in relation to water.
2. **Calibration reminder time:** sets the operation time until the digital burette Titrex reminds for a re-calibration process. After first installation and connection of the digital burette Titrex with Titrexsoft, the re-calibration reminder is set to 5.00hrs.

Re-Calibration process

In order to make a re-calibration, the following equipment is needed:

- Analytical balance with a resolution of 1mg, repeatability and linearity of 2mg and standard uncertainty of 3µl
- Temperature probe or thermometer with a standard uncertainty of maximum 0,2°C
- Barometer
- Vessel made of Borosilicate glass 3.3 or AR-glass
- Knowledge about the correct usage of digital burette Titrex and preparation for correct measurement

To support the recalibration process, you can find a spreadsheet in the installation path of your Titrex-software (Titrex-calibration-spreadsheet.xlsx).

Steps to Re-Calibrate the digital burette Titrex:

1. Prepare the digital burette Titrex according to the operation manual.
2. Make sure the vessel used for weighing is prepared for a correct result.
3. If you are using the Titrex-calibration-spreadsheet.xlsx fill out the fields for "Volume" (for Titrex calibration it is 50), "Room temperature", "Water temperature", "Air pressure" and choose "Type of glass" you are going to use for the weighing.
4. Switch the Titrex mode to "DOS"-mode and fill up until "50.00" is shown on the display.
5. Switch the Titrex mode to "TIT"-mode and dispense water until "50.00" is shown on the display into the previously prepared vessel.
6. Put the vessel onto the analytical balance and note the shown value in the Titrex-calibration-spreadsheet.xlsx or in your own existing water surcharge counting tool.
7. Tare the analytical balance and repeat the process from point 4 until five values have been weighed to create an average value.
8. If you have used the Titrex-calibration-spreadsheet.xlsx, you should now have the calculated CAL-value which can be stored into the Titrex memory (see "Write Cal values").
If you have used your own tool, please make sure, that you consider the surcharge for the liquid before changing the CAL value of the digital burette Titrex!

Please note: witeg Labortechnik GmbH isn't responsible for any changed values by the user! If you are uncertain about the calibration and/or re-calibration process, please contact us!

The following information is shown in the menu bar about the CAL values:

The current CAL-Value. **Standard value after delivery is 50.00.**

The operation time for the currently attached digital burette Titrex. **This value is set to 0hrs 0min after first installation and connection of the digital burette Titrex with the software!**

Countdown for next calibration time. **This value is set to 5hrs 0min after first installation and connection of the digital burette Titrex with the software!**

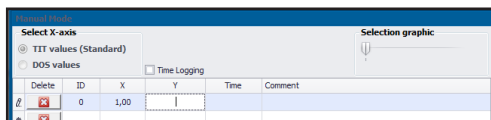
9. Working modes

With Titrexsoft, you have two working modes which support your whole titration application:


Manual mode

Automatic mode (a connected digital burette Titrex is needed)

Manual mode is indicated by a **blue** frame

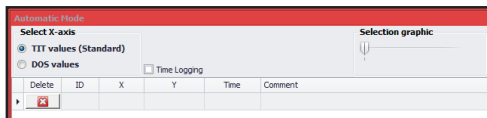Automatic mode is indicated by a **red** frame

a. Manual mode

The purpose of the manual mode is to type in existing hand-written titration documentations into a digital format for further usage. This could be printing out a pre-built protocol directly out of Titrexsoft or just to save your data on a hard-drive for digital storage (more information in chapter 6. Report functions & saving projects).

It is also possible to record data while doing the titration application without the need for a hand-written documentation.

After starting Titrexsoft, you automatically enter the **automatic mode** (indicated by a red framed value section). As soon as you enter the first value for the x- or y-axis you enter the **manual mode** (indicated by a blue framed value section).

Workflow:

- Type in the values for “X” and “Y” to automatically create the graph in the “graph section”.

b. Automatic mode

The purpose of the automatic mode is to transfer stored values from the digital burette Titrex (“MEM”-Button) directly into the value section of Titrexsoft. Therefore, a connected digital burette Titrex is needed (see chapter 6. **Device connection**).

Preparation:

- Connect your digital burette Titrex with Titrexsoft (explained in chapter 7).
- Click on “New” to create a new project and to enter the automatic mode.
- You'll be asked to save the current project.
- After saving/aborting the current project, Titrex memory will be entered by Titrexsoft. If there are any existing values in the memory, you'll get asked if they should be deleted or not.
 - **Option 1:** Don't delete the stored values if you want to work with a previously made titration application within Titrexsoft
 - **Option 2:** Delete the stored values to have a clear memory for a new titration application which should be directly transferred to Titrexsoft by pressing the “MEM”-Button

Workflow Option 1 - working with existing values in the Titrex memory:

- Click on “Readout” to enter the memory and transfer all stored values into the value section for further editing.
- After the values have been transferred you can enter the corresponding y-axis values. The graph will be created automatically.
- If you activate “Time logging”, the current computer time will be added to all rows
- If you change any x-axis value, the **automatic mode** will be left and you get access to the **manual mode** with all readout values.

Workflow Option 2 - working instantly with stored values in Titrexsoft:

- Click on “Continuous Readout” to gain permanent access to the memory of the digital burette Titrex.
- If you didn't delete the memory while creating a new project, all stored values will be transferred into the value section of Titrexsoft. You have now the possibility to add new values by pressing the “MEM”-Button on the digital burette Titrex.
- If you deleted the memory, you can now add new values by pressing the “MEM”-Button on the digital burette Titrex.
- If you activate “Time logging”, the current computer time while storing the value will be added to the row.
- After the values have been transferred, you can enter the corresponding y-axis values. You can also set a waiting time for Titrexsoft when the next reading cycle starts. This gives you the possibility to add y-axis values directly into Titrexsoft during the titration process.
Please note: The cursor will jump into the first row when a reading cycle starts! You have to determine the speed according to your needs and working speed if you want to add corresponding y-axis values during a titration process.

Features and best practice for both modes:

- Navigate through the cells with the arrow key on your keyboard or click in the cell with your cursor.
- If “Time logging” is activated, the computer time will be added automatically while typing in the values.
- If needed fill out the “Comment” column for additional information in a specific row.
- Wrong rows can be deleted with the “X” in the first column (Please note: a deleted row can't be retrieved afterwards and must be added manually).
- Sort the values according to different columns by pressing on the top line.
- The graph will always be created increasingly according to the values from the x-axis.
- Create a more detailed view on interesting sections of the graph by changing the first and last shown value with the “Selection graphic” slider.
- If you're using Titrexsoft in “English” language mode, please use a **decimal point** for typed-in values.

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