WAC-R

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Contents

1 Introduction	18
2 Safety Instructions	18
2.1 Description of symbols.....	18
2.2 Instructions	18
3 Packing content	20
4 Product description	20
4.1 Key features.....	20
4.2 General survey	21
4.3 Operating surface.....	22
4.4 Term description	22
5 Positioning the device and commissioning.....	22
5.1 Positioning the device	22
5.2 Commissioning the device	23
5.3 Emergency Lid unlock method	23
6 Operation	24
6.1 Menu structure	24
6.2 Set Operation Timer	24
6.3 Set Delay Timer	24
6.4 Set temperature.....	25
6.5 Temperature Offset	25
6.6 Set Printer	25
6.6.1 Set Date & Time	26
6.6.2 Set Name.....	26
6.7 Locking mode.....	26
6.8 Set Operation Cycle Mode	26
6.9 Storage function.....	27
7 Maintenance and cleaning.....	27
7.1 Change Printer Paper Roll	27
7.2 Battery Replacement Method (3V batteries for Date and Time Output).....	28
8 Troubleshooting Guide	29
9 Technical data.....	30
10 Certificate.....	31

1 Introduction

Thank you for purchasing our **WAC-R Steam Sterilizer**. You did a good choice buying a product from witeg Labortechnik.

This operation manual describes performance, usage and cautions for use of the purchased product. Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you specially have to follow the safety instructions of the second chapter imperatively.

Please keep this operating manual in a save place for future reference and hand it out if you consign the device to a third person.

Modifications of the capsule or amendments for the improvement of quality and efficiency can be made by **witeg** without prior notice.

2 Safety Instructions

2.1 Description of symbols

**WARNING**

Information for avoiding injuries or fatal accidents.

**ATTENTION**

Safety instruction to avoid a damage of the product.

**ATTENTION**

Information about technical parameters.

**Environment**

Information about the operating environment.

**WARNING of explosions**

Safety instruction to avoid explosions.

**ATTENTION to high temperatures**

Instructions to avoid injuries caused by high temperatures.

2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Connect the device only to the suitable power supply. Please have a look to the identification plate.

- Use the device only in electric circuits with ground fault circuit interrupter.
- Check the power supply cord before you stick it into the electrical outlet.
- Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.
- Do not touch the device with wet hands during the usage because this might cause an electrical shock.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.
- Do never try to open the lid during the operation.

- Do not use this device in area with potentially explosive atmosphere.

- Repairs and amendments should only be made by **witeg** authorized people. Contraventions exclude the warranty.

- Do not put heat-sensitive items at the device's housing.
- The lid and other parts could become very hot during the stirring process.
- Wear protection gloves.
- Warn all attendees if you leave the device unattended.
- Let the pressure fall to "0" and the temperature under 80°C before you open the lid.
- Pay attention when you open the lid, hot condensation water may drop down and hot steam could come out.
- Do not clean the scorching device.
- Ensure that the resins steam backwards of the device can not cause injuries or material damages.

- Use this device only if you read the whole safety instructions.
- This device should only be used by persons who are familiarized with safety instructions made for laboratories.
- The lid can open only if the pressure in the vessel is fallen to "0" and temperature is under 80 °C.
- Pay attention that no liquid reaches the controller or inward of the capsule.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective work clothing during the usage.
- Switch the device off before you disconnect it from the power grid.
- All circuits and electro technical components are **witeg** owned technology. In fact of this it is necessary that only **witeg** authorized personal repair the device.

- Pull out the power cable if you didn't want to use the device for a longer time.
- Do not cover the ventilating openings.
- Use only distilled water.
- Only sterilize autoclave able items otherwise they could lose their function and safety features.

- It is important that the device stands up on a horizontal, stable and firm subsurface to avoid vibrations.
- Make sure that the device cannot slip away, use therefore the caster's brakes.
- To avoid heat congestion or fires make sure that there is enough space for air circulations.
- Avoid direct solar radiation.
- Do not use the device near highly combustible materials.
- Do not set the device outdoors.
- It is superiorly to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- To set the device in wet or dusty places could cause overheating, short circuits or fires.

3 Packing content

Quantum	DH.WACR2047	DH.WACR2060	DH.WACR7080
1	WAC-R47	WAC-R60	WAC-R80
2	Wire basked		
1	Bottom strainer		
1	Printer Paper Roll		
1	Operating manual		

4 Product description


4.1 Key features

The **WAC-R Steam Sterilizer** is designed for chemical, biological, medical and pharmaceutical applications.


Key features are as follows:

- Long life by simply and solid construction.
- CE\MDD
- For Steam sterilisation up to 132°C, 2 bar.
- Simplest operating by patented Jog-Dial control system.
- The sterilization results printed out by built-in panel print are analyzed and reviewed, and the confirmation of the cycle execution record is provided.
- Solid/Liquid Mode can be chosen considering a Sample Phase.
- Fast commissioning and easy loading with top-loader.
- Highest temperature accuracy by digital Fuzzy controller.
- Temperature and delay time showed on illuminated Display.
- Storage functions for time and temperature values.
- Turn-off delay from one minute up to over 99 hours in one minute steps adjustable.
- Signal sound after end of turn-off delay.
- Casters for excellent mobility.

4.2 General survey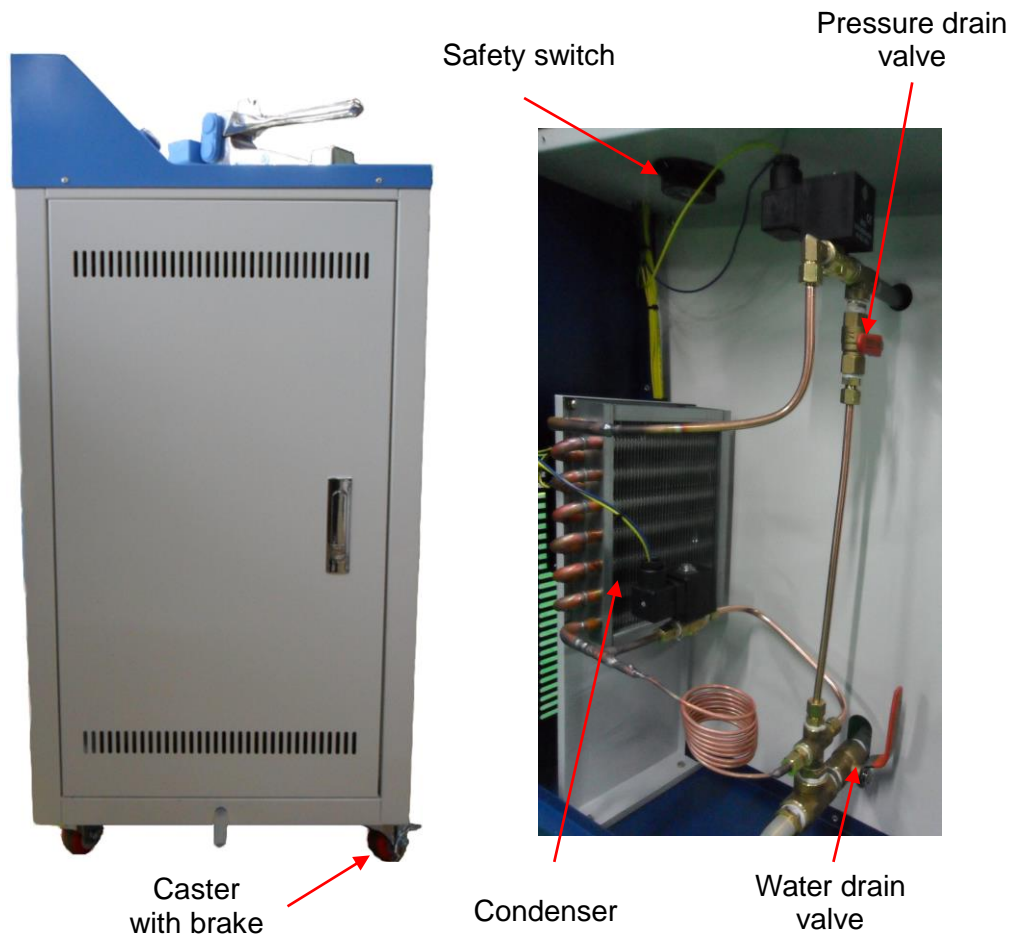

4.3 Operating surface

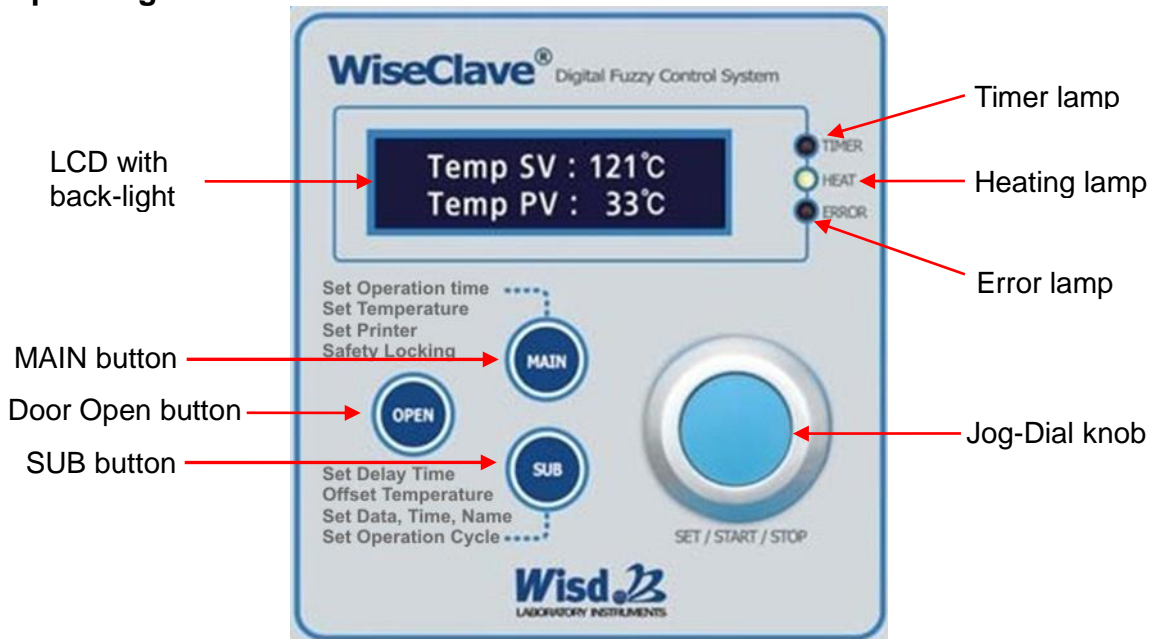4.4 Term description

SV: set value (upper line, in key lock menu no SV)

PV: present value, measured by the device (lower line, in key lock menu also upper line)

Timer lamp

- Is blinking in menu items for timer modulation
- Shines continuous in the key lock menu item when timer activated.

Heating lamp

- Shines if the heater is activated.

Error lamp

- Shines when an Error message is present.

Jog-Dial knob

- Changes values by turning, stores, confirm, stops and starts by pushing.

MAIN button

- For changing between the different menu items and the key combination.

SUB button

- To reach the sub menu Items.

OPEN button

- For unlocking the Opening mechanism

Safety switch

- The safety switch turns the heating off if your set maximal temperature is overshoot.
- Set the maximal temperature rough 20 – 30 °C over the desired maximal temperature, otherwise the device maybe not reaches your set value.

5 Positioning the device and commissioning

5.1 Positioning the device

1. Follow especially the safety instructions marked with following symbol:
2. Remove the package and perhaps soiling carefully.
3. If there are any unexpected transport damages, please inform your forwarder and your customer service.

4. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatized.
5. Connect the water outlet with a caisson or catch tank. Please recognize that hot water is flowing out.
6. Fix the brakes of the casters.

5.2 Commissioning the device


1. Before start, proceed as described in **5.3** before
2. Fill distilled water in and make sure the water is not topping the bottom strainer.
3. Make sure that you followed all safety instructions.
4. Connect the device with the power grid.
5. Load the vessel.
6. Pull the lid over the vessel and make sure it is exactly placed over the vessels opening. Otherwise may hot steam emit.
7. Close the lid with the hand lever.
8. Switch the device on and start the process (look at job description).

5.3 Emergency Lid unlock method

The autoclave has electronic safety lock system.

If power failure or something wrong in power supplier of units or commissioning the device, customer couldn't open the lid by push the "OPEN" button.


Please proceed as follows: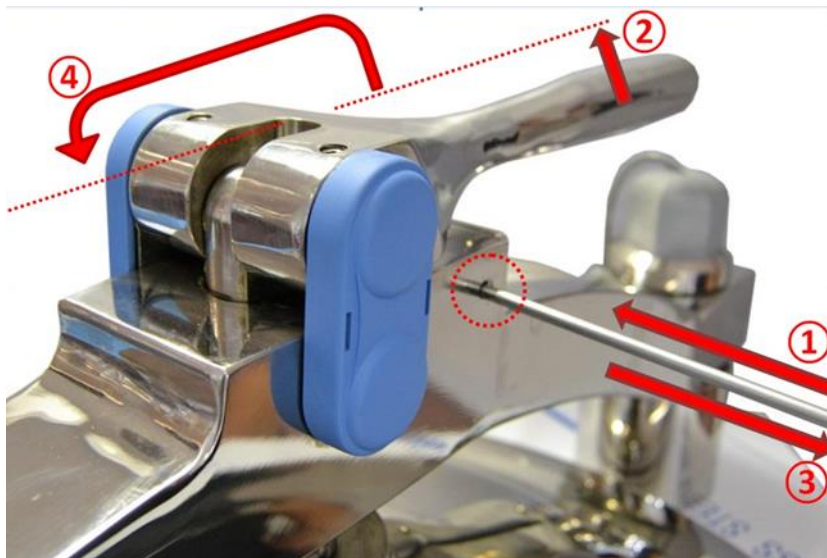

1. Perform an appropriate object (eg screwdriver) into the provided Opening, until you encounter resistance.
2. Pull lever in the opening direction.
3. Remove the object from the aperture.
4. Place the lever completely.

6 Operation

6.1 Menu structure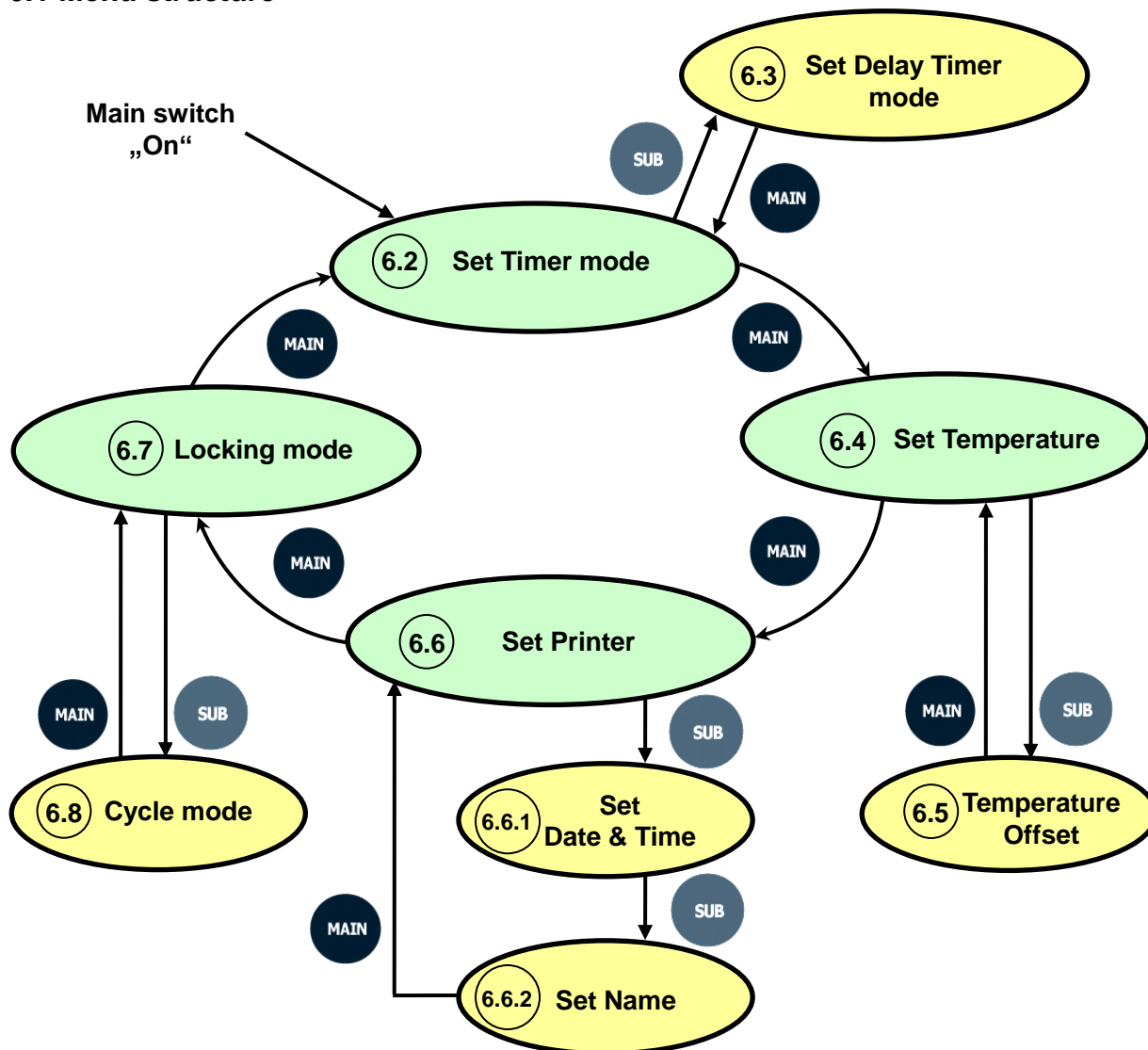

6.2 Set Operation Timer

The Operation Timer serves as automatically stop function and opens the steam outlet after the time is elapsed.

- Use MAIN button to reach next menu without setting a delay time.
- The delay time is adjustable from one minute till over 99 hours (00:01 - 99:59) in one minute steps.
- Turn the Jog-Dial knob to set the delay time. Push the Jog-Dial if you reached the desired value to confirm it. With the confirmation the timeframe starts elapsing.
- If you want to change the value after confirmation, you have to push the Jog-Dial once again. The PV initiates blinking and you can modify the value again.
- To confirm the value and start the delay time push the Jog-Dial yet again.
- **With “00:00” value of Timer SV, the machine cannot start operation.**

6.3 Set Delay Timer

The Delay Timer serves as automatically start function. It turns on the heating after the timeframe you set is elapsed.

To reach this sub menu item you have to use the SUB button in the menu “set turn-off delay”

- The delay time is adjustable from one minute till over 99 hours in one minute steps.
- Turn the Jog-Dial knob to set your delay time. Push the Jog-Dial if you reached the desired value to confirm it. With the confirmation the timeframe starts elapsing.
- If you want to change the value after confirmation, you have to push the Jog-Dial once again. The PV initiates blinking and you can modify the value again.
- To confirm the value and start the delay time push the Jog-Dial yet again.

6.4 Set temperature

- When you push the MAIN button in the menu item “set turn-off delay” you reach the menu item “set temperature”. You can recognize this menu at “°C” in SV and PV line.
- The temperature is adjustable from 25°C till 132°C in 0,1°C-steps.
- Push the MAIN button once again if you want to work without the heating function. Then you reach the next menu without collimated temperature value.
- Choose your desired temperature by turning the Jog-Dial and confirm the set value by pushing the Jog-Dial.
- For changing the confirmed value again, you first have to push the Jog-Dial knob again. “HEATING” starts blinking at the LCD and the temperature is adjustable again.
- When the heating lamp is shining, the heater is activated.

6.5 Temperature Offset

If the temperature measured by you differs from the temperature the device is showing (PV) you have the possibility to compensate the divergence by the Offset. So you can balance differences of about ± 10 °C.

The Offset is calculated as follows:

Temperature measured by you – showed temperature (PV) = Offset

- To reach this menu you have to use the SUB button in the menu item “set temperature”.
- Set the calculated value by turning the Jog-Dial. The Offset will be added to the PV immediately.
- Save the offset by holding the Jog-Dial pushed until the values in PV and SV line starts blinking.
- Push the MAIN button to leave the justification menu.

6.6 Set Printer

- When you push the MAIN button in the menu item “set temperature” you reach the menu item “set printer”.
- Turn ON/OFF the output functions of the printer that can record the sterilization operation process.
- If pressing the PUSH of Jog Dial Knob, the Display of ON/OFF of the printer function is changed. "Set to Print" and "Set not to Print" of the Display are alternately changed every time when the button is pressed.
- If pressing and holding the PUSH button, the set screen is stored, fixing while blinking the words of the Display.
- Printing is automatically carried out at 15 minutes (end of exhaust stage) after completing 1 Cycle (80°C basis) of normal sterilization.
- If pressing the Door Open button at the exhaust stage, the sterilization operation process is printed every time you press the button. However, if the user prints using the button, the automatic printing stage for 15 minutes is omitted.
- If an Error is occurred during operating, print the Data till immediately before the Error is

occurred, and then the Error message is printed. However, if an Error is occurred while proceeding till the sterilization stage, the Error details is printed out, and if an Error is occurred in the initial process, it is not printed. ex) Error_#1 : 2014.03.07_12:34:56

6.6.1 Set Date & Time

- To reach this menu you have to use the SUB button in the menu item “set printer”.
- Set up the current date and time to record on the printout paper of the printer.
- If indicating the standby mode while the cursor is blinking, the user can change the setting value by turning the Jog-Dial Knob to the right/left.
- If pressing the PUSH button, the Displayed value is stored, and the cursor will be moved to the next position.
- If pressing and holding the PUSH button after setting both DATE and TIME, the values are stored, fixing while blinking the screen of Display.

6.6.2 Set Name

- To reach this menu you have to use the SUB button in the menu item “Set Date & Time”.
- The title of experiment or the sterilization information to record on the printout paper of the printer can be entered.
- If indicating the standby mode while the cursor is blinking, the user can change the setting character and numbers by turning the Jog-Dial Knob to the right/left.
- If pressing the PUSH button, the Displayed value is stored, and the cursor will be moved to the next position.
- If pressing and holding the PUSH button after setting the Name, the values are stored, fixing while blinking the screen of Display.
- If pressing the MAIN button after completing all the settings, the screen is converted to the Set Printer mode.
- Press the MAIN button to leave the “set name” menu. Than you are in the menu item “set printer”.

6.7 Locking mode

- To avoid accidental changing of the set values, push the MAIN button in the menu item “set printer”.
- In this menu item can you see the actually temperature in the SV-line and the remaining time of turn-off delay in the PV-line. The function of the Jog-Dial knob is deactivated.
- Press the MAIN button once more to leave the key lock. Than you are in the menu item “set turn-off delay”.

6.8 Set Operation Cycle Mode

WAC-R Steam Sterilizer offer two different operation cycles named SOLID cycle mode and LIQUID cycle mode. As the name implies, LIQUID cycle is for autoclaving liquids while the SOLID cycle is for autoclaving solid or dry materials. The following table shows you the difference of the two cycles and proper use:

Type of Cycle	Appropriate Material	Description
SOLID or “fast exhaust”	Dry goods, glassware, etc.	This cycle charges the chamber with steam and holds it at a set pressure and temperature for a set period of time. At the end of the cycle, a valve opens and the chamber rapidly returns to atmospheric pressure.
LIQUID of “slow exhaust”	Liquids	This cycle prevents sterilized liquids from boiling. Steam is exhausted slowly at the end of the cycle, allowing the liquids (which will be super-heated) to cool.

- To reach this menu you have to use the SUB button in the menu item “locking mode”.
- Push the Jog-Dial knob to change the operation mode.
- Push the MAIN button to leave the operation cycle mode.

6.9 Storage function


The **WAC-R Steam Sterilizer** includes a storage function for temperature and time values. That is practical if you want to work with the same parameters often. Then you did not need to set them in every usage.

- For storing values you only need to hold the Jog-Dial knob pushed for two seconds. Successful storing is displayed by blinking of the SV for three times.
- Any stored parameters are erasable. Turn the value with the Jug-Dial till “0” and then hold the knob pressed for two seconds.
- If you want to overwrite a value it is not necessary to erase the old one. Set the desired value and use the Jog-Dial knob for storing.

7 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device.
- Pull the power plug out of the electrical outlet.
- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- Only use original spare parts.
- Please control every three month the condenser. If it is to much wasted with dust, please clean it with compressed air or a strong vacuum cleaner. Please handle the lamellas with care.
- The device’s pressure vessel has to been checked to ensure your safety regularly.

7.1 Change Printer Paper Roll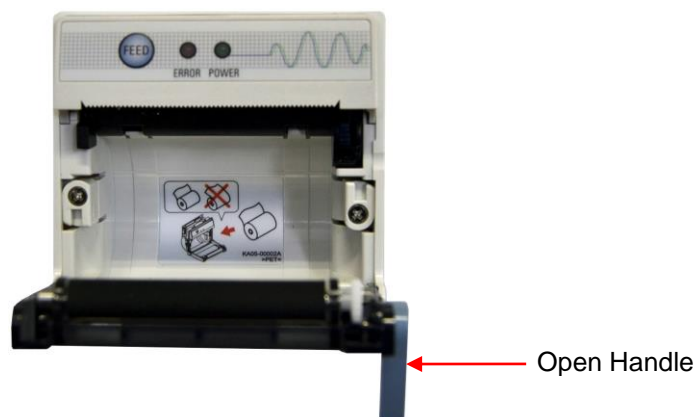

- Please check the amount of the paper before the sterilization operation, and when the printout paper is lack, replace it to new printout paper.
- Open the place where installs the paper of printer by lifting the handle located in the right of the printer, and remove the spent core of the paper.
- As shown in above Figure, insert the paper with facing up in the state that the starting part of new paper is gently come out of the printer, and close the door of printer and



complete the installation of the paper.

- If the end of the paper is hanging to a roller, or the margin needs back and forth of the outputs, the user can take the paper out a little bit by pressing the “FEED” button.

7.2 Battery Replacement Method (3V batteries for Date and Time Output)

- When the power is turned ON, need to replace the battery when the Date and Time stored disappear and initializes. (initialization state: DATE : 2000/00/00 , TIME : 00:00:00)
- **Shut off the power** of the product and open the Door on the Left Side. Remove the distribution board panel as shown in the figure.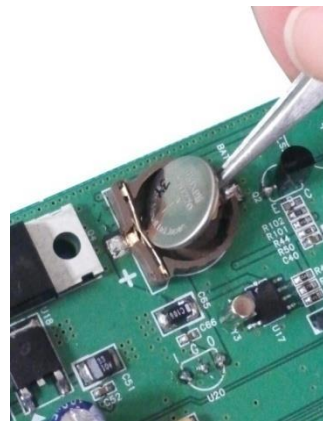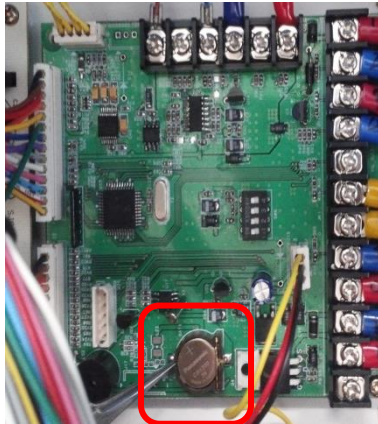

- After panel removal Check the Main Controller Board located in the power distribution board.
- After checking the with 3V battery of Main Controller Board, put a tweezers to gap and remove the battery.
- Replace a new battery and turn ON the power of product and then reset the Date and Time
- The battery replacement period is usually two years, please be sure to use the same specification of shipped factory battery.
- May be shortened replacement cycles depending on the initial condition and environment of the battery.

8 Troubleshooting Guide

Situation	Confirmations and Solutions
Power Failure	<ul style="list-style-type: none"> ✓ Check the supplied voltage is present. ✓ Check the power connection and the power cord for proper connection and integrity. ✓ Check if the MAIN Power Switch is "ON". ✓ Fuse is blown.
Unit Not Heating	<ul style="list-style-type: none"> ✓ Check if the MAIN Power Switch is "ON". ✓ Check if the set temperature (SV) confirmed with the Jog-Dial knob (by turn the Jog-Dial knob cannot change the set temperature – see point 6.4) ✓ Check the Safety switch is set higher (min. 10°C-20°C) than the set temperature (SV).
Heating, but unit not reaching temperature or temperature is fluctuated	<ul style="list-style-type: none"> ✓ Check the Safety switch is set higher (min. 10°C-20°C) than the set temperature (SV). ✓ Problem in the sensor or main controller.
ERROR 1	✓ Temperature Sensor Error.
ERROR 3	✓ Low level error. Please fill water in.

If other problems arise, or one of the above not solves problems, please contact your official agent or the manufacturer.

Reparations and amendments should only made by [witeg](http://www.witeg.de) authorized people. Contraventions exclude the warranty.

9 Technical data

		WAC-47	WAC-60	WAC-80
Dimension (B x T x H)	Inner (mm)	Ø 300 x h 670	Ø 350 x h 650	Ø 400 x h 650
	Capsule (mm)	705 x 475 x 1070	705 x 475 x 1070	815 x 585 x 1170
Capacity (litre)		47	60	80
Heating power (kW)		2.0	3.0	4.5
Temp.	Range	Up to 132 °C		
	Accuracy	± 0,5°C at + 121°C		
	Sensor	PT100		
Vessel Pressure		max. 2 bar		
Controller		Digital Fuzzy-controller with Jog-Dial knob (Turn + Push)		
Display		Digital LCD with Back-Light Function		
Timer		99 h 59 min (for Turn-off and Turn-on delay)		
Recorder				
Material	Vessel	Stainless Steel		
	Capsule	Powder coated Steel, polished stainless steel		
	Insulation	Polystyrol (20 mm)		
Safety devices		<ul style="list-style-type: none"> - over temp. and over current protector - sensor control - over pressure valve 		
Others		<ul style="list-style-type: none"> - Storage function (Temp., Timer) - Key lock (Jog-Dial deactivated) - Safety electronic door closing system <ul style="list-style-type: none"> - Easy to justify - Signal sound 		
Packing Size (mm) & Gross Weight		800 x 600 x 1.250 116 kg	800 x 600 x 1.250 120 kg	930 x 700 x 1.350 173 kg
Power supply		AC 230 V, 1N~, 50/60 Hz		AC 400 V, 3N~, 50/60 Hz

10 Certificate

Certificate

for

The quality and all features were checked by the manufacturer before the shipment.

We grant from date of purchase

two years guaranty.

This certificate excludes damages by natural disasters or incorrect usages by the costumer.

Please look on your account and complete following table:

Article	WAC-R Steam Sterilizer
Typ	
Serialno.	
Date	

witeg Labortechnik GmbH

Am Bildacker 16
D-97877 Wertheim
TEL: +49-9342-9301-0
FAX: +49-9342-9301-77
Email: info@witeg.de
www.witeg.de