WHM

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www.witeg.de

Content

1 Introduction	8
2 Safety Instructions	8
2.1 Description of symbols	8
2.2 Instructions	8
3 Product description	10
3.1 Key features	10
3.2 Restrictions	10
4 Package content.....	10
5 Commissioning	10
6 Maintenance and cleaning.....	10
7 Technical data	11
9 Certificate.....	12

1 Introduction

Thank you for purchasing our **WHM Heating Tape without Controller**. You did a good choice buying a **Wisd** product.

This operation manual describes performance, usage and cautions for use of the purchased product. Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you specially have to follow the safety instructions of the second chapter imperatively.

Please keep this operating manual in a save place for future reference and hand it out if you consign the device to a third person.

Modifications of the capsule or amendments for the improvement of quality and efficiency can be made by **witeg** without prior notice.

2 Safety Instructions

2.1 Description of symbols

**WARNING**

Information for avoiding injuries or fatal accidents.

**ATTENTION**

Safety instruction to avoid a damage of the product.

**ATTENTION**

Information about technical parameters.

**Environment**

Information about the operating environment.

**WARNING of explosions**

Safety instruction to avoid explosions.

**ATTENTION to high temperatures**

Instructions to avoid injuries caused by high temperatures.

2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Only use our controller **WHM-C10D** (Art.-No.: DH.WHM12500) or **WHM-C10A** (Art.-No.: DH.WHM13500) to heat up the Heating Tapes.
- Connect the device only to the suitable power supply. Please have a look to the identification plate.
- Due to its functions the Heating Tapes is not protected against incoming water. Only a professional handling can avoid hazardous accidents.
- Only persons with sufficient safety knowledge are allowed to use this device.
- Use the device only in electric circuits with ground fault circuit interrupter.
- Do not touch the device if any liquid is flowed into the device (e.g. broken glass vessel) and firstly switch the power supply off. After that please pull out the power plug.
- Check the power supply cord before you stick it into the electrical outlet.
- Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.
- Do not touch the device with wet hands during the usage because this might cause an electrical shock.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.

- Do not put heat-sensitive items at the device.
- Only use adequate sized vessels and avoid an over boil of the liquids.
- Do not clean the scorching device.

- Use this device only if you read the whole safety instructions.
- This device should only use by persons who are familiarized with safety instructions made for laboratories.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective work clothing during the usage.
- Pull out the power cable if you didn't want to use the device for a longer time.
- Do not heat up the heating tape without any vessel to avoid an overheating.
- Do not heat the up the device more than it is necessary - for lower energy consumption and your own safety.
- Make sure that the device cannot slip away.
- To avoid heat congestions or fires make sure that there is enough space for air circulations.
- Avoid direct solar radiation.
- Do not use the device near highly combustible materials.
- Do not set the device outdoors.

- It is superiorly to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- To set the device in wet or dusty places could cause overheating, short circuits or fires.

- Do not use this device in areas with potentially explosive atmosphere.

- Reparations and amendments should only made by **witeg** authorized people. Contraventions exclude the warranty.

3 Product description

3.1 Key features

The **Wisdom WHM Heating Tape without Controller** is designed for chemical, biological, medical and pharmaceutical heating and extraction applications up to 250 °C. The device was developed for simplest usage and a very long life. This Heating Tapes has to be used only with our **WHM-C10D** or **WHM-C10A** external controller.

3.2 Restrictions

- Please only use glass vessels that are made for heating applications. A minimum heat resistance like borosilicate glass 3.3 has is needed.
- Usage is only allowed with controller **WHM-C10D** controller (Art.-No.: DH.WHM12500) or **WHM-C10A** controller (Art.-No.: DH.WHM13500).

4 Package content

Quantity	Description
1	WHM Heating Tape without Controller
1	Operating manual

5 Commissioning

1. Follow the safety instructions.
2. Remove the package carefully.
3. If there are any unexpected transport damages, please inform your supplier.
4. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatized.
5. Connect the heating tapes power cord to the controllers power socked for heating tapes.
6. Switch the device on.
7. Please follow the job description of **WHM-C10 #** controller.

6 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device.
- Pull the power plug out of the electrical outlet.

- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- Only use original spare parts and accessories.
- Make sure that no liquid gets into the device.
- Do not clean the hot device.

7 Technical data

Heating Tapes, 250°C, 230V		
Article number	Width x Length	Watts
DH.WHM12311	13 mm x 0,6 m	200
DH.WHM12312	13 mm x 1,0 m	200
DH.WHM12313	13 mm x 1,5 m	250
DH.WHM12314	13 mm x 2,0 m	300
DH.WHM12315	13 mm x 2,5 m	350
DH.WHM12316	13 mm x 3,0 m	400
DH.WHM12321	25 mm x 0,6 m	200
DH.WHM12322	25 mm x 1,0 m	200
DH.WHM12323	25 mm x 1,5 m	250
DH.WHM12324	25 mm x 2,0 m	300
DH.WHM12325	25 mm x 2,5 m	350
DH.WHM12326	25 mm x 3,0 m	400
DH.WHM12327	25 mm x 4,0 m	550
DH.WHM12328	25 mm x 6,0 m	700

9 Certificate

Certificate

for

The quality and all features were checked by the manufacturer before the shipment.

We grant from date of purchase
two years guaranty.

This certificate excludes damages by natural disasters or incorrect usages by the customer.

Please look on your account and complete following table

Article	WHM Heating Tape without Controller
Type	
Serial no.	
Date	

witeg Labortechnik GmbH

Am Bildacker 16

D-97877 Wertheim

TEL: +49-9342-9301-0

FAX: +49-9342-9301-77

Email: info@witeg.de

www.witeg.de