## **WHM-C10A**

# **Laborregler für Heizmäntel**

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# **Controller for Heating**

# **Mantles**

**Operating Manual**

**witeg Labortechnik GmbH**

(English manual at page 6)

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# Contents

|  |           |
|--|-----------|
| <b>1 Introduction .....</b>                            | <b>7</b>  |
| <b>2 Safety Instructions .....</b>                     | <b>7</b>  |
| <b>2.1 Description of Symbols .....</b>                | <b>7</b>  |
| <b>2.2 Instructions .....</b>                          | <b>7</b>  |
| <b>3 Product description .....</b>                     | <b>8</b>  |
| <b>3.1 Key features .....</b>                          | <b>8</b>  |
| <b>3.2 General Survey.....</b>                         | <b>9</b>  |
| <b>3.3 Term description .....</b>                      | <b>9</b>  |
| <b>4 Packing contents.....</b>                         | <b>9</b>  |
| <b>5 Positioning the device and commissioning.....</b> | <b>9</b>  |
| <b>7 Maintenance and cleaning.....</b>                 | <b>10</b> |
| <b>8 Technical data.....</b>                           | <b>10</b> |
| <b>9 Certificate.....</b>                              | <b>11</b> |

# 1 Introduction

Thank you for purchasing our **WHM-C10A Controller for Heating Mantles**. You did a good choice buying a **Wisd** product.

Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you have to follow the safety instructions of the second chapter.

Please keep this operating manual in a save place for future reference and hand it out if you consign the device to a third person.

**Witeg** reserves to modify the capsule or make amendments for the improvement of quality and efficiency without prior notice.

## 2 Safety Instructions

### 2.1 Description of Symbols

**WARNING**

Information for avoiding injuries or fatal accidents.

**ATTENTION**

Safety instruction to avoid a damage of the product.

**ATTENTION**

Information about technical parameters.

**Environment**

Information about the operating environment.

**WARNING of explosions**

Safety instruction to avoid explosions.

### 2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Connect the device only to the suitable power supply. Please have a look to the identification plate.
- Use the device only in electric circuits with ground fault circuit interrupter (RCD).
- Check the power supply cord before you stick it into the electrical outlet.
- Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.

- Do not touch the device with wet hands during the usage because this might cause an electrical shock.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.

- Reparations and amendments should only be made by **witeg** authorized people. Contraventions exclude the warranty and may cause fatal injuries.

- Use this device only if you read the whole safety instructions.
- This device should only be used by persons who are familiarized with safety instructions made for laboratories.
- Pay attention that no liquid reaches the controller or inward of the capsule.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective work clothing during the usage.
- All circuits and electrotechnical components are **witeg** owned technology. In fact of this it is necessary that only **witeg** authorized personnel repair the device.
- Pull out the power cable if you didn't want to use the device for a longer time.

- It is important that the device stands up on a horizontal and firm subsurface.
- Make sure that the device can not slip away.
- To avoid heat congestion or fires make sure that there is enough space for air circulations.
- Avoid direct solar radiation.
- Do not use the device near highly combustible materials.
- Do not set the device outdoors.
- It is superiorly to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- To set the device in wet or dusty places could cause overheating, short circuits or fires.

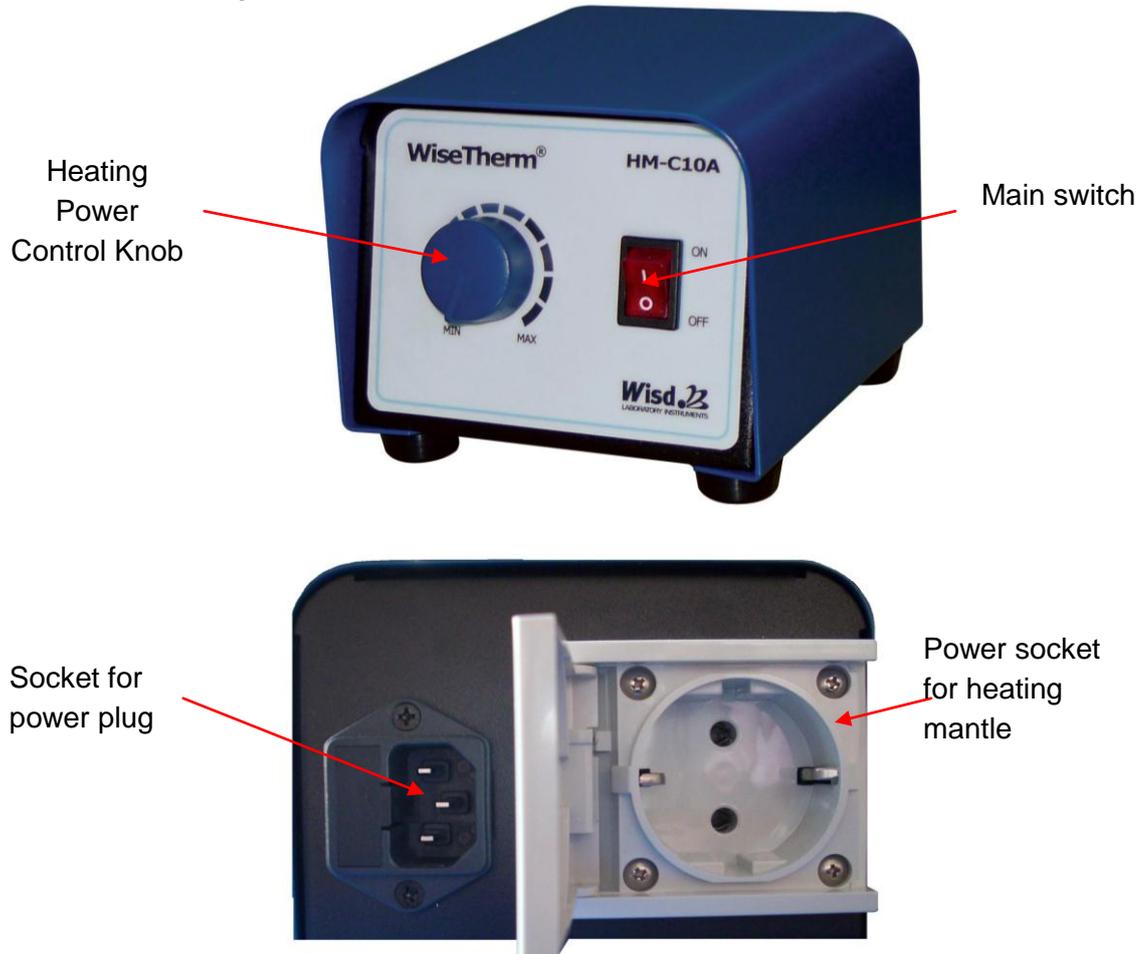
- Do not use this device in an area with a potentially explosive atmosphere.

## 3 Product description

### 3.1 Key features

The **Wisdom WHM-C10A Controller for Heating Mantles** is designed for all Heating Mantles without controller up to 1.200 Watts overall power consumption.

### 3.2 General Survey

### 3.3 Term description

#### Heating Power Control Knob

- Adjustment of heating power

#### Main switch

- On / off switch.

#### Power socket for heating mantle

- Power connection for heating mantle.

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## 5 Positioning the device and commissioning

1. Follow especially the safety instructions marked with following symbol: 
2. Remove the package carefully.
3. If there are any unexpected transport damages, please inform your forwarder and your customer service.
4. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatised.
5. Make sure that you follow all safety instructions.

6. Connect the device with the power outlet.
7. Switch the device on.
8. Set your values and start the process (look at job description).

## 6 Job description

### Temperatur Control

- When you turn the control knob to clockwise, click sounds will on. It means that heater is on for start heating in mantle. If you want more heating, have to turn the control knob to clockwise. On the other hand, if you want less heating, have to turn the control knob to other clockwise.
- Also if you want stop heating, have to turn the control knob to other clockwise until listen click sounds for control knob.

## 7 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device and pull the power plug out of the electrical outlet.
- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- Only use original spare parts.

## 8 Technical data

|                        | WHM-C10A                  |
|------------------------|---------------------------|
| Dimension (mm) (BxTxH) | 160 x 128 x 145           |
| Controller             | Analog Phase Controller   |
| Control Range          | 5 ~ 100% of Rated Voltage |
| Power outlet           | 0 – 1, 2 kW               |
| Weight                 | 1,0 kg                    |
| Others                 | CE Certified              |
| Power plug type        | Hybrid E+F                |
| Power supply           | AC 230 V, 1N~, 50/60 Hz   |

## 9 Certificate

# Certificate

for

**The quality and all features were checked by the manufacturer before the shipment.**

We grant from date of purchase

**two years guaranty.**

This certificate excludes damages by natural disasters or incorrect usages by the costumer.

Please look on your account and complete following table:

|                  |  |
|------------------|--|
| <b>Article</b>   | <b>WHM-C10A Controller for Heating Mantles</b> |
| <b>Type</b>      |  |
| <b>Serialno.</b> |  |
| <b>Date</b>      |  |

**witeg Labortechnik GmbH**  
Am Bildacker 16  
D-97877 Wertheim  
TEL: +49-9342-9301-0  
FAX: +49-9342-9301-77  
Email: info@witeg.de  
[www.witeg.de](http://www.witeg.de)