**WLF**

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# 1 Introduction

Thank you for purchasing our **WLF-420 Digital Laboratory Freezer**. You did a good choice buying a **Wisdom** product.

Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you have to follow the safety instructions of the second chapter.

Please keep this operating manual in a safe place for future reference and hand it out if you consign the device to a third person.

**Witeg** reserves to modify the capsule or make amendments for the improvement of quality and efficiency without prior notice.

## 2 Safety Instructions

### 2.1 Description of symbols

**WARNING**

Information for avoiding injuries or fatal accidents.

**ATTENTION**

Safety instruction to avoid a damage of the product.

**ATTENTION**

Information about technical parameters.

**Environment**

Information about the operating environment.

**WARNING of explosions**

Safety instruction to avoid explosions.

### 2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Connect the device only to the suitable power supply. Please have a look to the identification plate.
- Use the device only in electric circuits with ground fault circuit interrupter.
- Check the power supply cord before you stick it into the electrical outlet.
- Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.

- Do not touch the device with wet hands during the usage because this might cause an electrical shock.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.

- Reparations and amendments should only be made by **witeg** authorized people. Contraventions exclude the warranty.

- Use this device only if you read the whole safety instructions.
- This device should only be used by persons who are familiarized with safety instructions made for laboratories.
- Pay attention that no liquid reaches the controller or inward of the capsule.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective work clothing during the usage.
- All circuits and electrotechnical components are **witeg** owned technology. In fact of this it is necessary that only **witeg** authorized personnel repair the device.
- Pull out the power cable if you don't want to use the device for a longer time.

- It is important that the device stands up on a horizontal and firm subsurface to avoid vibrations.
- If you set the device in a site with vibrations, make sure that the device can not slip away.
- The device could get damaged if the vibrations are too strong.
- To avoid heat congestion or fires make sure that there is enough space for air circulations.
- Avoid direct solar radiation.
- Do not use the device near highly combustible materials.
- Do not set the device outdoors.
- It is superior to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- To set the device in wet or dusty places could cause overcooling, short circuits or fires.

- Do not use this device in areas with potentially explosive atmospheres.

### 3 Product description

#### 3.1 Key features

The **Wisdom WLF-420 Digital Laboratory Freezer** are useful for various kind of laboratory works such like Chemistry, Life Science, Medical Science, Pharmacology, etc. Suitable for Storage of Bio Reagent, Seed, Enzyme and Various Samples.

Convenience and safety in use is considered in developing and manufacturing this product. The device was developed for simplest usage and a very long life.

Key features are as follows:


- Simplest control by Jog-Shuttle knob.
- Fast commissioning.
- Time and temperature showed on illuminated LCD.
- Storage functions for adjusted time and temperature values.
- Turn-off and turn-on delay from one minute up to over 99 hours.
- Signal sound after end of delay time.
- Easy to justify.
- Key lock against inadvertent changing of set temperature and time values.
- Corrosion Resistant 304 Stainless Steel Chamber and Shelves.
- Intelligent Control of Refrigerator for Preventing Overload of Compressor.
- Inner room made o stainless steel.
- Automatic Defrost System.

#### 3.2 General survey

### 3.3 Operating surface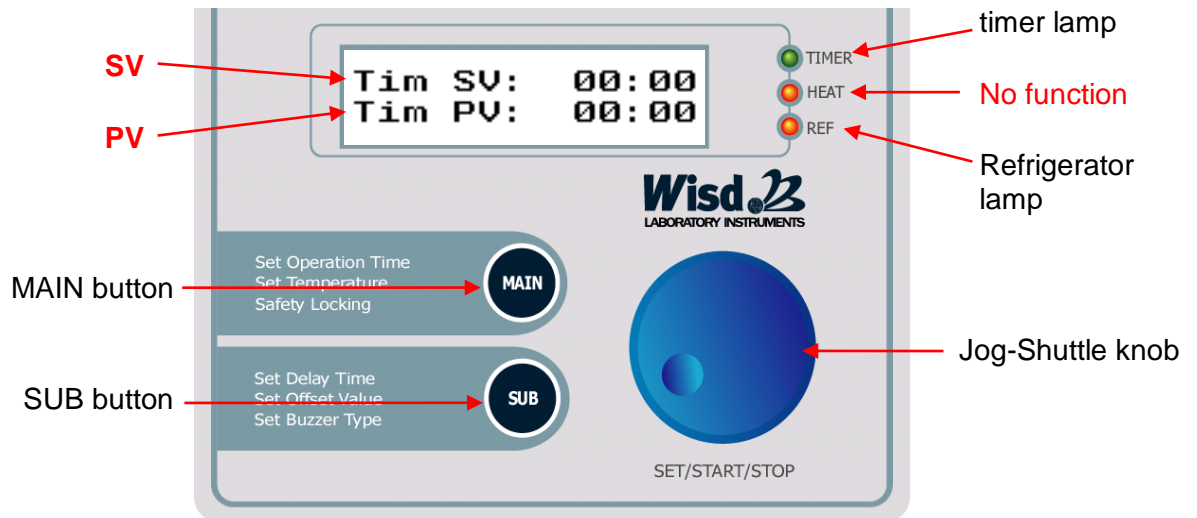

### 3.4 Term description

**SV:** set value (upper line, in key lock menu item no SV)

**PV:** present value, measured by the device (lower line, in key lock menu item also upper line)

#### Main switch

- On / off switch.

#### Timer lamp

- Is blinking in menu items for timer modulation
- Shines continuous in the key lock menu item.

#### Refrigerator lamp

- Shines continuous in the key lock menu item when refrigerator runs.

#### Jog-Shuttle knob

- Change SV by turning.
- Push the knob for start and stop or probate set values.

#### MAIN button

- For changing between the different menu items.

#### SUB button

- For reaching the sub menu Items.

## 4 Positioning the device and commissioning


1. Follow especially the safety instructions marked with following symbol:
2. Remove the package carefully.
3. If there are any unexpected transport damages, please inform your forwarder and your customer service.
4. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatised.
5. Make sure that you follow all safety instructions.
6. Connect the device with the power outlet.
7. Switch the device on.
8. Set your values and start the process (look at job description).

## 5 Job description

### 5.1 Menu structure

The **Wisdom WLF-420 Digital Laboratory Freezer** proffers you the genial combination of MODE and SUB button and the Jog-Shuttle knob. That allows a simple and comfortable usage. The menu structure is well-arranged and pictured as follows.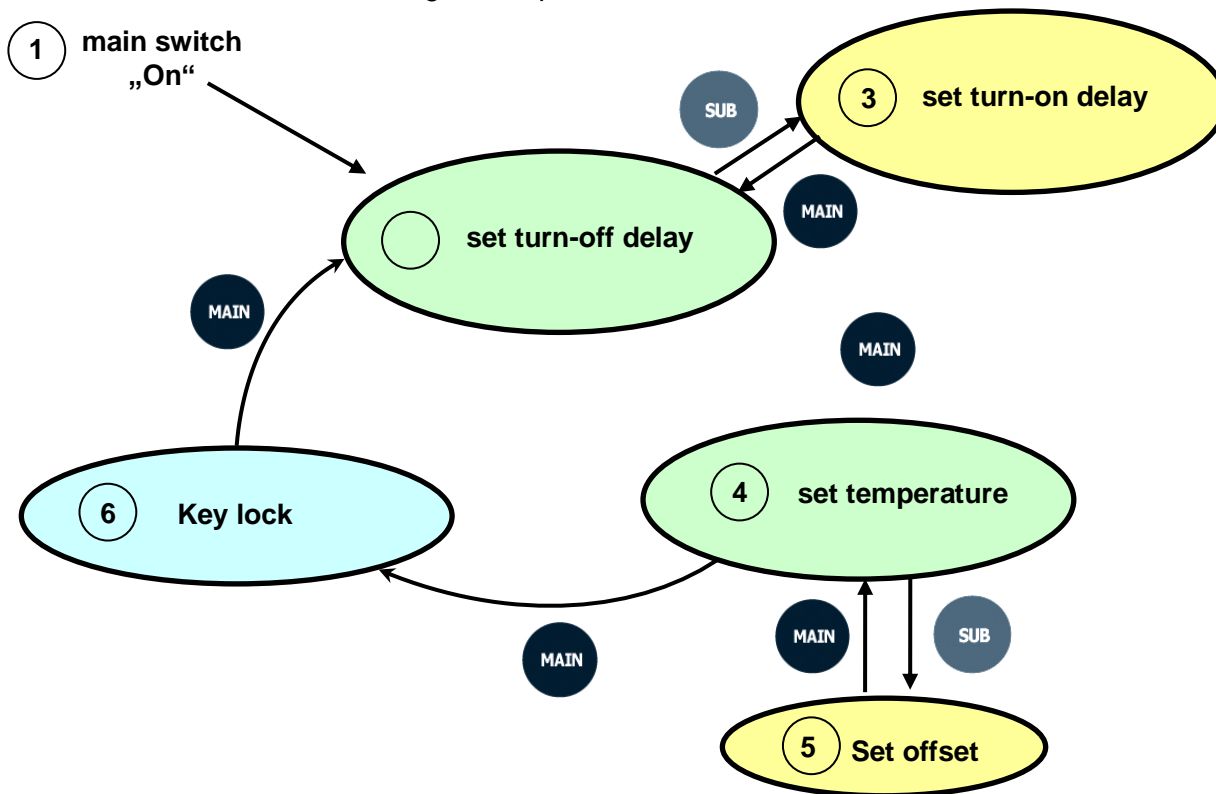

### 5.2 Set turn-off delay

The turn-off delay timer serves as automatically stop function. It turns off the cooling after elapsing of the timeframe you set.

- First menu item after switching the device on is the menu “set turn-off delay”. The timer lamp is blinking.
- The delay time is adjustable from one minute till over 99 hours (00:01 - 99:59) in one minute steps.
- “00:00” stands for endurance run.
- Turn the Jog-Shuttle to choose your desired value and probate it by pushing the Jog-Shuttle. The delay time starts to run by probating the value.
- If you want to change the confirmed value, you have to push the Jog-Shuttle once again.

### 5.3 Set turn-on delay

The turn-on delay timer serves as automatically start function. It turns on the cooling and turn-off delay timer after the timeframe you set is elapsed.

- To reach this sub menu item you have to push the SUB button in the menu “set turn-off delay”
- The delay time is adjustable from one minute till over 99 hours (00:01 - 99:59) in one minute steps.
- Turn the Jog-Shuttle to choose your desired value and probate it by pushing the Jog-Shuttle. The delay time starts to run by probating the value.

- If you want to change the confirmed value, you have to push the Jog-Shuttle once again.
- Turn the Shuttle knob clockwise to increase the time and counter-clockwise to reduce it.
- You only reach this menu item if there is no turn-off delay time running.

#### 5.4 Set temperature

- To reach this menu item you have to push the MAIN button in the menu "set turn-off delay"
- The temperature is adjustable from -35 °C to +10 °C.
- Turn the Jog-Shuttle to choose your desired value and probate it by pushing the Jog-Shuttle. The delay time starts to run by probating the value.
- If you want to change the confirmed value, you have to push the Jog-Shuttle once again.
- Turn the Shuttle knob clockwise to increase the temperature and counter-clockwise to reduce it.
- The refrigerator lamp shines, if the cooling is activated.

#### 5.5 Set Offset

If the temperature measured by you differs from the temperature the device is showing (PV) you have the possibility to compensate the divergence by the Offset. So you can balance differences of about  $\pm 10^{\circ}\text{C}$ . The device will show the same temperature as the thermometer you used.

The Offset is calculated as follows:

**Temperature measured by you – showed Temperature (PV) = Offset**

- To reach this menu you have to push the SUB button in the menu item "set temperature".
- You can recognize this menu at the PV showed in the SV-line.
- If you turn the Jog-Shuttle knob.
- Set the calculated value by turning the Jog-Shuttle. The Offset will be added to the PV immediately.
- Save and confirm the Offset by pushing the MODE button, you will also reach the previous menu.

#### 5.6 Key lock

- To avoid accidental changing of the set values, push the MODE button in the menu item "set temperature".
- In this menu item can you see the actually temperature in the SV-line and the remaining time of turn-off delay in the PV-line. The function of the Jog-Shuttle knob is deactivated.
- Press the MODE button once more to leave the key lock. Than you are in the menu item "set turn-off delay".

#### 5.7 Storage function

The **WLF-420 Digital Laboratory Freezer** includes a storage function for temperature and time values. That is practical if you want to work with the same parameters often. Then you didn't need to set them in every usage.

- For storing values you only need to hold the Jog-Shuttle knob pushed for two seconds. Successful storing is displayed by blinking of the SV for three times.
- Any stored parameters are erasable. Turn the value with the Jug-Shuttle till "0" and then hold the knob pressed for two seconds.
- If you want to overwrite a value it is not necessary to erase the old one. Set the desired value and use the Jog-Shuttle knob for storing.

## 6 Packing contents

quantum	denomination
1	<b>WLF-420 Digital Laboratory Freezer</b>
3	<b>Stainless Steel Shelves</b>
1	<b>operating manual</b>

## 7 Automatic Defrost System

If the unit runs for a long time, there could be generated “frost” inside the chamber owing to even a little humid existing in the chamber. Frost in the chamber could fall the refrigerating performance seriously. So, if the unit runs longer than a week continuously, Automatic Defrost function is executed for 10 minutes once in a week. It is an automatic / periodic operation for removing potential frost in the chamber.

## 8 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device.
- Pull the power plug out of the electrical outlet.
- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- Only use original spare parts.

## 9 Troubleshooting Guide

Situation	Confirmations and Solutions
Power Failure	<ul style="list-style-type: none"> <li>✓ Check the supplied voltage is present.</li> <li>✓ Check the power connection and the power cord for proper connection and integrity.</li> <li>✓ Check if the Switch for Refrigerating is “ON”.</li> <li>✓ Fuse is blown.</li> </ul>
Unit Not Refrigerating	<ul style="list-style-type: none"> <li>✓ Check if the Switch for Refrigerating is “ON”.</li> <li>✓ Check if the set temperature (SV) confirmed with the Jog-Shuttle knob (by turn the Jog-Shuttle knob cannot change the set temperature – see point 5.4).</li> </ul>
Refrigerating but it doesn't fall to the SV temperature	<ul style="list-style-type: none"> <li>✓ Problem with the cooling circuit.</li> </ul>
ERROR 1	<ul style="list-style-type: none"> <li>✓ Temperature Sensor or main controller Error.</li> </ul>

**If other problems arise, or one of the above not solves problems, please contact your official agent or the manufacturer.**

**Reparations and amendments should only made by [witeg](#) authorized people. Contraventions exclude the warranty.**

## 10 Technical data

		WLF-420
<b>Dimension (w×d×h)</b>	<b>Intern (mm)</b>	600 × 700 × 1.000
	<b>Extern (mm)</b>	800 × 800 × 1.700
<b>Capacity (Lit.)</b>		420
<b>Refrigerating Power</b>		750 W
<b>Temp.</b>	<b>Range</b>	-35 °C to +10 °C.
	<b>Sensor</b>	PT100
<b>Controller</b>		Digital Fuzzy Controller with Jog-Shuttle Switch (Turn + Push)
<b>Display</b>		Digital LCD Display with Back-Light Function
<b>Timer</b>		99 hr 59 min, (Delay & Continuous Function Included)
<b>Material</b>	<b>Internal</b>	Stainless Steel (#304)
	<b>External</b>	Powder Coated Steel
<b>Shelves</b>		3ea Stainless Steel Shelves(#304), 16 kg
<b>Refrigerant</b>		CFC-Free (134A) Refrigeration System
<b>Safety Device</b>		Over temp. and over current protector, sensor error detector
<b>Others</b>		Storage function (Temp., Timer), Locking mode (Jog-Shuttle input disabled)
<b>Power Supply</b>		AC 230 V, 1N~, 50/60 Hz

## 11 Certificate

# Certificate

for

**The quality and all features were checked by the manufacturer before the shipment.**

We grant from date of purchase

**two years guaranty.**

This certificate excludes damages by natural disasters or incorrect usages by the costumer.

Please look on your account and complete following table:

<b>Article</b>	<b>WLF-420 Digital Laboratory Freezer</b>
<b>Typ</b>	
<b>Serialno.</b>	
<b>Date</b>	

**witeg Labortechnik GmbH**

Am Bildacker 16  
D-97877 Wertheim  
TEL: +49-9342-9301-0  
FAX: +49-9342-9301-77  
Email: info@witeg.de  
[www.witeg.de](http://www.witeg.de)