WUC-A

Ultraschallreiniger Bedienungsanleitung Ultrasonic Cleaner Operating Manual

gilt für / written for
WUC-A 01H/ 02H/ 03H/ 06H/ 10H/ 22H

witeg Labortechnik GmbH

(English manual at page 7)

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Contents

1 Introduction	8
2 Safety Instructions.....	8
2.1 Description of symbols.....	8
2.2 Instructions	8
3 Product description	10
3.1 Key features	10
3.2 General survey.....	10
4 Packing content	11
5 Positioning the device and commissioning.....	11
6 Job description	11
7 Maintenance and cleaning	11
8 Technical data	12
9 Certificate	13

1 Introduction

Thank you for purchasing our **WUC-A Ultrasonic Cleaner**. You did a good choice buying a **Wisd** product.

Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you have to follow the safety instructions of the second chapter.

Please keep this operating manual in a save place for future reference and hand it out if you consign the device to a third person.

Witeg reserves to modify the capsule or make amendments for the improvement of quality and efficiency without prior notice.

2 Safety Instructions

2.1 Description of symbols

**WARNING**

Information for avoiding injuries or fatal accidents.

**ATTENTION**

Safety instruction to avoid a damage of the product.

**ATTENTION**

Information about technical parameters.

**Environment**

Information about the operating environment.

**WARNING of explosions**

Safety instruction to avoid explosions.

**ATTENTION to high temperatures**

Instructions to avoid injuries caused by high temperatures.

2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Connect the device only to the suitable power supply. Please have a look to the identification plate.

- Use the device only in electric circuits with ground fault circuit interrupter.
- Check the power supply cord before you stick it into the electrical outlet.
- Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.
- Do not touch the device with wet hands during the usage because this might cause an electrical shock.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.
- Switch the device off during when you load the device.
- Avoid injuries: never put your hands or other parts of your body in the working device.
- Do not clean living creatures in this device.

- It is important that the device stands up on a horizontal, stable and firm subsurface to avoid vibrations.
- Make sure that the device can not slip away.
- To avoid heat congestion or fires make sure that there is enough space for air circulations.
- Avoid direct solar radiation.
- Do not use the device near highly combustible materials.
- Do not set the device outdoors.
- It is superiorly to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- To set the device in wet or dusty places could cause overheating, short circuits or fires.

- Do not use this device in area with potentially explosive atmosphere.

- Reparations and amendments should only made by **witeg** authorized people. Contraventions exclude the warranty.

- Do not put heat-sensitive items at the device.
- Warn all attendees if you leave the device unattended.
- Pay attention when you lift up the lid, hot condensation water may drop down or hot steam could come out.
- Do not clean the hot device.

- Use this device only if you read the whole safety instructions.
- This device should only be used by persons who are familiarized with safety instructions made for laboratories.
- Pay attention that no liquid reaches the controller or inward of the capsule.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective work clothing during the usage.
- Pull out the power cable if you didn't want to use the device for a longer time.

3 Product description

3.1 Key features

The **Wisdom WUC-A Ultrasonic Cleaner** is designed for chemical, biological, medical and pharmaceutical applications. The device was developed for simplest usage and a very long life.

Key features are as follows:

- Long life by simply and solid construction.
- 6 different sizes (1,2; 1,8; 3,3; 6; 10; 22 liter).
- Analog timer and temperature adjustment.
- Frequencies: 28 kHz and 40 kHz (from up to 3 liter).
- Temperature for devices up to 2 liter capacity 75 °C (fix).
- Temperature for devices up to 3 liter capacity up to 90 °C adjustable.
- Stainless steel bath.
- Drain valve for devices up to 10 liter capacity.
- Stainless Steel lid inclusive.
- Over temperature protection by safety switch.
- A lot of laboratory devices are mountable.

3.2 General survey

4 Packing content

Article	A01H	A02H	A03H	A06H	A10H	A22H
1	Ultrasonic Cleaner					
1	Stainless steel lid					
1	Power cord					
1	Operating manual					

5 Positioning the device and commissioning

1. Follow especially the safety instructions marked with following symbol:
2. Remove the package and perhaps soiling carefully.
3. If there are any unexpected transport damages, please inform your forwarder and your customer service.
4. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatised.
5. Make sure that you follow all safety instructions.
6. Connect the device with the power outlet.
7. Fill enough operating liquid in.
8. Switch the device on and start the process (look at job description).

6 Job description

1. Position the device according the instructions.
2. Switch on the heating and wait until the water has reached the desired temperature.
3. Put the wasted object into the basket.
4. Putt he basket in the device. Make sure there is a small space between object and bath.
5. Switch on the device with the timer knob.
6. Remove the cleaned object out of the bath after elapsed time.

7 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device.
- Pull the power plug out of the electrical outlet.
- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- Only use original spare parts.
- There is no need to maintenance the device.

8 Technical data

	A01H	A02H	A03H	A06H	A10H	A22H
Capacity	1,2 l	1,8 l	3,3 l	6 l	10 l	22 l
Temperature	65 °C fix	65 °C fix	30-85 °C	30-85 °C	30-85 °C	30-85 °C
Capsule size	177x163x 168	177x163x 211	298x162x 250	359x176x 303	393x265x 370	605x325x 332
Bath size	150x135x 60	150x135x 100	240x140x 100	290x150x 150	290x240x 150	500x300x 150
Timer	-	0-30 min	0-30 min	0-30 min	0-30 min	0-30 min
Drain valve	-	-	-	-	½" NPT	½" NPT
HF-Power (output)	72 W	75 W	124 W	168 W	265 W	379 W
Heating power	86 W	86 W	172 W	172 W	258 W	275 W
Overall power	158 W	161 W	296 W	340 W	523 W	654 W
HF-Frequency	40 kHz					
Weight	1,8 kg	2,5 kg	4 kg	5,3 kg	7 kg	11 kg
Power Supply	AC 230 V, 1N~, 50/60 Hz					

9 Certificate

Certificate

for

The quality and all features were checked by the manufacturer before the shipment.

We grant from date of purchase

two years guaranty.

This certificate excludes damages by natural disasters or incorrect usages by the costumer.

Please look on your account and complete following table:

Article	WUC-A Ultrasonic Cleaner
Typ	
Serialno.	
Date	

witeg Labortechnik GmbH

Am Bildacker 16
D-97877 Wertheim
TEL: +49-9342-9301-0
FAX: +49-9342-9301-77
Email: info@witeg.de
www.witeg.de